

Assembly instructions Tripod

KERN PCD-A03

Version 1.0

08/2011

GB





KERN PCD-A03

Version 1.0 08/2011

Assembly instructions Tripod

Scope of delivery:



Qty.	Components
1	Tripod
4	Screws
1	Cable guidance

Technical data

Tripod height	25 cm
Dimensions	195 x 195 x 250 mm

Assembly:

There are several possibilities to use the tripod PCD-A03:


Variant 1

The tripod is placed on a base foot next to the balance.

Variant 2:


The tripod is assembled to the platform of the balance

- ⇒ Slacken 4 screws on the rear side of the display unit and remove the plastic covers [1]

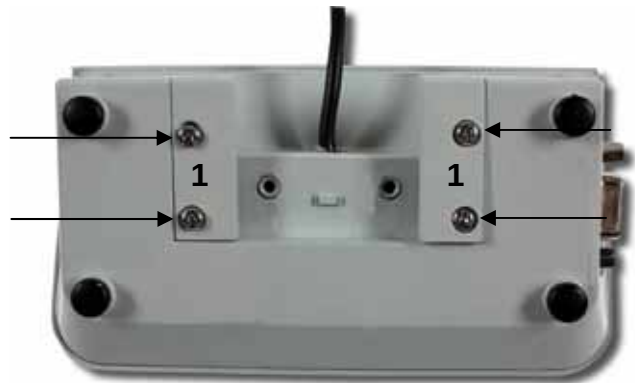


- ⇒ Plug together display unit and tripod acc. to illustration

Make sure that the cable is passed through the designated opening [2].

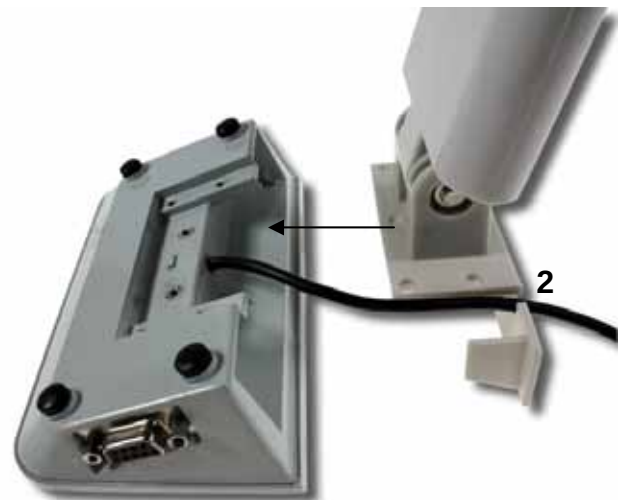


- ⇒ Tighten 4 screws

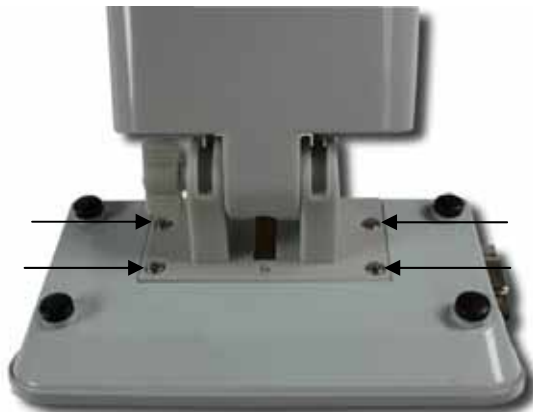

Variant 1

- ⇒ Put the cable on the rear side of the tripod through the cable guidance and fasten cable guidance by 4 screws

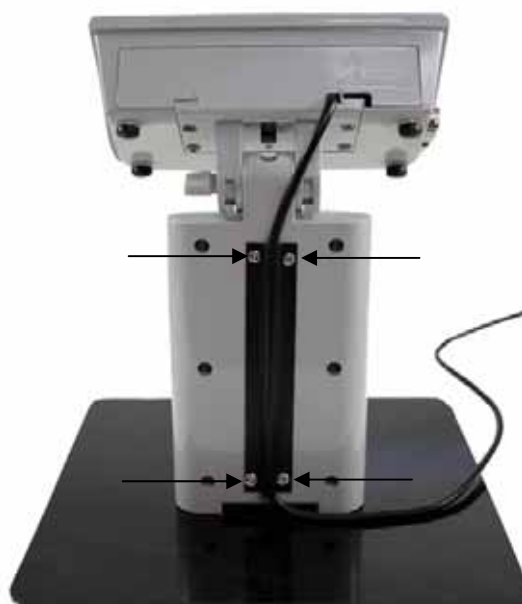

Variant 2:

- ⇒ Unscrew 2 screws at the lower side of the adjustment foot



- ⇒ Push together tripod and support of the platform



⇒ Fasten with screws



⇒ Put the cable on the rear side of the tripod through the cable guidance and fasten cable guidance by 4 screws



