



KERN & Sohn GmbH

Ziegelei 1

D-72336 Balingen

E-Mail: info@kern-sohn.com

Tel: +49-[0]7433- 9933-0

Fax: +49-[0]7433-9933-149

Internet: www.kern-sohn.com

Assembly instructions Tripods

KERN SFE-A01 / A02 / A03

Version 2.0

04/2014

GB



Example SFE-A02

SFE-A01/A02/A03-IA-e-1420



KERN SFE-A01 /A02 / A03

Version 2.0 04/2014

Assembly instructions Tripod

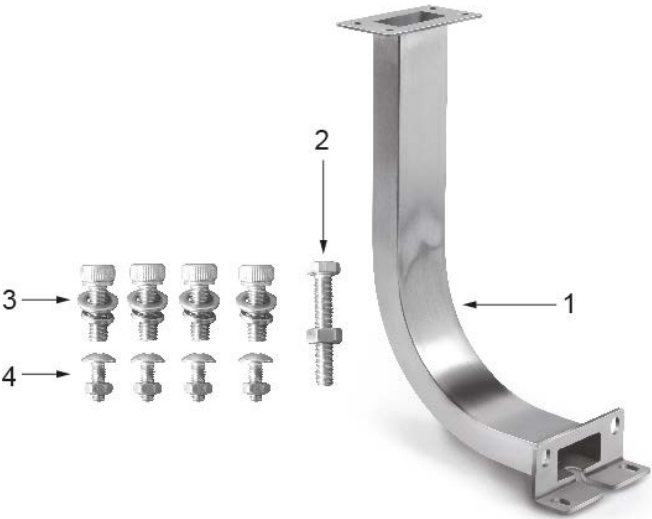
Contents

1	Technical data	2
2	Scope of delivery	3
3	Assembly	3
3.1	Assemble the tripod to the balance	3
3.2	Assemble the display unit to the tripod.....	4

1 Technical data

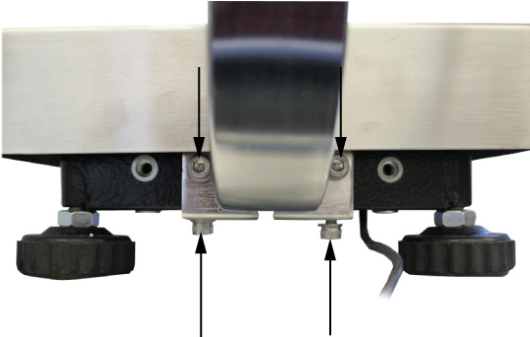
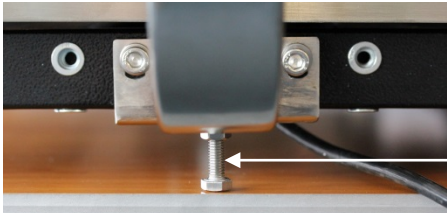
KERN	SFE-A01	SFE-A02	SFE-A03
Tripod height	20 cm	40 cm	60 cm
Net weight	800 g	1600 g	2400 g
For SFE models with weighing plate (mm)	all sizes	400 x 300 x 118 500 x 400 x 137 650 x 500 x 142	500 x 400 x 137 650 x 500 x 142

2 Scope of delivery

	1. Tripod 2. Support screw 3. 4 screws (large) with securing disks and washers for assembly to a balance 4. 4 screws (small) with nuts for assembly of the display unit
---	---

3 Assembly

3.1 Assemble the tripod to the balance

	Attach the tripod to the platform acc. to illustration using the 4 large screws, the securing disks and the washers.
	Screw-out support screw till it is safely fixed.

3.2 Assemble the display unit to the tripod

	Screw tight the display unit to the tripod using the 4 small screws and the nuts according to the illustration. Fix the cable using cable ties in a way that it does not disturb the weighing process.
---	--