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Ziegelei 1

D-72336 Balingen

E-Mail: [info@kern-sohn.com](mailto:info@kern-sohn.com)

Tel: +49-[0]7433- 9933-0

Fax: +49-[0]7433-9933-149

Internet: [www.kern-sohn.com](http://www.kern-sohn.com)

**Betriebsanleitung Temperatur-Kalibrierset  
Operating instruction temperature calibrating set  
Mode d'emploi jeu de calibrage de la température**

**KERN DAB-A01**

Version 1.0

07/2015


Deutsch

English

Français

DAB\_A01-BA-def-1510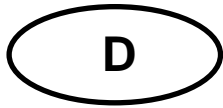

# KERN DAB-A01

Version 1.0 07/2015

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---

---

## **Calibrate / adjust temperature of the heating module**

### **1. Temperature calibration**

We recommend performing occasional tests of the instrument's temperature value, using the optional temperature calibration set **KERN DAB-A01**.

Before you do this, allow the device to cool down for at least 3 hours after the last heating phase.


#### **Preparation:**

- ⇒ Switch off moisture analyser
- ⇒ Install temperature calibration set as per illustration


- ⇒ Turn on moisture analyser


**Set test parameters:**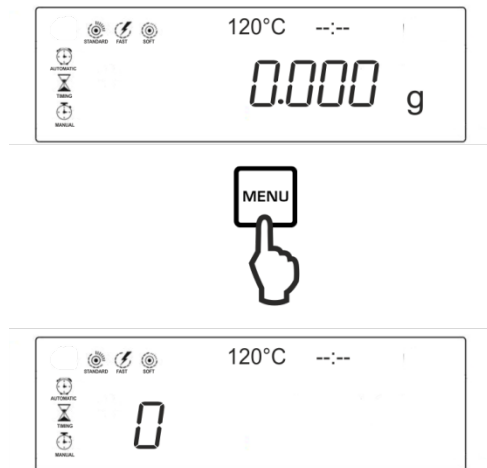

- ⇒ After pressing **MENU** the first ident number <0> will be displayed.
- ⇒ Using arrow keys **↓** **↑** select e.g. ident number <1> and confirm by **ENTER**. The currently set heating profile flashes.

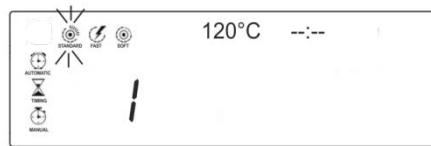
- ⇒ Using arrow keys **↓** **↑** select heating profile <SOFT> and confirm by **ENTER**.

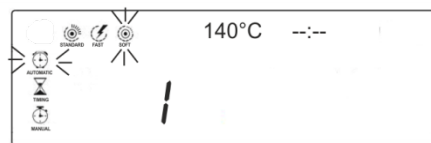
The currently set heating profile and switch-off criterion flashes.

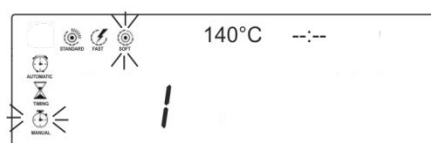
- ⇒ Use the arrow keys **↓** **↑** to select switch-off criterion <MANUAL>.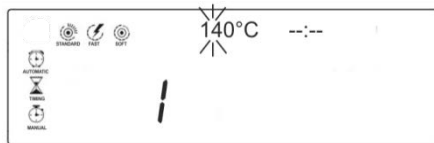


⇒ Confirm by , the currently set drying temperature flashes.

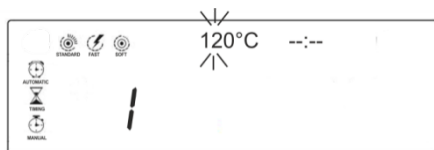⇒ Use the arrow keys   to enter the desired test temperature. For the first or the first two digits select a value and confirm by .

⇒ Using arrow keys   select the desired value for the last digit.

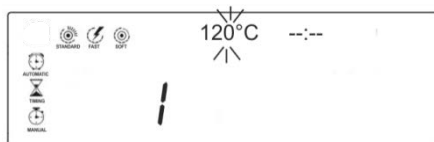⇒ Confirm input by .

### Begin temperature calibration:

⇒ Switch on temperature calibration set.

⇒ Close heating hood and press . The instrument will be automatically heated to the set temperature. The current temperature and the elapsed time will be indicated in the display.

⇒ After approx. 15 minutes finish by . Compare the temperature value of the moisture analyser with that of the temperature calibration set. At a difference of  $\pm 5^{\circ}\text{C}$  we recommend to adjust the temperature, see following chapter.

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
If the admissible deviation is exceeded / not reached during temperature calibration the temperature of the instrument can be adjusted as described below.

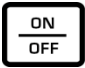1. Switch off instrument.

2. Press , quickly release, then press  and  at the same time.


„8.1x“ will be displayed. If not, separate device from power supply and restart step two.

3. Adjust temperature using arrow keys  . Each time you press the button the temperature will be increased or reduced by 1°C.
4. Confirm input by .

For further information on how to operate the moisture meter please refer to the operating instructions enclosed with each model.

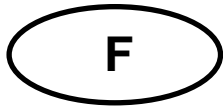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