Bedienungsanleitung User Manual

PC Software PCE-RVI 2/8/10

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English Contents

1	Safety notes	9
2	System requirements	9
3	Installation	9
4	Description of the user interface	10
5	Meaning of the individual icons on the toolbar	11
6	Operation	12
6.1	The first use of the software	12
6.2	Connect to the PCE-RVI 2/8/10	12
6.3	Disconnect from the PCE-RVI 2/8/10	12
6.4	Performing the measurement	13
6.5	Loading and saving series of measurements	13
6.6	Export of series of measurements	13
6.7	List of measurement series	14
6.8	Numerical display	14
6.9	Statistical data	15
6.10	Graphical representation	15

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1 Safety notes

Acquaint yourself with the operation of the measuring device before using it in combination with the PC software. For this purpose, use the instruction manual which comes with your PCE product. All safety notes from the device's manual have to be observed as well when the device is used with this software.

The device and the software may only be used by qualified personnel. Damage or injuries caused by non-observance of the manuals are excluded from our liability and not covered by our warranty.

We do not assume liability for printing errors or any other mistakes in this manual.

We expressly point to our general warranty terms which can be found in our general terms of business.

If you have any questions, please contact PCE Instruments. The contact details can be found at the end of this manual.


2 System requirements

- Operating system Windows 7 or higher
- USB port (2.0 or higher)
- An installed .NET framework 4.0
- A minimum resolution of 800x600 pixels
- Optional: a printer
- Processor with 2 GHz
- 4 GB RAM recommended

3 Installation

Please run the "Setup PCE-RVI 2-8-10.exe" and follow the instructions of the setup.

4 Description of the user interface

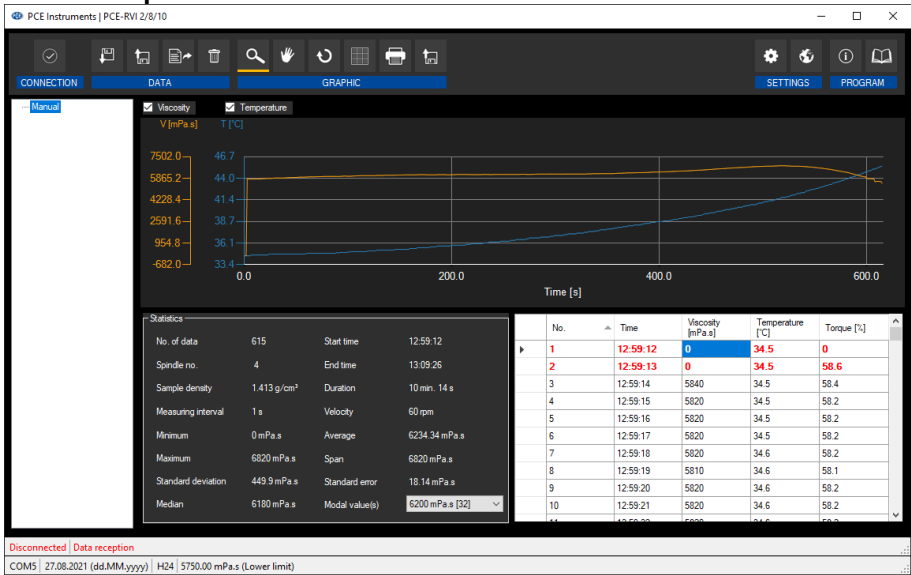The main window consists of several areas:

Below the title bar, there is a toolbar the icons of which are functionally grouped.

Below this toolbar, there is a list of available measurement series in the left part of the window.


The right part of the window is the area for the graphical representation of the readings of a series of measurements.

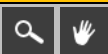Below the graphical view, you can see some statistics and also a tabular listing of the measured values.

At the bottom of the main window there are two status bars. The lower of the two shows the static information: the COM port, the limit values for the selected series of measurements if applicable, and the date and time format.

The upper status bar shows the dynamic data like the connection status.


5 Meaning of the individual icons on the toolbar

Group "Connection"	
	Connect to the PCE-RVI 2/8/10
	Disconnect from the PCE-RVI 2/8/10
Group "Data"	
	Start a measurement
	Complete an ongoing measurement
	Load readings from a file
	Save the readings to a file
	Export readings (CSV format)
	Delete currently selected series of measurements
Group "Graph"	
	Enlarge the graph area ("zooming") or move the enlarged graph
	Restore the original graph
	Change the background and representation of the graph
	Print the currently visible graph
	Save the currently visible graph
Group "Settings"	
	Open settings dialogue
	Select one of the languages supported by the programme
Group "Programme"	
	Display an information dialogue
	Exit the programme

6 Operation

6.1 The first use of the software

Before the PCE-RVI 2/8/10 can work with the software, the assigned COM port must be set in the software once. It can be set via the "settings dialogue" .

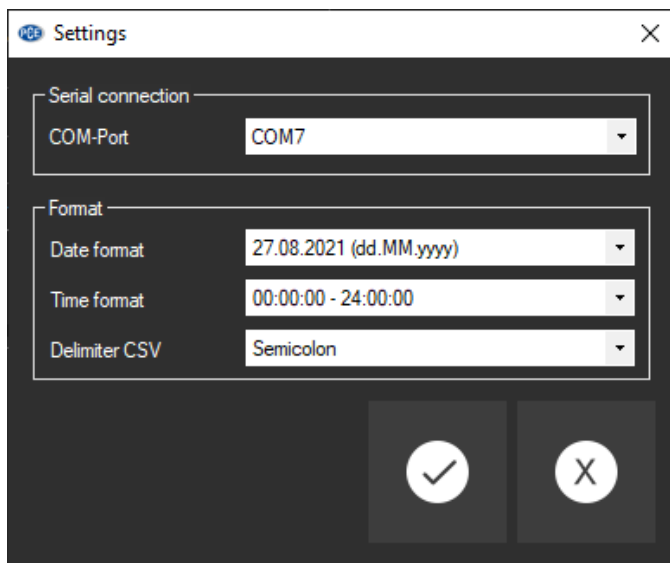In addition to the connection data, further settings for the date and time format and the delimiter for the export of the series of measurements can be made here.

6.2 Connect to the PCE-RVI 2/8/10

After the desired settings have been made and the settings dialogue has been closed with a click on the "Apply" button, the connection to the device can be established by clicking on the  icon.

If the connection could be successfully established, the term "Connected" is displayed in **green** on the status bar for the dynamic data.

6.3 Disconnect from the PCE-RVI 2/8/10

By clicking on the  icon, an active connection to the meter can be terminated.

If you exit the software while the connection is active, the connection will also be terminated.

When the software is not connected, "Disconnected" will be displayed in **red** in the status bar for the dynamic data.

6.4 Performing the measurement

After clicking on the  icon, a new dialogue will be displayed where the name of the new series of measurements can be determined. After confirming via the “Apply” button, the measurement is started.

6.5 Loading and saving series of measurements

To save a series of measurements, it must first be selected in the overview of the series of measurements currently available in the software (left-hand side of the programme).

Now click on the  button to name the file and determine a folder for this series of measurements and to save the name and folder.

A series of measurements can be loaded by clicking on the  icon and selecting a folder and determining a file name.

Alternatively, you can “drag & drop” the file from the Windows Explorer to the list of series of measurements in the software.


6.6 Exporting series of measurements

Using the  icon in the toolbar, the complete content of a single series of measurements can be exported in CSV format.

If desired, the delimiter can be changed with the help of the settings dialogue .


6.7 List of measurement series

In the left part of the main window, there is an overview of all series of measurements currently available in the software.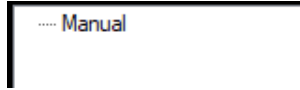

Depending on what series of measurements has been selected, the graph and also the statistical data are updated.

A right click allows you to set limit values for the viscosity of this series of measurements, to rename or to delete a series of measurements. A series of measurements can also be renamed by pressing the “F2” key or deleted by pressing the “Del” key or clicking on the  icon.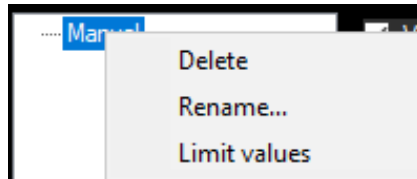

6.8 Numerical display

	No.	Time	Viscosity [mPa.s]	Temperature [°C]	Torque [%]
▶	1	12:59:12	0	34.5	0
	2	12:59:13	0	34.5	58.6
	3	12:59:14	5840	34.5	58.4
	4	12:59:15	5820	34.5	58.2
	5	12:59:16	5820	34.5	58.2
	6	12:59:17	5820	34.5	58.2
	7	12:59:18	5820	34.6	58.2
	8	12:59:19	5810	34.6	58.1
	9	12:59:20	5820	34.6	58.2
	10	12:59:21	5820	34.6	58.2

In the lower part of the main window, all the measured values are tabulated.


6.9 Statistical data

Statistics			
No. of data	615	Start time	12:59:12
Spindle no.	4	End time	13:09:26
Sample density	1.413 g/cm ³	Duration	10 min. 14 s
Measuring interval	1 s	Velocity	60 rpm
Minimum	0 mPa.s	Average	6234.34 mPa.s
Maximum	6820 mPa.s	Span	6820 mPa.s
Standard deviation	449.9 mPa.s	Standard error	18.14 mPa.s
Median	6180 mPa.s	Modal value(s)	6200 mPa.s [32]

To the left of the chart, you can see the statistics of the currently selected measurement.

6.10 Graphical representation

The graph shows the course of the temperature and the viscosity. The x axis shows the time in seconds.

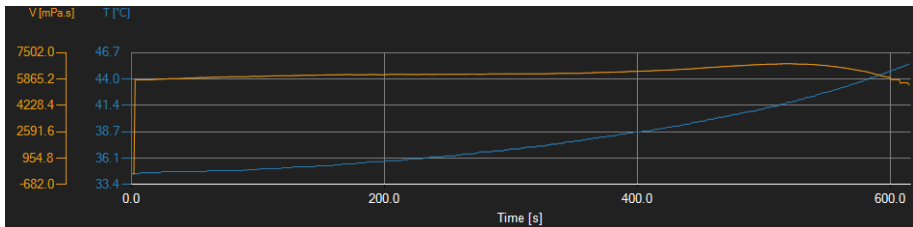Information window

Moving the mouse cursor to a point in the displayed line makes a small information window with the data (time and unit) of the currently selected measurement value pop up after a short time.

Viscosity: 6170 [mPa.s]
 Temperature: 36.3 [°C]
 Time: 9/11/2020 1:03:27 PM

Enlargement/zooming

A freely selectable part of the displayed graph can be enlarged.

For this purpose, the corresponding icon in the toolbar ("Enlarge the graph area ("Zooming") or move the enlarged graph) must be a magnifying glass .

Then, a rectangle can be drawn over a part of the graph by holding the mouse button down. When the mouse button is released, the selected area appears as a new graph.

As soon as at least one enlargement has been made, it is possible to switch from enlargement mode to shift mode by clicking the icon ("Enlarge the graph area ("Zooming") or move the enlarged graph) with the magnifying glass icon.

Shift mode is represented by the  icon.

If the mouse is now placed over the graph area and then the left mouse button is pressed, the depicted section can be moved by holding the mouse button down.

Another click on the  icon changes back to enlargement mode, which is recognizable by the  icon.

Restore complete view

The original graph can be restored at any time by clicking on the  icon next to the magnifying glass or hand.

Activating the background

The background of the graph and its representation can be changed via the  icon to the right. A click on the icon works like a switch:

A single click makes the division of the background finer and adds some more dots to the graph. A further click on the icon changes back to standard view.

Printing the graph

The currently displayed graph can also be printed.

A click on the  icon opens the "Print" dialogue.

Saving the graph

To save the currently displayed graph, just click on the  icon. These formats are possible for the graph: BMP, JPEG, PNG, GIF, TIFF and VMF.

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User manuals in various languages (français, italiano, español, português, nederlands, türk, polski, русский, 中文) can be found by using our product search on: www.pce-instruments.com

Specifications are subject to change without notice.

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PCE Americas Inc.
PCE Teknik Cihazları Ltd.Şti.