

USER MANUAL

SPECTROPHOTOMETER

PCE-CHM 700

ENGLISH



User manuals in various languages (français, italiano, español, português, nederlands, türk, polski) can be found via our product search on:
www.pce-instruments.com

PRODUCT INTRODUCTION

The instrument is a multifunctional color haze meter. It adopts full spectrum LED light source and spectral sensor design, which is used for the measurement of haze, transmittance and color of various diffuse transmittance materials such as milky white, matte, frosted surface and various transparent and semi-transparent regular transmittance materials, and it can also display the spectral transmittance curve. The instrument is equipped with 7-inch color screen and capacitive touch screen, which provides excellent operation experience.

Standards for the product

JJF 1303-2011 Calibration Specification for Hazemeter

GBT 2410-2008 Determination of the luminous transmittance and haze of transparent plastics

GBT 36142-2018 Test methods of color and color difference for architectural glass

JJG 595-2002 Colorimeters and Color Difference Meters

GB/T 3978-2008 Standard illuminants and geometric conditions

GBT 7921-2008 Uniform color space and color difference formula

GB 3143-1982 Colour determination method of liquid chemicals (Hazen unit--platinum-cobalt scale)

JJF 1947-2021 Calibration Specification for Platinum-Cobalt Colorimeter

JJG 880-2006 Verification Regulation of Turbidimeters

PARAMETERS

Illumination geometry	D/0°
Illumination light source	Full spectrum LED light source
Spectral Range	400 ... 700nm
Spectral Interval	10nm
Measuring aperture	5mm/14mm
Measurement conditions	Light source D65, field of view 10°; Light source A, field of view 2°; Light source C, field of view 2°
Measuring time	About 3s
Measuring parameter	Haze, Transmittance, CIE Lab, LCh, Yxy, Spectral transmittance, Clarity, Turbidity, Pt-Co(Hazen)
Transmittance resolution	0.01%
Transmittance accuracy	Better than ±1%
Transmittance repeatability	≤0.03 (The standard deviation of 30 times measurements of a haze standard with haze measurement of approximately 30 at 3s intervals after calibration)

Haze/Clarity resolution	0.01%
Haze/Clarity accuracy	Better than $\pm 2\%$
Haze/Clarity repeatability	≤ 0.03 (The standard deviation of 30 times measurements of a haze standard with haze measurement of approximately 30 at 3s intervals after calibration)
Color resolution	0.01
Color inter-instrument agreement	$\Delta E^*_{ab} \leq 0.4$ (Measurement of a haze standard with haze measurement of approximately 30)
Color repeatability	Standard deviation ΔE^*_{ab} is within 0.03 (Average of 30 measurements of 30 times of a haze standard at an interval of 3s after calibration)
Turbidity resolution	0.1NTU
Turbidity range	0 ... 1000NTU
Turbidity accuracy	$\leq \pm(5\%H+1\text{NTU})$ H is standard
Pt-Co resolution	0.1
Pt-Co range	0 ... 1000
Pt-Co accuracy	$\leq \pm(5\%H+1)$ H is standard
Power supply	100 ... 277V 0.4A 50/60HZ
Display	7-inch 1024*600 dot matrix IPS color screen
Language	Simplified Chinese, English
Data transmission	USB (Type-C)
Operating temperature range	0 ... 45°C, 0 ... 85%RH (no condensation)
Storage temperature range	-25 ... 55°C, 0 ... 85%RH (no condensation)
Size	21.1cm×26.4cm×36.0cm(LWH)
Weight	5.3kg

FEATURES

- » The instrument is fashioned with a full-spectrum LED light source and spectral sensor, and it is used for haze, transmittance, color difference multiple parameters measurement, and can display the spectral transmittance curve.
- » Adopt a 7-inch capacity touch panel with human-machine interaction interface for convenient operation.
- » The instrument has a built-in compensation optical path with stable values, allowing long-period calibration.
- » QC testing function: the threshold can be set.
- » Can store 1000 measurement records with large storage space.
- » A powerful function of color difference analysis as well as data import and export with PC software

OPERATION

Power on/off

Power on: Plug in the power supply, short press the right side button to turn on the instrument.

Power off: Long press the right side button to turn off or press "Power off" in the menu bar.

Calibration

After entering „calibration“ interface, you can perform calibration operation as the calibration animation, or skip the calibration. It is recommended to calibrate after a long time of no usage.

For failed calibration, the possible reasons are as below:

- » The dust cover has not been removed or the test holes are covered with samples;
- » Tested in a bright light environment;
- » The attenuation of the light source leads to failure of normal use and it needs to be returned to the factory for inspection and repair.

Measurement

The default measurement parameter of the instrument is „Haze/Transmittance“. Users can click the „Param“ at the top of the screen or the „Param Setting“ in menu bar to enter the setting interface to select the parameters to be measured (H/T, H, T, CIE_Lab, CIE_LCh, Yxy, Spectra, Clarity, Turbidity, Pt-Co). Entering the measurement interface, you can click the „ Compare“ on the top of the screen to switch the compare measurement mode.

Compare off

In the measurement interface, click on „Scan“ or short press the button on the right side of the instrument to measure. The following steps can be performed after completing measurement:

- » Save - To save the comparison result, press „Save“ on the upper right corner. You can choose default or name the record manually.
- » History - If users need to view the saved results, click „History“ to view the records, and records can be paged, deleted, renamed and searched.

Compare on

In the measurement interface, press the „Scan Standard“ button to set current value as standard. Press the „Scan Sample“ button to use current value as sample. If short press the button on the right side, another measurement will also be taken and the current value will replace the previous one. When measuring or adding a standard and measuring a sample, the difference or color difference between the two measured values will be calculated with the following results:

- » **The difference value** - The difference Δ is the sample data minus the standard data;
- » **Color Bias** - When the measurement parameter is selected as Lab, the color bias will be judged by ΔL , Δa , and Δb ;
- » **ΔE value** - When the measurement parameter is selected as Lab, LCh, or Yxy, ΔE is calculated using the ΔE^*_{ab} color difference formula;
- » **QC test** - Suitable for difference inspection in production process. If ΔE is higher than the set threshold, it will read “NG” and shown in red; if ΔE is less than or equal to the set threshold, it will read “Pass” and display in green.

The following steps can be performed after completing compare measurement:

- » Save - To save the comparison result, press „Save“ on the upper right corner. You can choose default or name the record manually.
- » History - If users need to view the saved results, click „History“ to view the records, and records can be paged, deleted, renamed and searched.
- » Save standard - After the user has finished measuring the standard value, if need to save it for later calling, click the „ Save standard“ to save it, and the saved value can be queried in the add standard interface.
- » Add standard - After clicking „Add standard“, the interface will jump to „Add standard“, click the value you want to add, and the selected value will be used as the standard, and will automatically return to the measurement interface to compare with the current sample or the upcoming sample.
- » Input standard - The user can manually input the Haze, Transmittance, Lab, LCh, Yxy value as standard.
- » Tolerance Setting - Setting the threshold for QC inspection.

Menu bar

In the measurement interface, click the icon on the upper left corner to pop up the menu bar, with the following options: Language, Calibration, Standard, Param Settings, Device info, Reset to defaults and Power off.

Calibration

Can enter the calibration interface to perform calibration operation of the instrument.

Language

Support Simplified Chinese and English.

Standard

Selection of ASTM/ISO standards, switching of measurement aperture 14mm/5mm, and setting of light source and observer's field of view.

Param Settings

Select the parameter desired to be measured and displayed. Shortcut keys for the same functions are available at the top of the measurement interface.

Device Info

„Device Info“ displays the basic information of the instrument, including: Device Model, Firmware Version, and Device Status.

Reset to defaults

To restore parameters to the factory settings and clear the data in the measurement interface.

Power Off

Press „Power off“ to power off the instrument.

PC SOFTWARE

You can use the PC software to connect the computer via USB. The software has functions of compare measurement, reading compare record, export compare data to Excel, qualified number, unqualified number, total number statistics, report generation and printing, etc. For specific operation, please refer to the „Color Haze Meter Software Operation Manual“.

PRECAUTION

- » When the colorimeter has not been used for a long time, it is recommended to perform calibration before using.
- » Please ensure that the sample is evenly colored with a flat and clean surface, otherwise it will affect the measurement accuracy.
- » Avoid liquids entering the instrument through the measurement aperture, as this will damage the instrument and affect measurement accuracy and operational safety.
- » When the instrument is not in use, put on the dust cover to prevent dust from entering or prolonged humidity, which may affect the measurement accuracy.
- » It's recommended to calibrate the instrument once each year. And we offer calibration service.
- » Due to color difference in the LCD display, the color displayed on the instrument screen is for reference only.

DISPOSAL

For the disposal of batteries in the EU, the (EU) 2023/1542 directive of the European Parliament applies. Due to the contained pollutants, batteries must not be disposed of as household waste. They must be given to collection points designed for that purpose. In order to comply with the EU directive 2012/19/EU we take our devices back. We either re-use them or give them to a recycling company which disposes of the devices in line with law. For countries outside the EU, batteries and devices should be disposed of in accordance with your local waste regulations. If you have any questions, please contact PCE Instruments.

PCE INSTRUMENTS CONTACT INFORMATION

Germany

PCE Deutschland GmbH
Im Langel 26
D-59872 Meschede
Deutschland
Tel.: +49 (0) 2903 976 99 0
Fax: +49 (0) 2903 976 99 29
info@pce-instruments.com
www.pce-instruments.com/deutsch

France

PCE Instruments France EURL
2, rue Georges Kuhnmmunch
67250 Soultz-sous-Forêts
France
Tel.: +33 (0) 972 35 37 17
Fax: +33 (0) 972 35 37 18
info@pce-france.fr
www.pce-instruments.com/french

Spain

PCE Ibérica S.L.
Calle Mula, 8
02500 Tobarra (Albacete)
España
Tel.: +34 967 543 548
info@pce-iberica.es
www.pce-instruments.com/espanol

United Kingdom

PCE Instruments UK Ltd
Trafford House
Chester Rd, Old Trafford
Manchester M32 0RS
United Kingdom
Tel: +44 (0) 161 464902 0
Fax: +44 (0) 161 464902 9
info@pce-instruments.co.uk
www.pce-instruments.com/english

Italy

PCE Italia s.r.l.
Via Pesciatina 878 / B-Interno 6
55010 Loc. Gragnano
Capannori (Lucca)
Italia
Telefono: +39 0583 975 114
Fax: +39 0583 974 824
info@pce-italia.it
www.pce-instruments.com/italiano

Turkey

PCE Teknik Cihazları Ltd.Şti.
Halkalı Merkez Mah.
Pehlivan Sok. No.6/C
34303 Küçükçekmece - İstanbul
Türkiye
Tel: 0212 471 11 47
Faks: 0212 705 53 93
info@pce-cihazlari.com.tr
www.pce-instruments.com/turkish

The Netherlands

PCE Brookhuis B.V.
Twentepoort West 17
7609 RD Almelo
Nederland
Telefoon: +31 (0)53 737 01 92
info@pcebenelux.nl
www.pce-instruments.com/dutch

United States of America

PCE Americas Inc.
1201 Jupiter Park Drive, Suite 8
Jupiter / Palm Beach
33458 FL
USA
Tel: +1 (561) 320-9162
Fax: +1 (561) 320-9176
info@pce-americas.com
www.pce-instruments.com/us

Denmark

PCE Instruments Denmark ApS
Birk Centerpark 40
7400 Herning
Denmark
Tel.: +45 70 30 53 08
kontakt@pce-instruments.com
www.pce-instruments.com/dansk