


Leica A60 S – Assembly Sheet

This assembly sheet does **not** replace the complete manual on provided CD (i.e. LED illumination installation).


Note: For safety reasons, it is highly recommend that **two people** assemble the column to the base plate. One person should hold the base plate securely on a non-skid surface, while the second person attaches the column to the base.

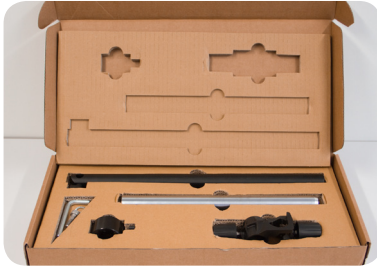

One box contains horizontal arm, cross joint, safety clamp & tools




One box contains the base plate.

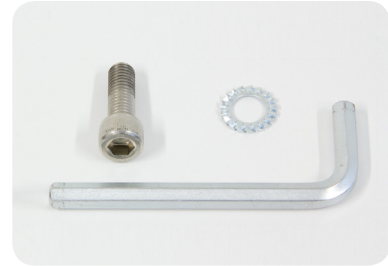


Allen key, washer and allen screw



Take out parts from boxes. Place base plate on non skidding surface.

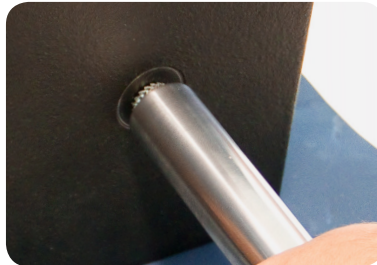

Tighten allen screw connecting column to base. **Attention:** washer must be placed between base and column.

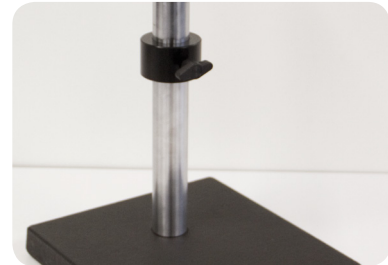

Mount the safety clamp to the column.



Insert horizontal arm into the cross joint observing the flattening of the arm and orientation of joint according to image.



Place the cross joint on top of the safety clamp on the column and lock arm with referring knob.



Remove allen screw and washer from the Leica A60 instrument as shown.



Mount the Leica A60 to the horizontal arm from below...



... and immediately secure the instrument with previously removed washer and screw.



If necessary connect base plate to grounding cable with suitable plug.

Leica A60 F – Assembly Sheet

This assembly sheet does **not** replace the complete manual on provided CD (i.e. LED illumination installation).




Box contains flex arm and table clamp.



Take out table clamp from box.



Pick one of three positions of the clamp according to table thickness.

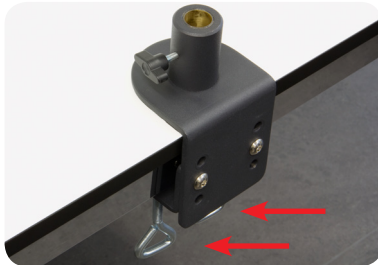

Choose position on table edge and tighten well the two screws.



Insert arm into table clamp and tighten screw.



Open the two screws on the focus arm of the Leica A60...



...and bring pin into horizontal position. Tighten screws again.



Attach Leica A60 to flex arm and tighten knob to secure the instrument.



Your instrument is now ready for usage!