

Digital Eyepiece Quick Start Guide

EN

Thank you for purchasing this product – **Digital Eyepiece**.

Please read these instructions carefully.

The Digital Eyepiece you purchased is provided with CD Instruction Manual.

For detailed information on all features and warranty provision of this Digital Eyepiece, please refer to the CD. PDF reader is required to view the User Manual.

Preparations


Connect Digital Eyepiece to computer USB interface

Plug the USB Digital eyepiece into computer USB interface, the driver will install automatically by the operating system.

Insert the CD to install the software

- The supplied CD and installation software is available for Microsoft Windows System only.
1. Put your CD into the CD-ROM driver.
 2. Run the Digital micro Eyepiece program and follow the prompt information to finish installation

How to fit the digital eyepiece to microscopes

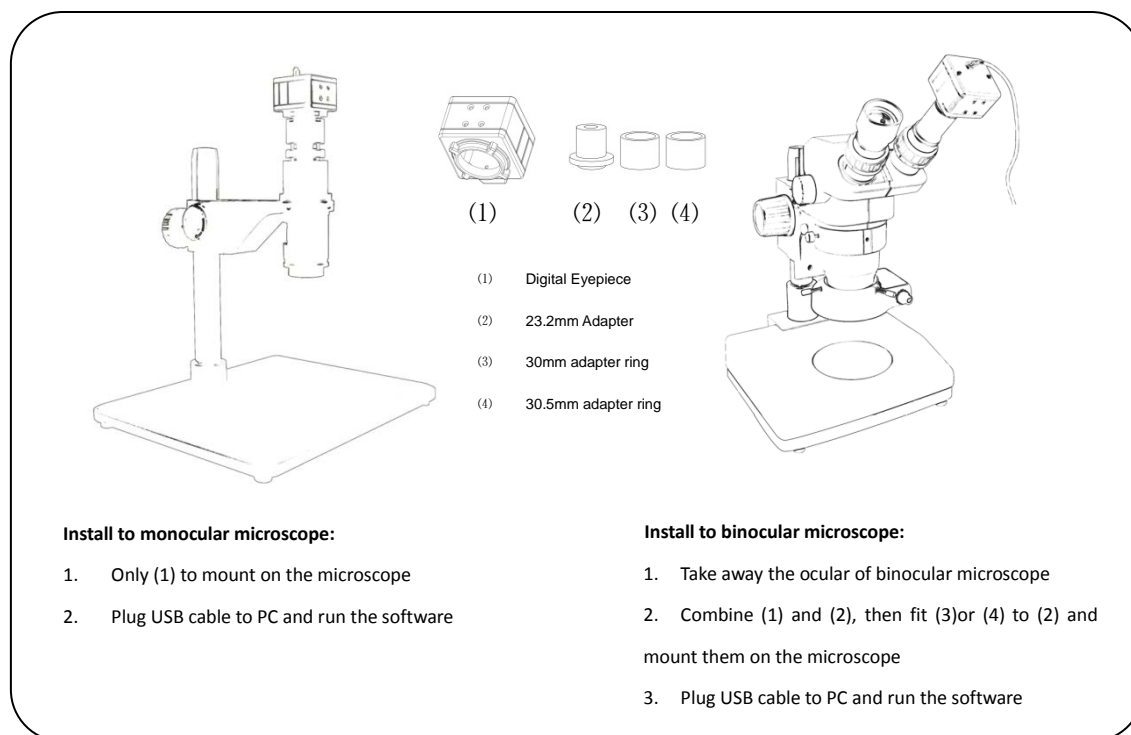

EN

Safety Precautions

- To avoid danger or damage incurred to the sensor, do not touch it directly with your fingers.
- To avoid failure or electric hazard and so on, do not disassemble or modify the internal structure of the device.
- Do not plug in or unplug the USB connectors when the hands are wet
- Do not use alcohol and other organic solvents to clean the device.
- If the sensor is dirty or damp, you should better use the dry and clean cotton swab tissue to clean it. To avoid scratches on the surface, do not touch the sensor with your fingers and wipe the sensor lightly.
- The products are not specifically designed for an outdoor use. Do not expose it to outdoor environment without any protection, because the excessive temperature and humidity will damage the sensor
- Please store and use in the following environment: operation: temp= -20~60 °C HR=10%~85% ;

storage: temp= -40 ~ +8°C HR=5~95%

- If there is any foreign matter / water or liquid enter into the device by accident, disconnect the USB cable immediately. Please send it to the maintenance center and do not use the hair dryer to dry it by yourself.
- To prevent being tripped over or dropping the digital eyepiece, please put away the USB cable of digital eyepiece.
- To avoid electric shock by accident, please turn off and power off the digital eyepiece, before you move the PC.

Warning

1. Do not use the digital eyepiece near the flammable or explosive objects.
2. Keep the young children and infants away from the digital eyepiece.
Always use and store the digital eyepiece out of the reach of young children and infants to prevent the following dangerous situations which might cause serious injury:
 - 1) Becoming entangled in the digital eyepiece strap, causing strangulation.
 - 2) Accidentally being injured by the moving parts of the digital eyepiece

Trademarks

1. Microsoft Windows are registered trademarks of Microsoft Corporation
2. All other company and product names are registered trademarks and/or trademarks of their respective owners.