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Ziegelei 1

D-72336 Balingen

E-Mail: [info@kern-sohn.com](mailto:info@kern-sohn.com)

Tel: +49-[0]7433- 9933-0

Fax: +49-[0]7433-9933-149

Internet: [www.kern-sohn.com](http://www.kern-sohn.com)

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## 11 Preload, Deadload and Overload settings

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

















## 12 Anhang

| ZONE | REV. | DESCRIPTION                 | DATE      | APPROVED |
|------|------|-----------------------------|-----------|----------|
| 1    | B    | REDRAWN FROM 021450321_REVA | 9/19/2023 | ZHUANG   |
|      |      |                             |           |          |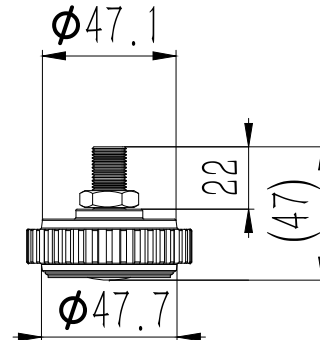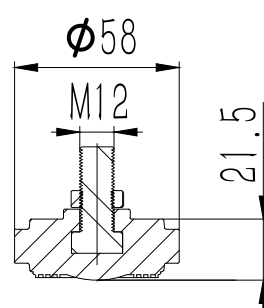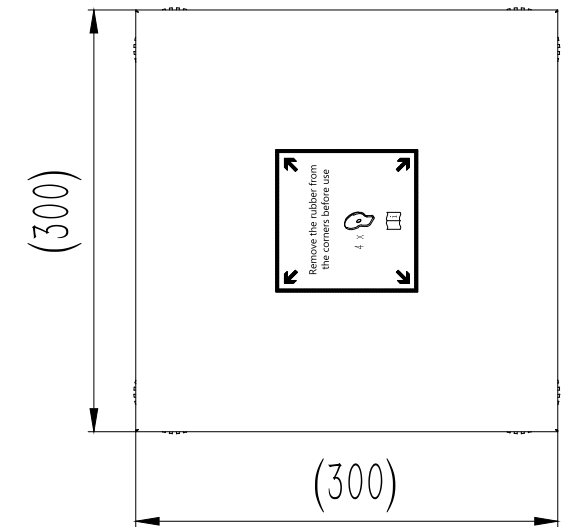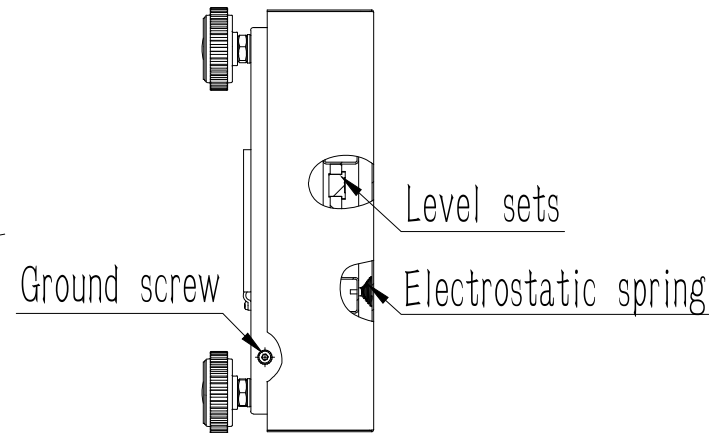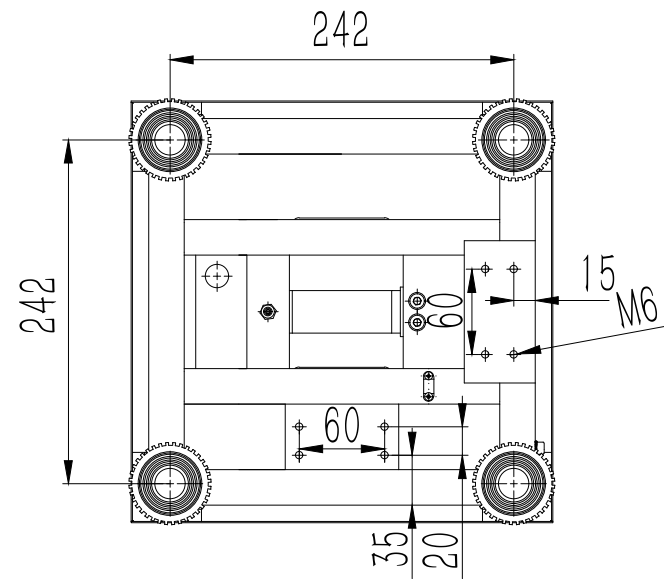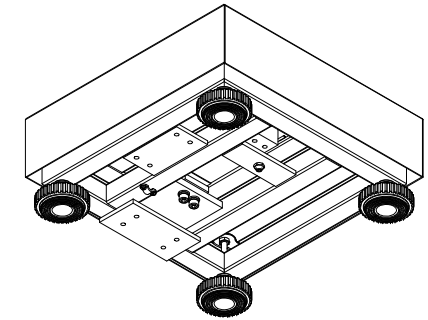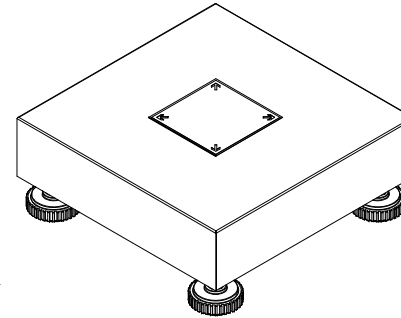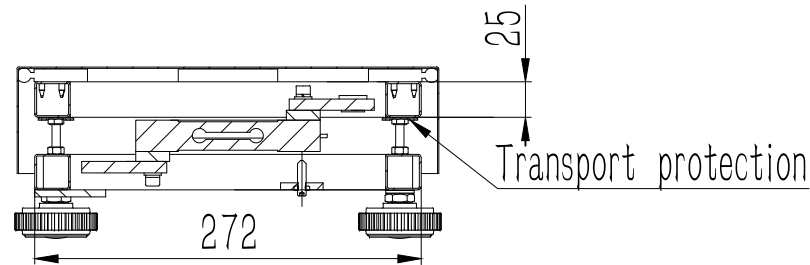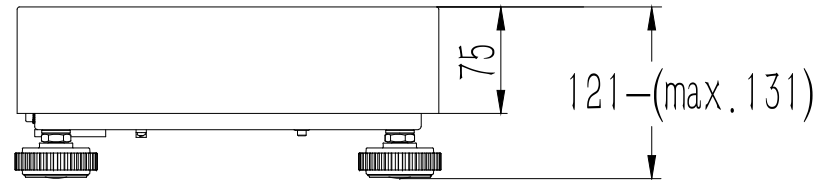









|                                                                                       |     |                      |                                                              |                                          |                           |
|---------------------------------------------------------------------------------------|-----|----------------------|--------------------------------------------------------------|------------------------------------------|---------------------------|
| TOLERANCE UNLESS OTHERWISE SPECIFIED<br>DEC ANGLES<br>±0.1 FOR .X<br>±0.05 FOR .XX    |     | ±0.5                 |                                                              | <b><i>KERN® KERN &amp; Sohn GmbH</i></b> |                           |
|  |     | 3RD ANGLE PROJECTION |                                                              | TITLE: TKBP 6V20LM-A                     |                           |
| DRAWN                                                                                 |     | ZHUANG               |                                                              |                                          |                           |
| CHECKED                                                                               | LIN | MODEL NO.            | B                                                            | PART NO.                                 | 021450321                 |
| APPROVED                                                                              | LIN | MATERIAL             | Weighing plate, stainless steel,<br>Platform, steel, coated, | FINISH                                   | Powder coat "pantone 445" |
| UNIT                                                                                  | mm  | SCALE                |                                                              | DO NOT SCALE DRAWING                     | SHEET OF                  |









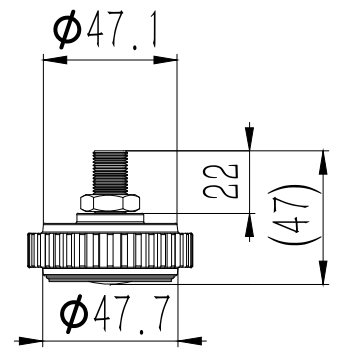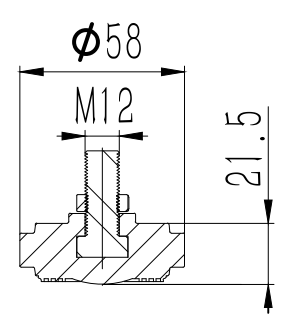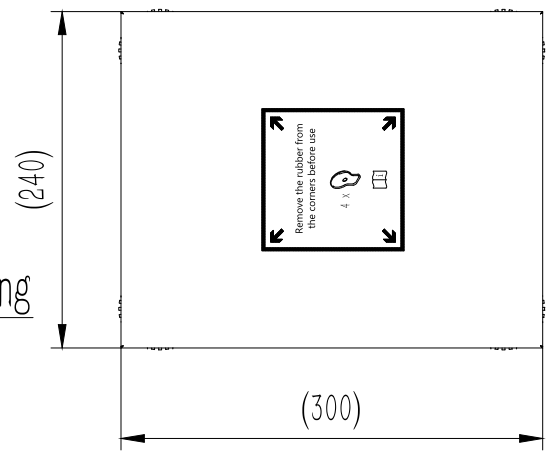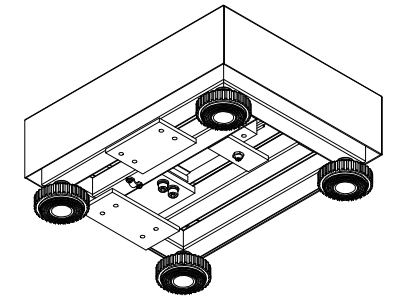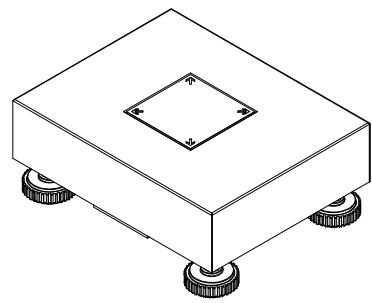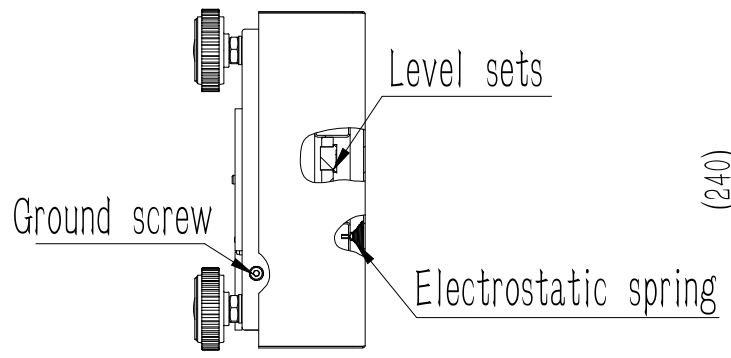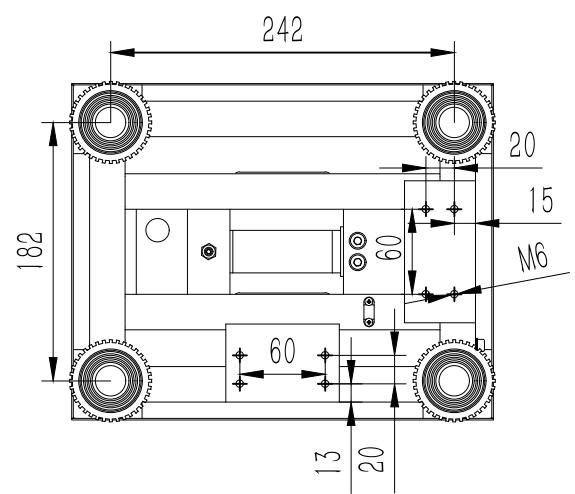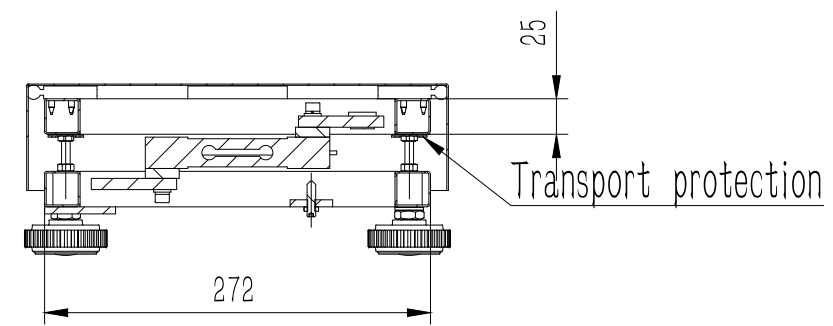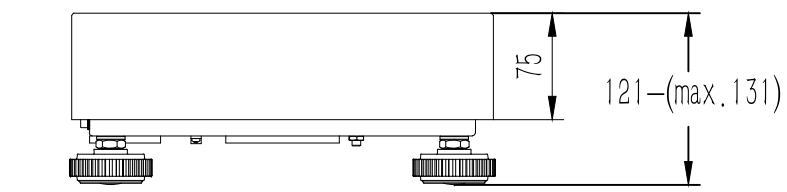
| REVISIONS |      |                             |           |          |
|-----------|------|-----------------------------|-----------|----------|
| ZONE      | REV. | DESCRIPTION                 | DATE      | APPROVED |
| 1         | B    | REDRAWN FROM 021450322_REVA | 9/19/2023 | ZHUANG   |
|           |      |                             |           |          |

|                                                                                                            |  |                                                                       |  |                                  |                                          |          |  |
|------------------------------------------------------------------------------------------------------------|--|-----------------------------------------------------------------------|--|----------------------------------|------------------------------------------|----------|--|
| TOLERANCE UNLESS OTHERWISE SPECIFIED<br>DEC ANGLES<br>±0.1 FOR .X<br>±0.05 FOR .XX                         |  | ±0.5                                                                  |  |                                  | <b><u>KERN</u>® KERN &amp; Sohn GmbH</b> |          |  |
|  3RD ANGLE PROJECTION |  | TITLE: TKBP 15V20M-A                                                  |  |                                  |                                          |          |  |
| DRAWN ZHUANG                                                                                               |  | MODEL NO. B                                                           |  | PART NO. 021450322               |                                          |          |  |
| CHECKED LIN                                                                                                |  | MATERIAL Weighing plate, stainless steel,<br>Platform, steel, coated, |  | FINISH Powder coat "pantone 445" |                                          |          |  |
| UNIT mm                                                                                                    |  | SCALE                                                                 |  | DO NOT SCALE DRAWING             |                                          | SHEET OF |  |

| ZONE | REV. | DESCRIPTION                 | DATE      | APPROVED |
|------|------|-----------------------------|-----------|----------|
| 1    | B    | REDRAWN FROM 021450323_REVA | 9/19/2023 | ZHUANG   |
|      |      |                             |           |          |


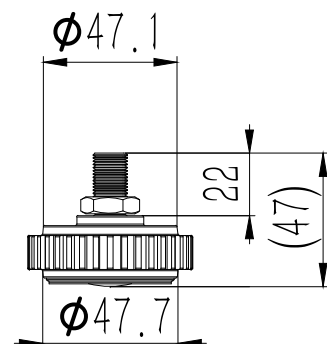
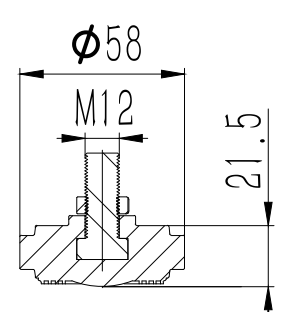
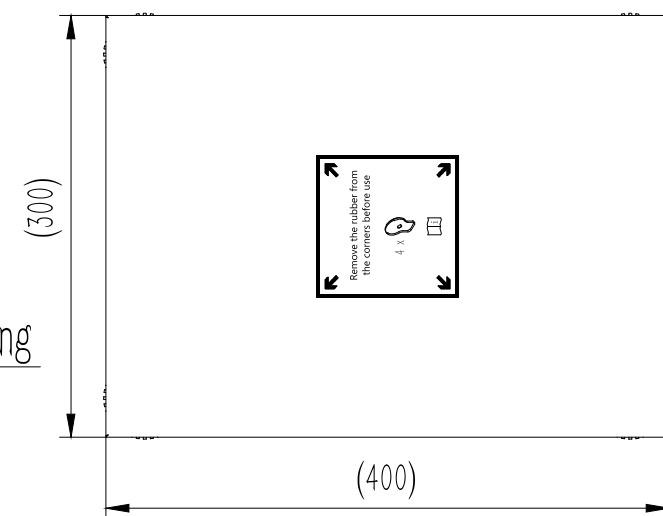
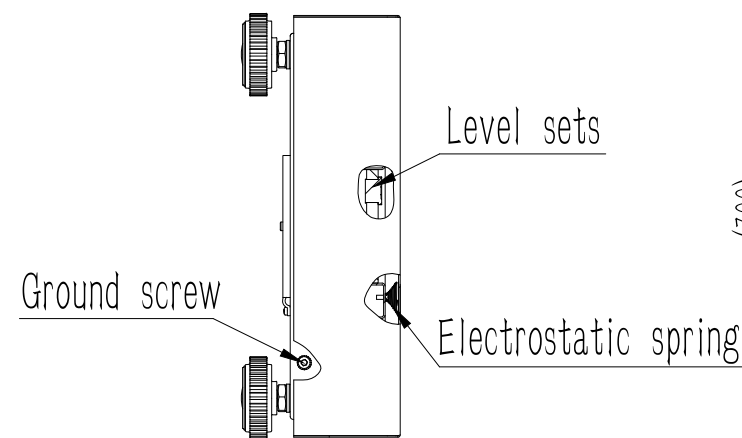
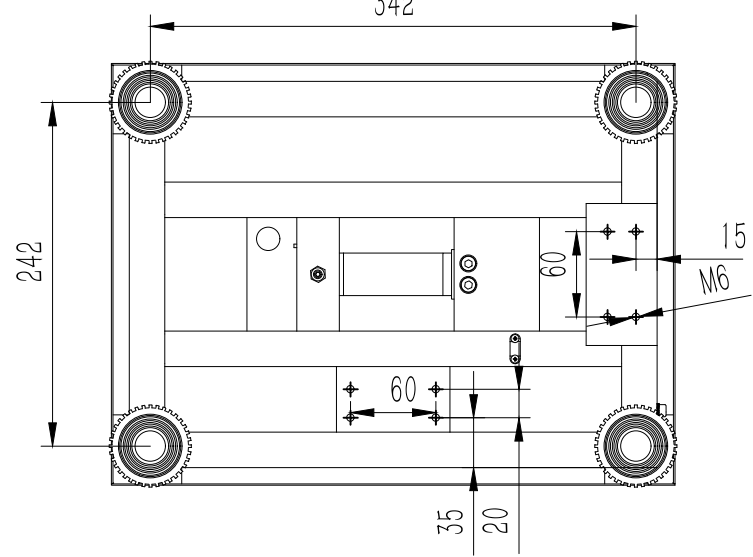
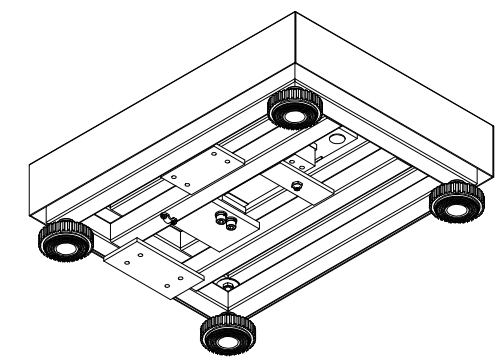
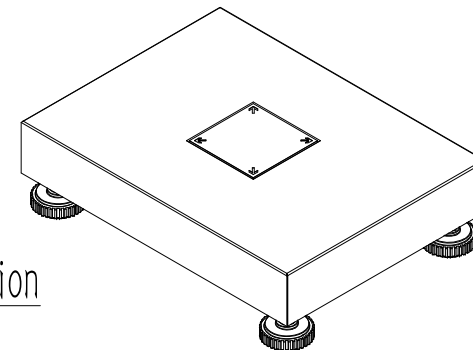
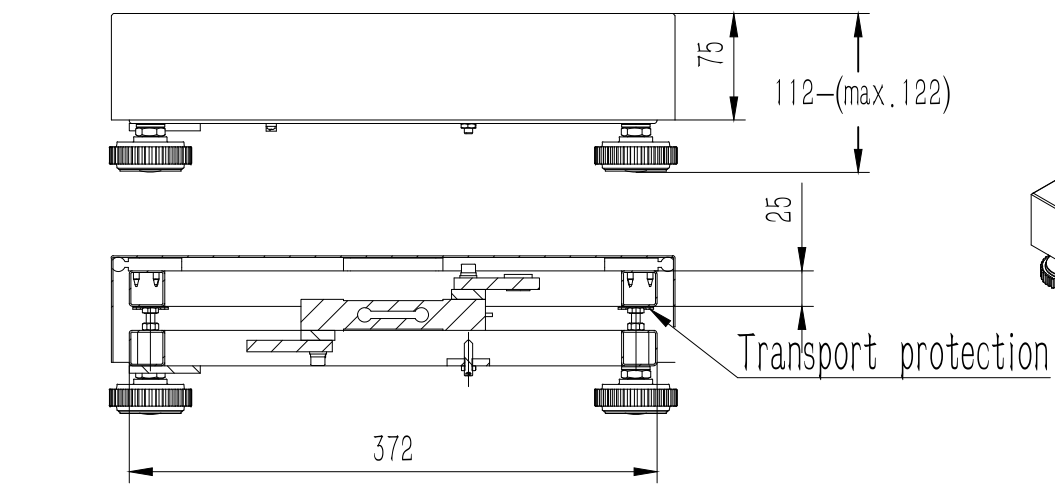
|                                                                                                            |        |                       |                                                             |                                          |                           |
|------------------------------------------------------------------------------------------------------------|--------|-----------------------|-------------------------------------------------------------|------------------------------------------|---------------------------|
| TOLERANCE UNLESS OTHERWISE SPECIFIED<br>DEC ANGLES<br>±0.1 FOR .X<br>±0.05 FOR .XX                         |        | ±0.5                  |                                                             | <b><u>KERN</u>® KERN &amp; Sohn GmbH</b> |                           |
|  3RD ANGLE PROJECTION |        | TITLE: TKBP 15V20LM-A |                                                             |                                          |                           |
| DRAWN                                                                                                      | ZHUANG |                       |                                                             |                                          |                           |
| CHECKED                                                                                                    | LJAN   | MODEL NO.             | B                                                           | PART NO.                                 | 021450323                 |
| APPROVED                                                                                                   | LIN    | MATERIAL              | Weighing plate; stainless steel;<br>Platform; steel coated; | FINISH                                   | Powder coat "pantone 445" |
| UNIT                                                                                                       | mm     | SCALE                 | DO NOT SCALE DRAWING                                        |                                          | SHEET OF                  |
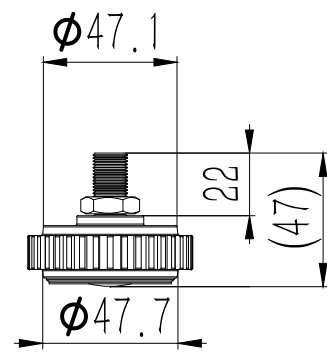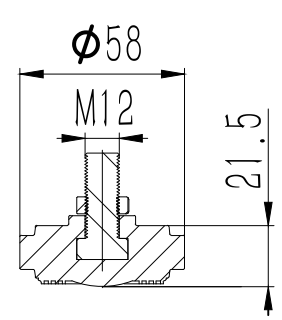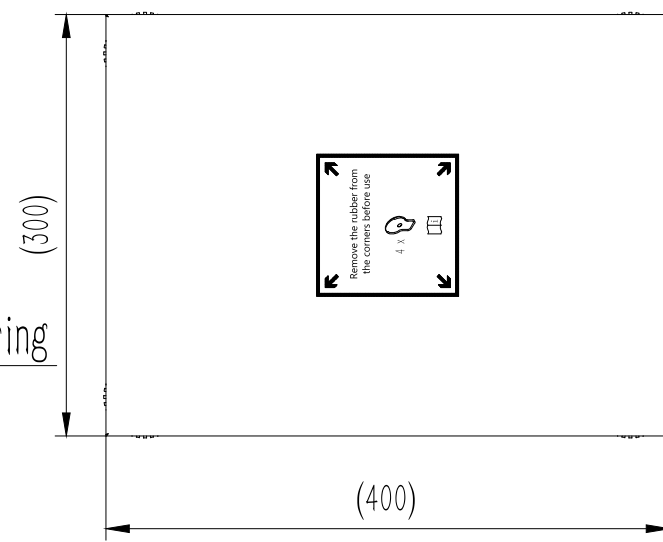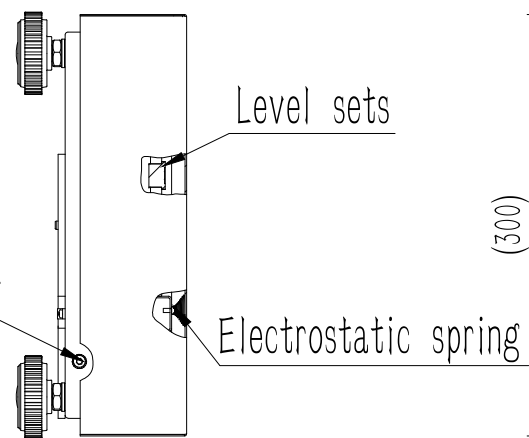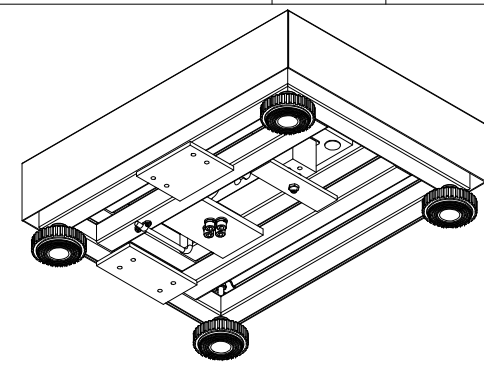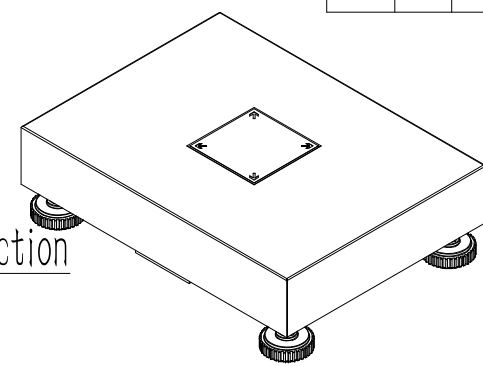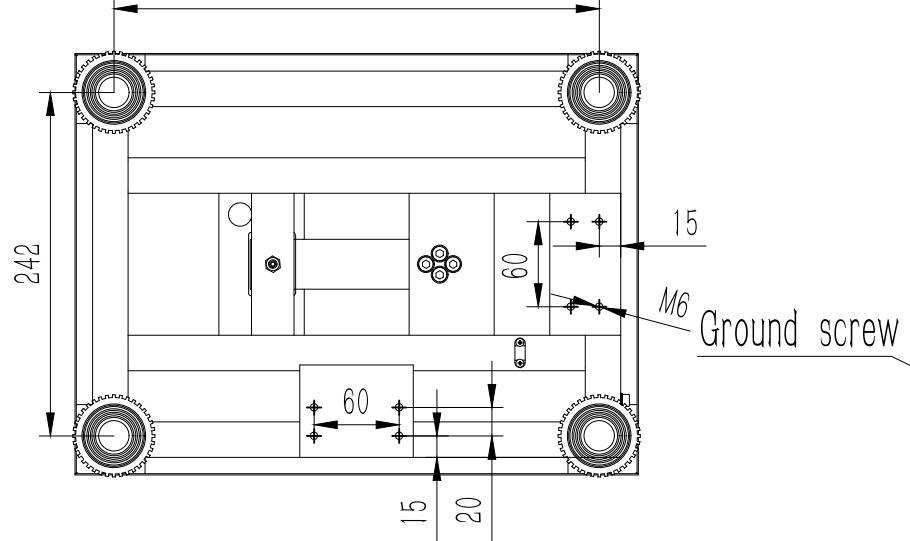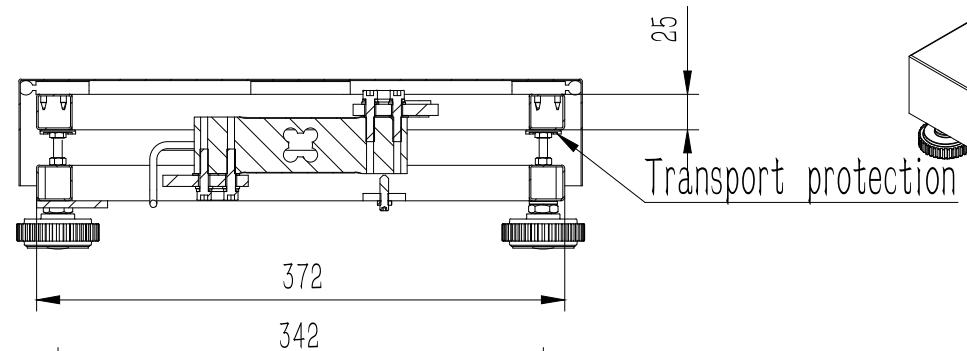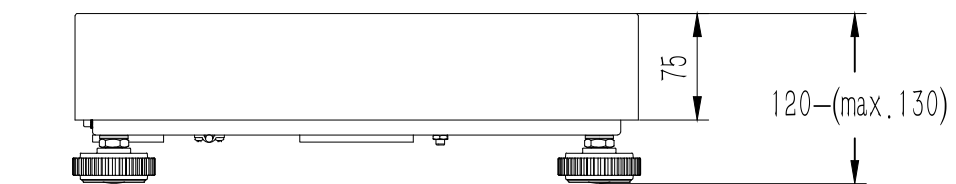
| REVISIONS |      |                             |           |          |
|-----------|------|-----------------------------|-----------|----------|
| ZONE      | REV. | DESCRIPTION                 | DATE      | APPROVED |
| 1         | B    | REDRAWN FROM 021450324_REVA | 9/20/2023 | ZHUANG   |
|           |      |                             |           |          |




|                                                                                                            |        |                                          |                                                             |          |                                  |
|------------------------------------------------------------------------------------------------------------|--------|------------------------------------------|-------------------------------------------------------------|----------|----------------------------------|
| TOLERANCE UNLESS OTHERWISE SPECIFIED<br>DEC ANGLES<br>±0.1 FOR .X<br>±0.05 FOR .XX                         |        | <b><i>KERN® KERN &amp; Sohn GmbH</i></b> |                                                             |          |                                  |
| ±0.5                                                                                                       |        | TITLE: TKBP 30V20M-A                     |                                                             |          |                                  |
|  3RD ANGLE PROJECTION |        |                                          |                                                             |          |                                  |
| DRAWN                                                                                                      | ZHUANG | MODEL NO.                                | B                                                           | PART NO. | 021450324                        |
| CHECKED                                                                                                    | LJAN   | MATERIAL                                 | Weighing plate, stainless steel,<br>Platform; steel coated, |          | FINISH Powder coat "pantone 445" |
| APPROVED                                                                                                   | LIN    | UNIT                                     | mm                                                          | SCALE    | DO NOT SCALE DRAWING             |
|                                                                                                            |        |                                          |                                                             | SHEET OF |                                  |
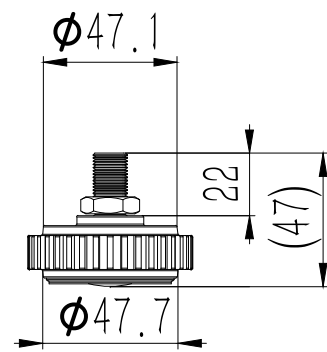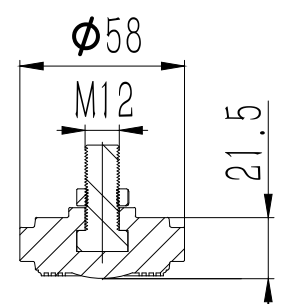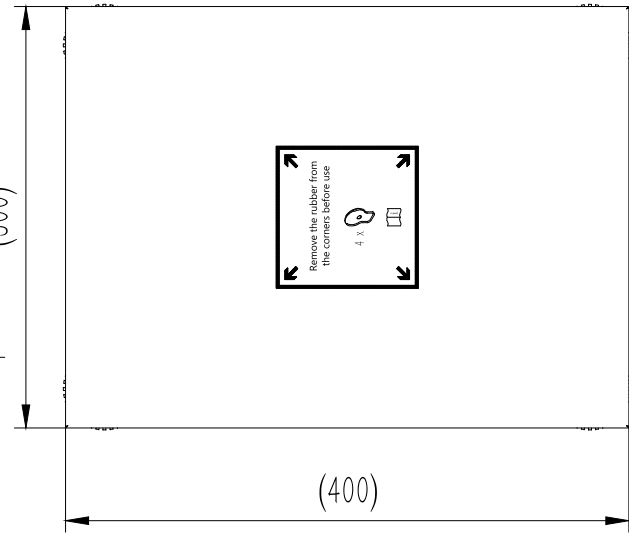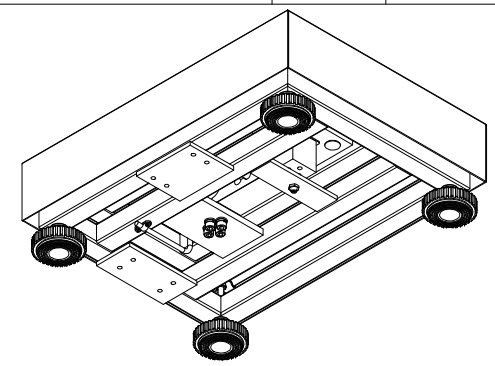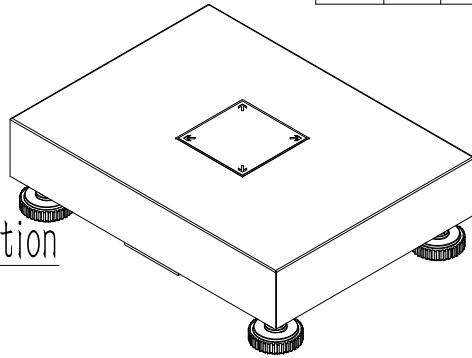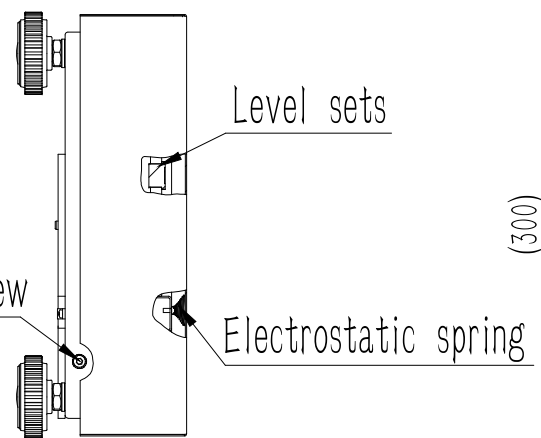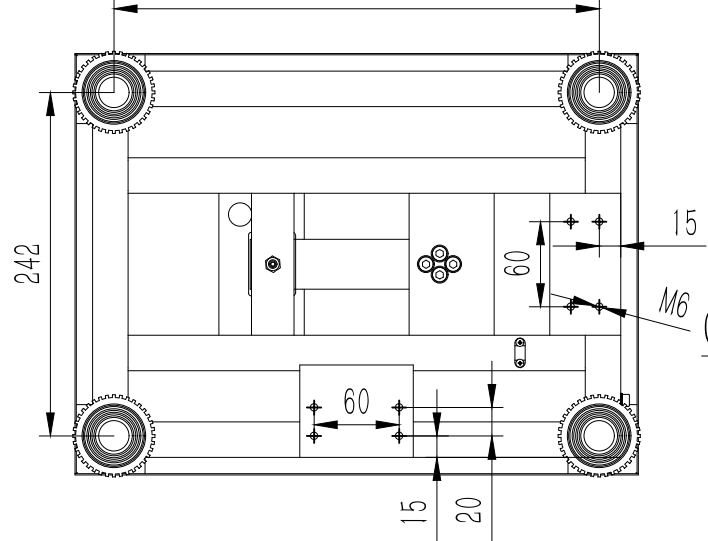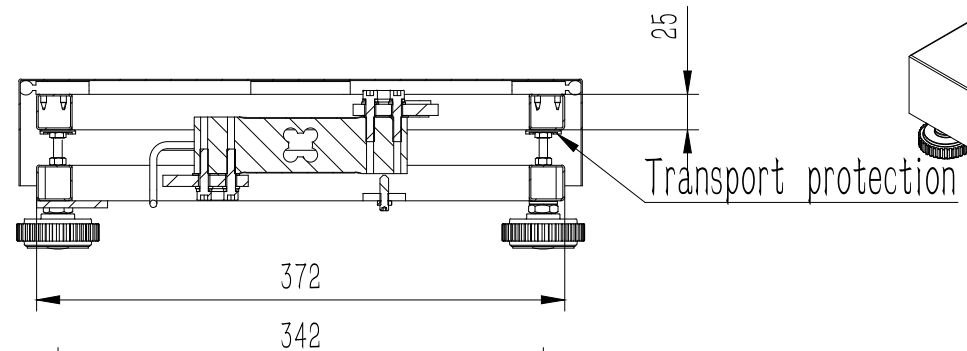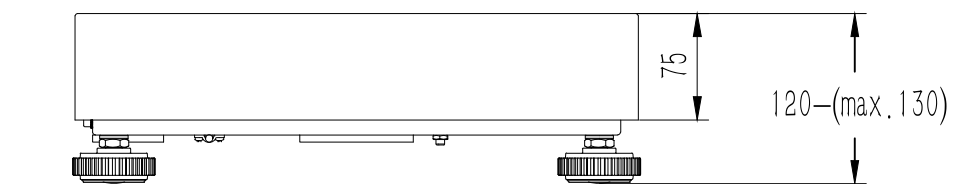
| REVISIONS |      |                             |           |          |
|-----------|------|-----------------------------|-----------|----------|
| ZONE      | REV. | DESCRIPTION                 | DATE      | APPROVED |
| 1         | B    | REDRAWN FROM 021450325_REVA | 9/20/2023 | ZHUANG   |
|           |      |                             |           |          |








|                                                                                       |        |                      |                                                              |                                          |                           |          |
|---------------------------------------------------------------------------------------|--------|----------------------|--------------------------------------------------------------|------------------------------------------|---------------------------|----------|
| TOLERANCE UNLESS OTHERWISE SPECIFIED<br>DEC ANGLES<br>±0.1 FOR .X<br>±0.05 FOR .XX    |        | ±0.5                 |                                                              | <b><u>KERN</u>® KERN &amp; Sohn GmbH</b> |                           |          |
|  |        | 3RD ANGLE PROJECTION |                                                              | TITLE: TKBP 60V20M-A                     |                           |          |
| DRAWN                                                                                 | ZHUANG |                      |                                                              |                                          |                           |          |
| CHECKED                                                                               | LIAN   | MODEL NO.            | B                                                            | PART NO.                                 | 021450325                 |          |
| APPROVED                                                                              | LIN    | MATERIAL             | Weighing plate, stainless steel,<br>Platform, steel, coated, | FINISH                                   | Powder coat "pantone 445" |          |
| UNIT                                                                                  | mm     | SCALE                |                                                              | DO NOT SCALE DRAWING                     |                           | SHEET OF |

| ZONE | REV. | DESCRIPTION                 | DATE      | APPROVED |
|------|------|-----------------------------|-----------|----------|
| 1    | B    | REDRAWN FROM 021450326_REVA | 9/20/2023 | ZHUANG   |
|      |      |                             |           |          |








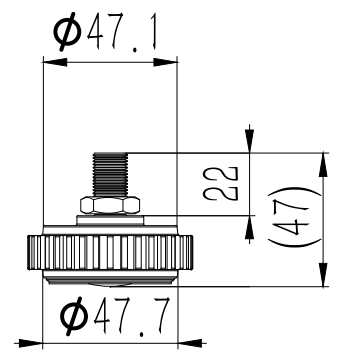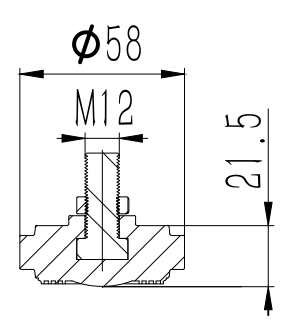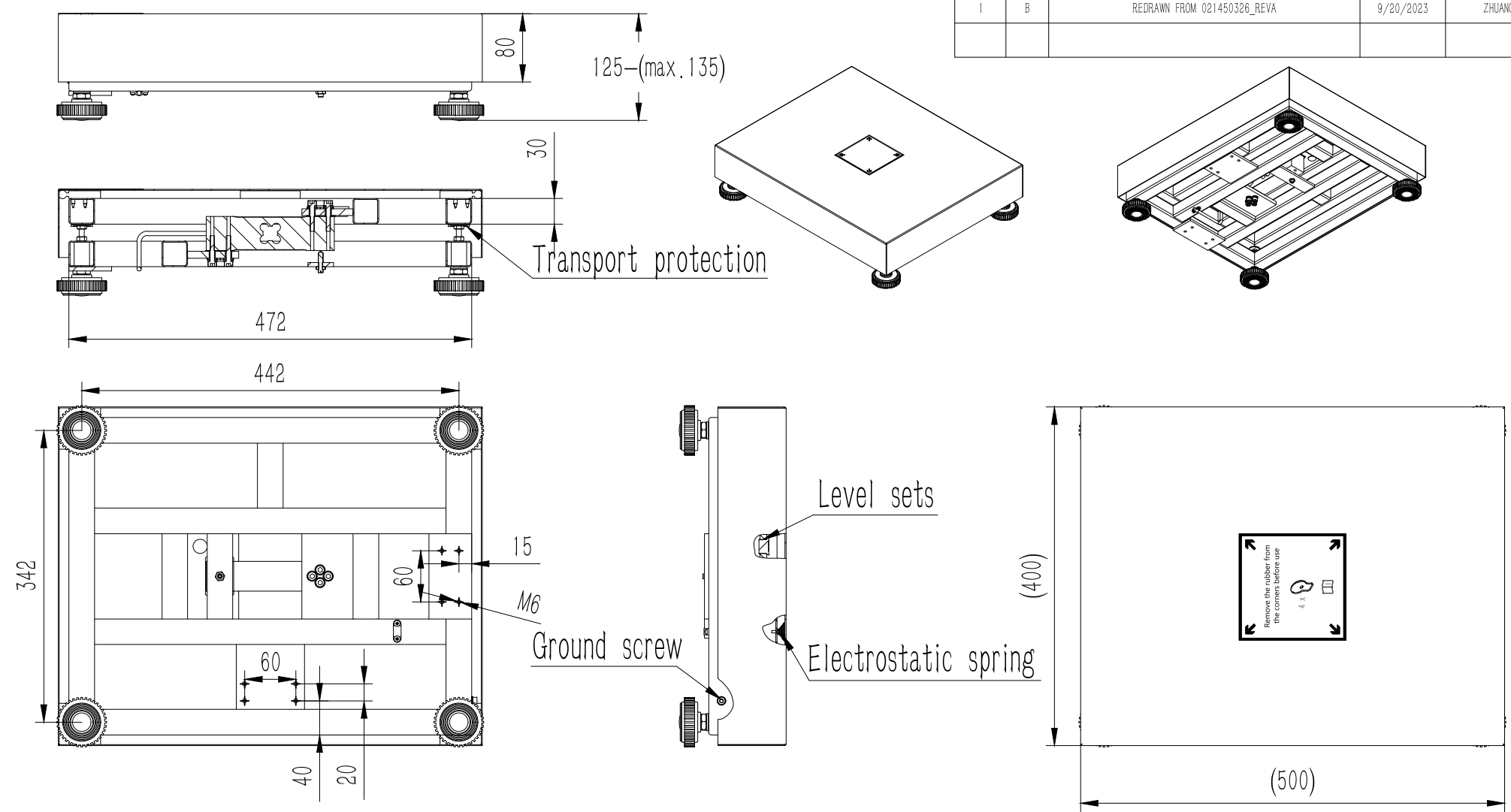
|                                                                                                            |        |                       |                                                             |                                          |                           |
|------------------------------------------------------------------------------------------------------------|--------|-----------------------|-------------------------------------------------------------|------------------------------------------|---------------------------|
| TOLERANCE UNLESS OTHERWISE SPECIFIED<br>DEC ANGLES<br>±0.1 FOR .X<br>±0.05 FOR .XX                         |        | ±0.5                  |                                                             | <b><u>KERN</u>® KERN &amp; Sohn GmbH</b> |                           |
|  3RD ANGLE PROJECTION |        | TITLE: TKBP 60V20LM-A |                                                             |                                          |                           |
| DRAWN                                                                                                      | ZHUANG | MODEL NO.             | B                                                           | PART NO.                                 | 021450326                 |
| CHECKED                                                                                                    | LIAN   | MATERIAL              | Weighing plate, stainless steel,<br>Platform, steel coated, | FINISH                                   | Powder coat "pantone 445" |
| APPROVED                                                                                                   | LIN    | SCALE                 |                                                             | DO NOT SCALE DRAWING                     | SHEET OF                  |
| UNIT                                                                                                       | mm     |                       |                                                             |                                          |                           |

| ZONE | REV. | DESCRIPTION                 | DATE      | APPROVED |
|------|------|-----------------------------|-----------|----------|
| 1    | B    | REDRAWN FROM 021450327_REVA | 9/20/2023 | ZHUANG   |
|      |      |                             |           |          |
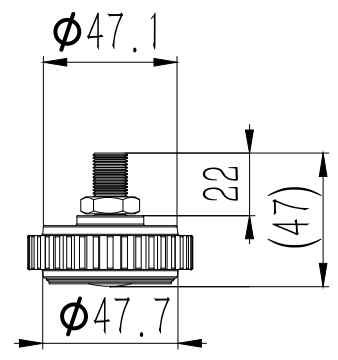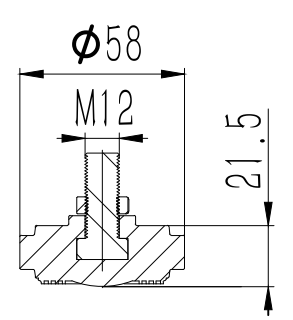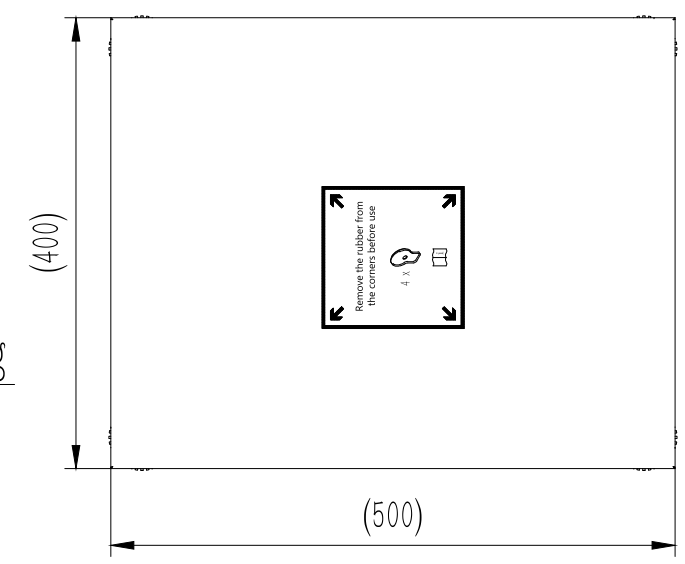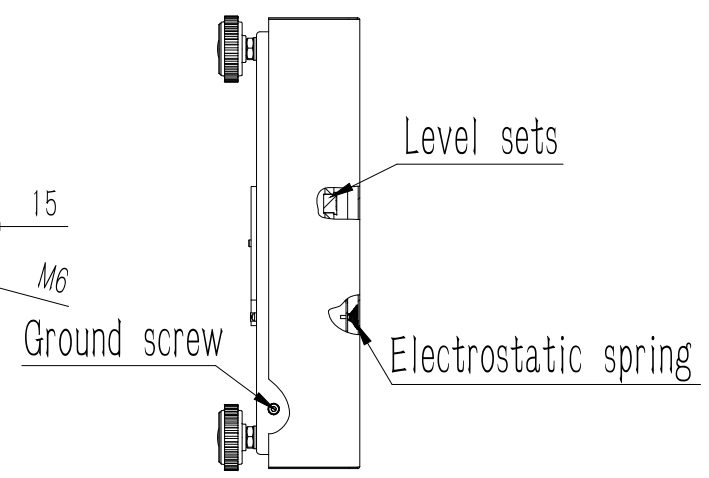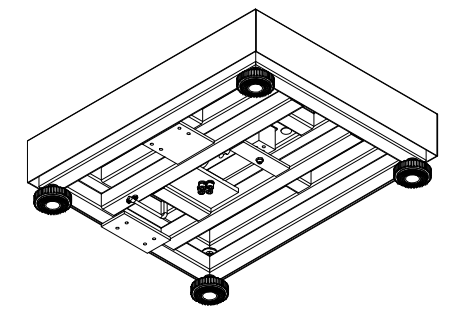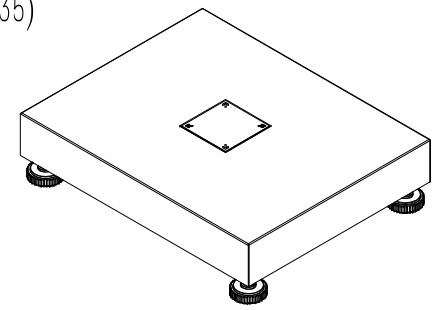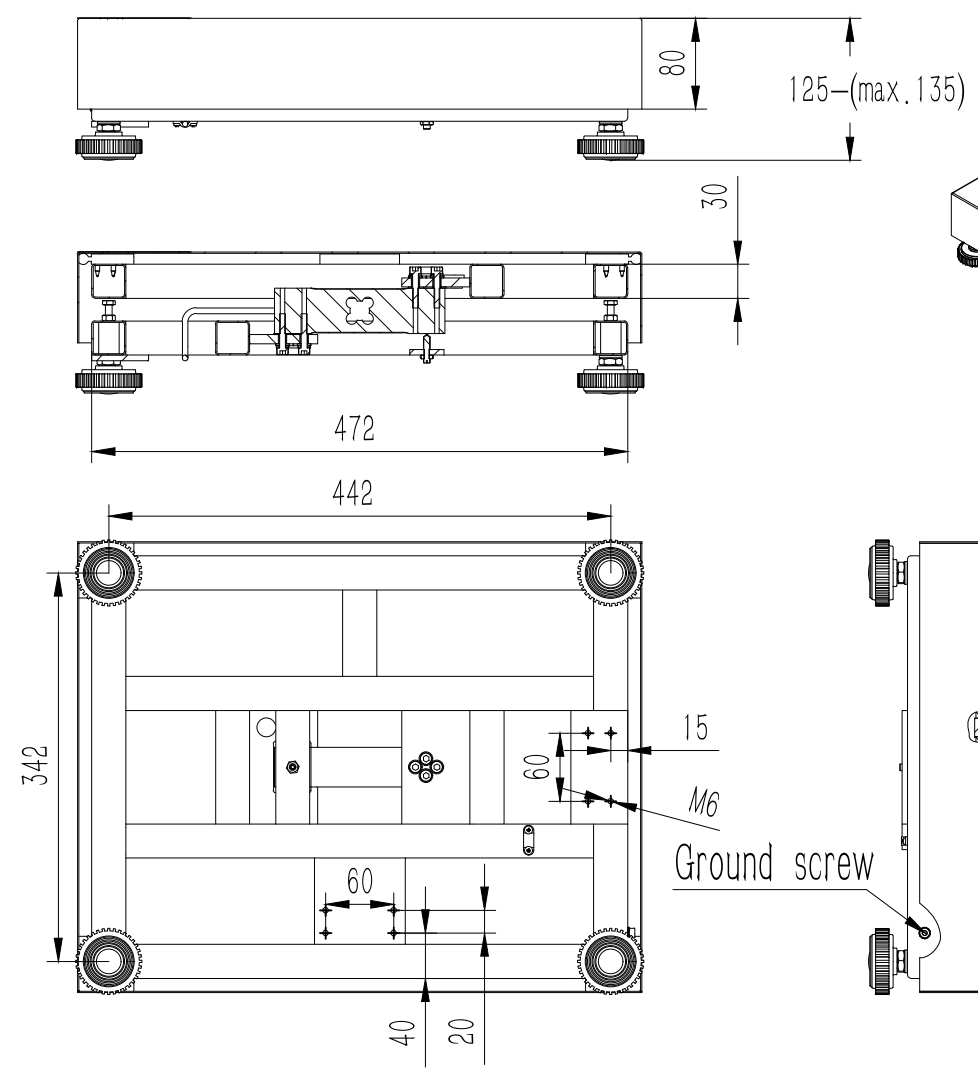
|                                                                                       |        |                      |                                                             |                                          |                                  |
|---------------------------------------------------------------------------------------|--------|----------------------|-------------------------------------------------------------|------------------------------------------|----------------------------------|
| TOLERANCE UNLESS OTHERWISE SPECIFIED<br>DEC ANGLES<br>±0.1 FOR .X<br>±0.05 FOR .XX    |        | ±0.5                 |                                                             | <b><u>KERN</u>® KERN &amp; Sohn GmbH</b> |                                  |
|  |        | 3RD ANGLE PROJECTION |                                                             | TITLE: TKBP 150V20M-A                    |                                  |
| DRAWN                                                                                 | ZHUANG |                      |                                                             |                                          |                                  |
| CHECKED                                                                               | LIAN   | MODEL NO.            | B                                                           | PART NO.                                 | 021450327                        |
| APPROVED                                                                              | LIN    | MATERIAL             | Weighing plate; stainless steel,<br>Platform; steel coated, |                                          | FINISH Powder coat "pantone 445" |
| UNIT                                                                                  | mm     | SCALE                |                                                             | DO NOT SCALE DRAWING                     | SHEET OF                         |









| REVISIONS |      |                             |           |          |
|-----------|------|-----------------------------|-----------|----------|
| ZONE      | REV. | DESCRIPTION                 | DATE      | APPROVED |
| 1         | B    | REDRAWN FROM 021450328_REVA | 9/20/2023 | ZHUANG   |
|           |      |                             |           |          |
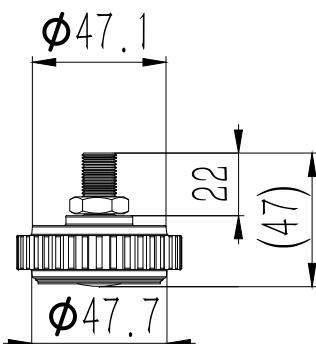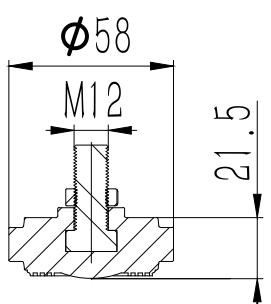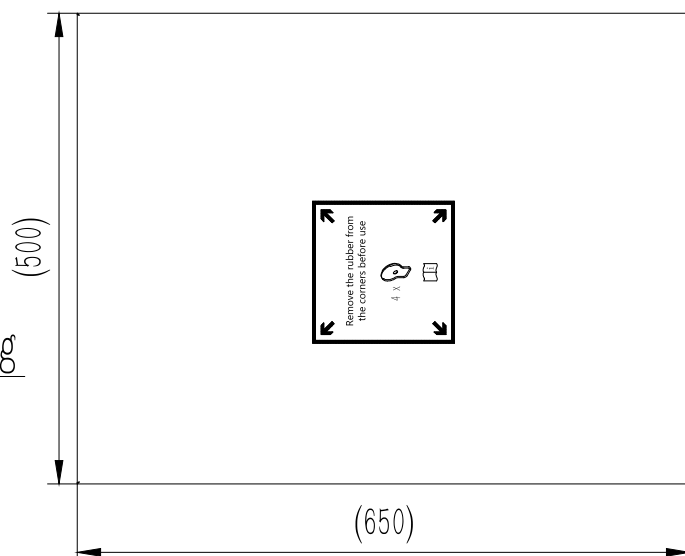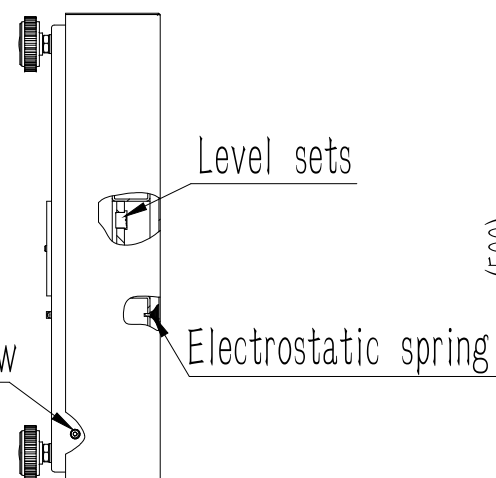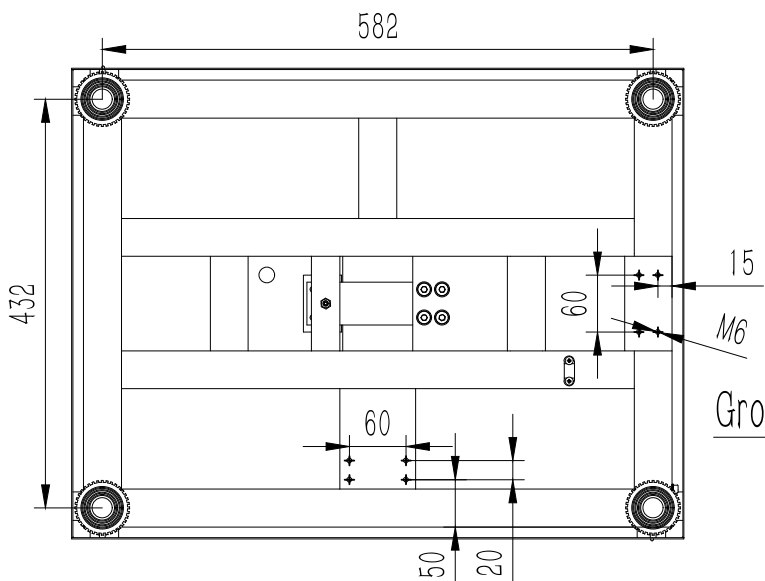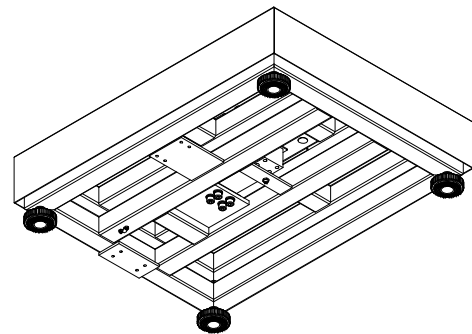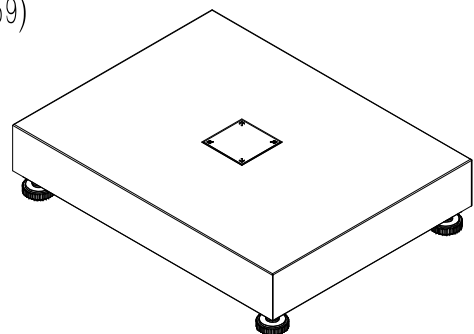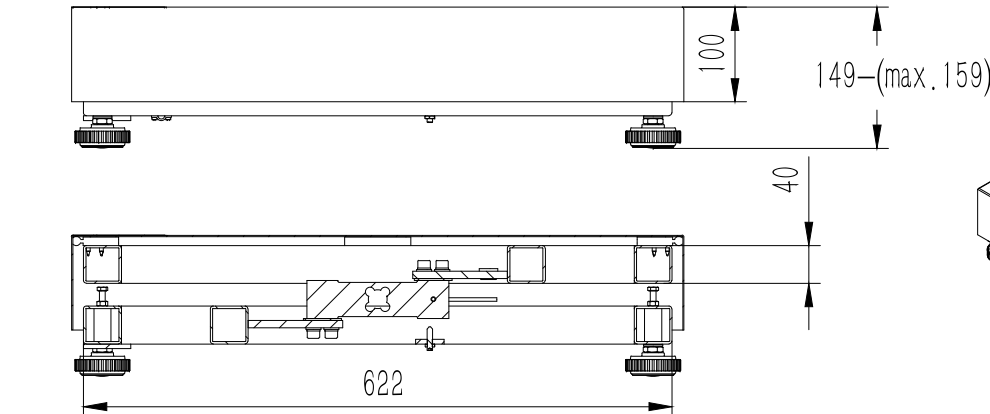
|                                                                                       |    |                      |  |                                          |                                                             |          |           |                           |
|---------------------------------------------------------------------------------------|----|----------------------|--|------------------------------------------|-------------------------------------------------------------|----------|-----------|---------------------------|
| TOLERANCE UNLESS OTHERWISE SPECIFIED<br>DEC ANGLES<br>±0.1 FOR .X<br>±0.05 FOR .XX    |    | ±0.5                 |  | <b><u>KERN®</u> KERN &amp; Sohn GmbH</b> |                                                             |          |           |                           |
|  |    | 3RD ANGLE PROJECTION |  | TITLE: TKBP 150V20LM-A                   |                                                             |          |           |                           |
| DRAWN                                                                                 |    | ZHUANG               |  |                                          |                                                             |          |           |                           |
| CHECKED                                                                               |    | LIAN                 |  | MODEL NO.                                | B                                                           | PART NO. | 021450328 |                           |
| APPROVED                                                                              |    | LIN                  |  | MATERIAL                                 | Weighing plate, stainless steel,<br>Platform, steel coated, |          | FINISH    | Powder coat "pantone 445" |
| UNIT                                                                                  | mm | SCALE                |  | DO NOT SCALE DRAWING                     |                                                             |          | SHEET OF  |                           |








| REVISIONS |      |                             |           |          |
|-----------|------|-----------------------------|-----------|----------|
| ZONE      | REV. | DESCRIPTION                 | DATE      | APPROVED |
| 1         | B    | REDRAWN FROM 021450329_REVA | 9/20/2023 | ZHUANG   |
|           |      |                             |           |          |
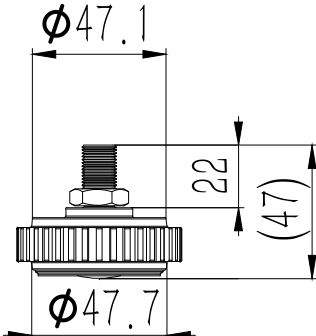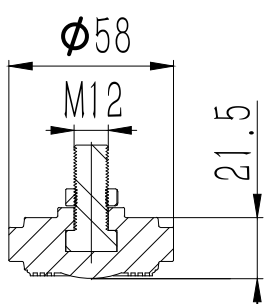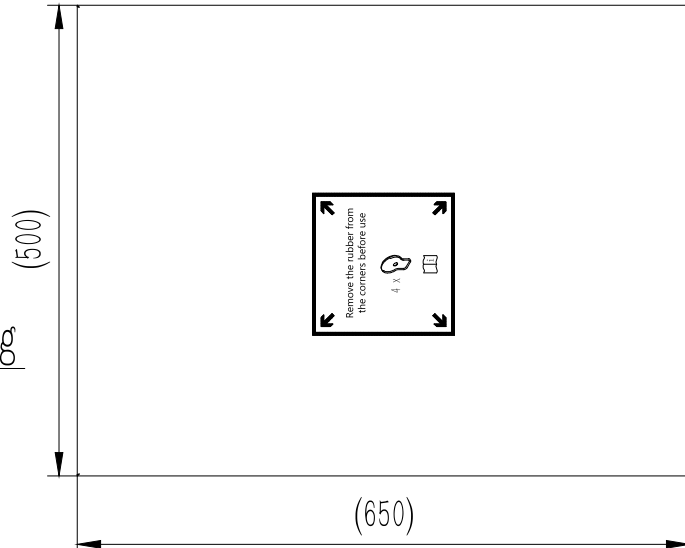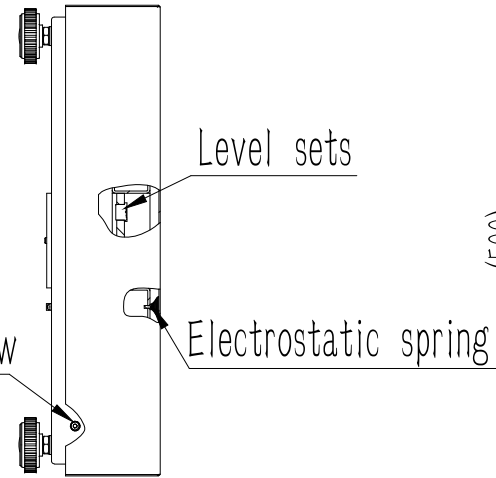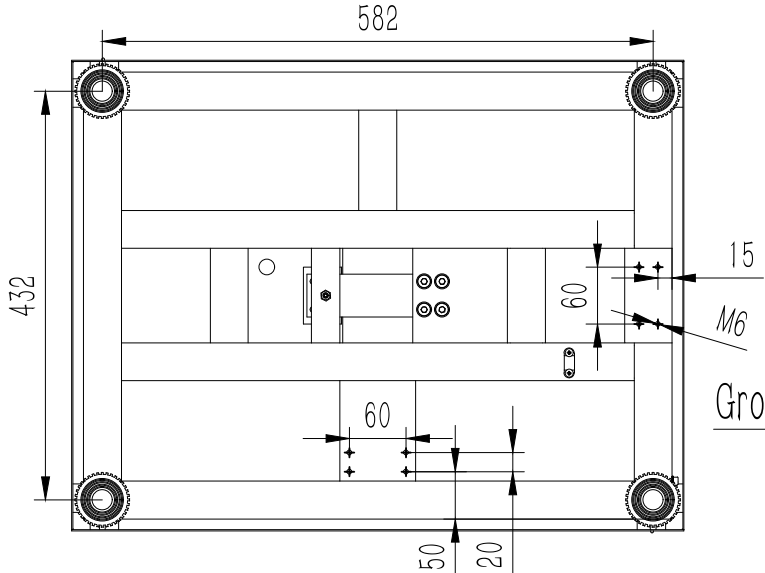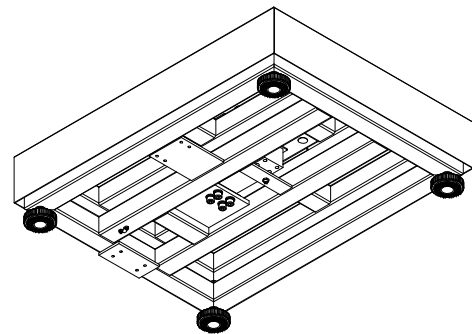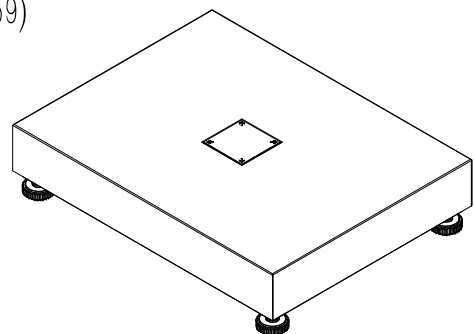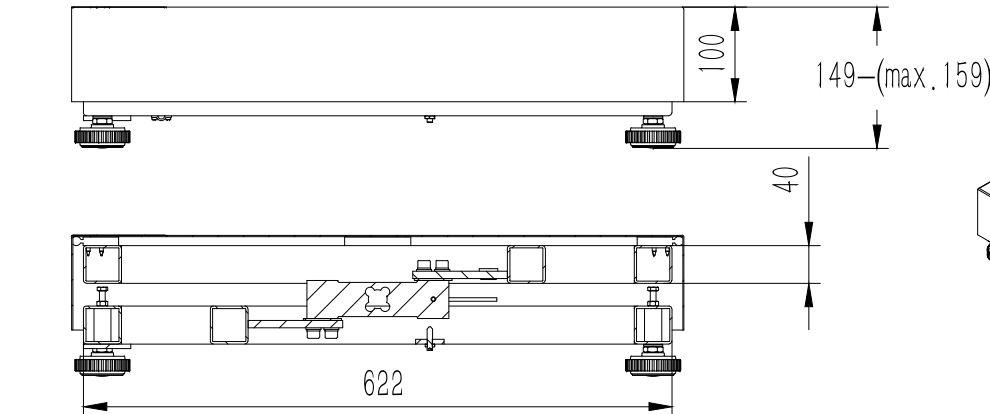
|                                                                                       |        |                      |           |                                                              |                                  |
|---------------------------------------------------------------------------------------|--------|----------------------|-----------|--------------------------------------------------------------|----------------------------------|
| TOLERANCE UNLESS OTHERWISE SPECIFIED<br>DEC ANGLES<br>±0.1 FOR .X<br>±0.05 FOR .XX    |        | ±0.5                 |           | <b><u>KERN®</u> KERN &amp; Sohn GmbH</b>                     |                                  |
|  |        | 3RD ANGLE PROJECTION |           | TITLE: TKBP 300V20M-A                                        |                                  |
| DRAWN                                                                                 | ZHUANG |                      | MODEL NO. | B                                                            | PART NO. 021450329               |
| CHECKED                                                                               | LIAN   |                      | MATERIAL  | Weighing plate, stainless steel,<br>Platform, steel, coated, | FINISH Powder coat "pantone 445" |
| APPROVED                                                                              | LIN    |                      | UNIT      | mm                                                           | SCALE DO NOT SCALE DRAWING       |
|                                                                                       |        |                      |           | SHEET OF                                                     |                                  |

| REVISIONS |      |                             |           |          |
|-----------|------|-----------------------------|-----------|----------|
| ZONE      | REV. | DESCRIPTION                 | DATE      | APPROVED |
| 1         | B    | REDRAWN FROM 021450330_REVA | 9/20/2023 | ZHUANG   |
|           |      |                             |           |          |
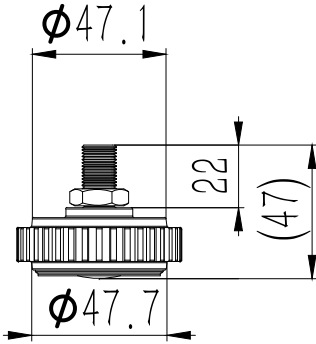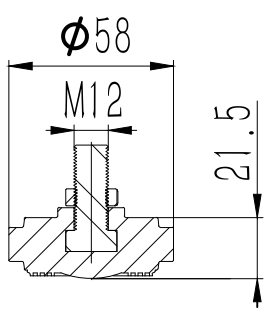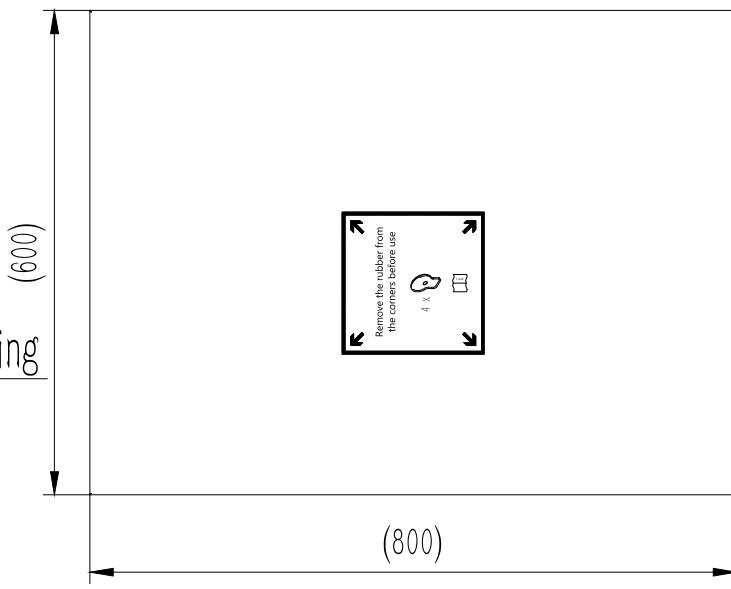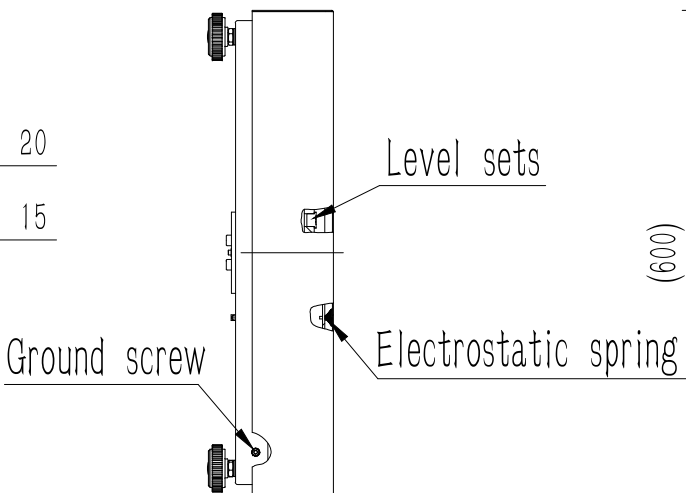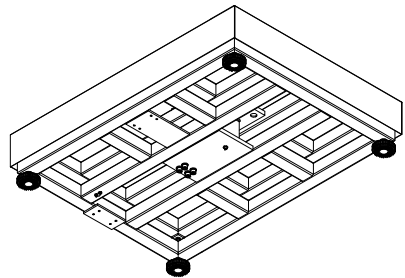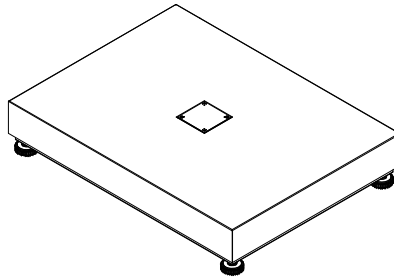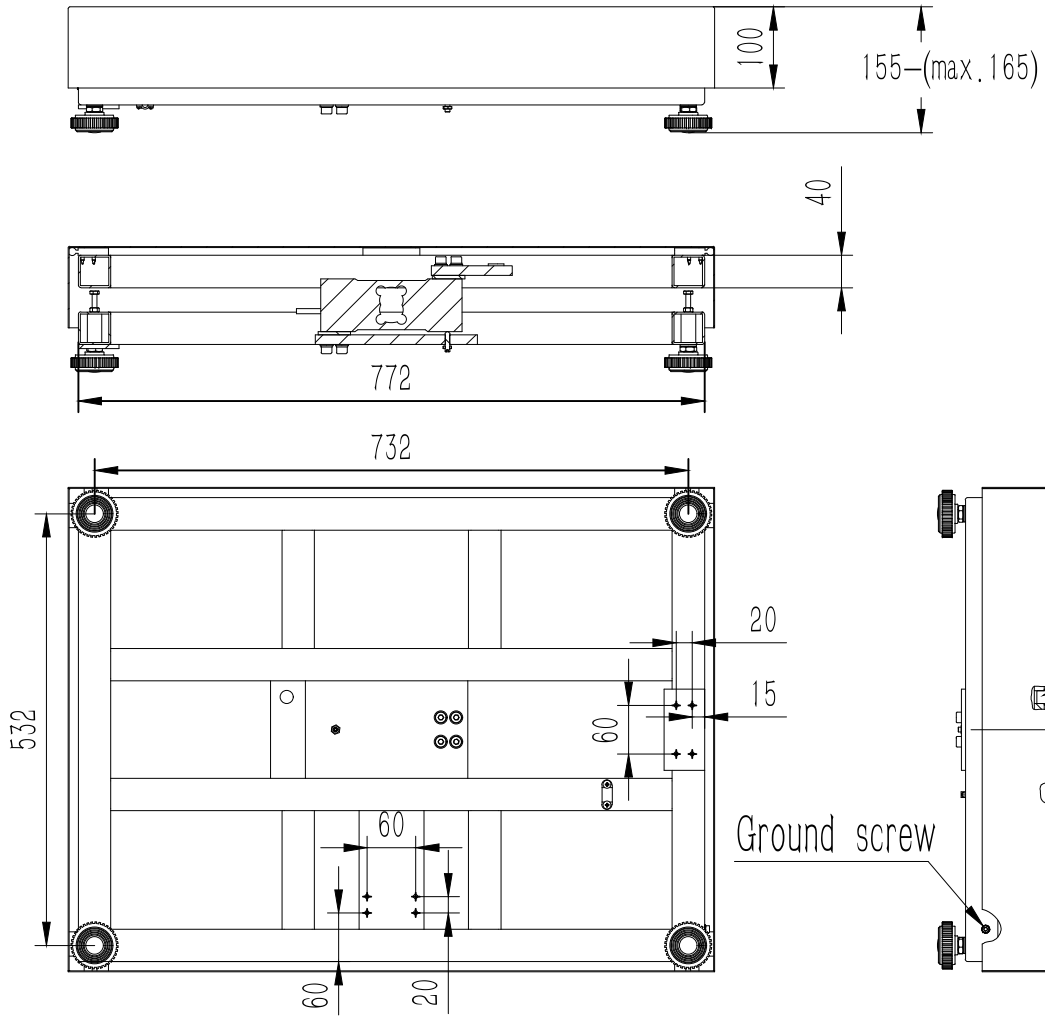
|                                                                                       |        |                      |           |                                                             |                      |                                  |
|---------------------------------------------------------------------------------------|--------|----------------------|-----------|-------------------------------------------------------------|----------------------|----------------------------------|
| TOLERANCE UNLESS OTHERWISE SPECIFIED<br>DEC ANGLES<br>±0.1 FOR .X<br>±0.05 FOR .XX    |        | ±0.5                 |           | <b><u>KERN</u>® KERN &amp; Sohn GmbH</b>                    |                      |                                  |
|  |        | 3RD ANGLE PROJECTION |           | TITLE: TKBP 600V20M-A                                       |                      |                                  |
| DRAWN                                                                                 | ZHUANG |                      |           |                                                             |                      |                                  |
| CHECKED                                                                               | LIAN   |                      | MODEL NO. | B                                                           | PART NO.             | 021450330                        |
| APPROVED                                                                              | LIN    |                      | MATERIAL  | Weighing plate, stainless steel,<br>Platform, steel coated, |                      | FINISH Powder coat "pantone 445" |
| UNIT                                                                                  | mm     |                      | SCALE     |                                                             | DO NOT SCALE DRAWING | SHEET OF                         |