

USER MANUAL

INFRARED STERILIZER

PCE-IRS 150 / PCE-IRS 240



ENGLISH



User manuals in various languages (français, italiano, español, português, nederlands, türk, polski) can be found via our product search on:
www.pce-instruments.com

BRIEF INTRODUCTION

This brand infrared sterilizer adopts international advanced design concepts and manufacturing technology, with smooth and beautiful appearance, compact and firm structure, high versatility, simple operation and other advantages. It is a rare instrument of scientific research and experiment. This instrument is mainly used for high-temperature sterilization and disinfection of small items such as inoculation loops and inoculation needles. It can replace alcohol lamps and is convenient and quick.

PRODUCT INTRODUCTION

- 1 - Desinfection hole
- 2 - Switch/gear
- 3 - Plum socket



TECHNICAL DATA

Model	PCE-IRS 150	PCE-IRS 240
Voltage	AC220V 50Hz	
Power	150	240
Max. temperature inside the heating hole	825°C ±50°C	
Heating temperature	Two temperatures available: 400°C / 850°C	
Max. diameter of sterilizer goods	ø 14	ø 35
Length of heating zone	150	100
Warm-up time	10 ... 20 min	
Fuse	250V 3A 5x20	

Model	PCE-IRS 150	PCE-IRS 240
Ambient temperature	5 ... 40°C	
Relative humidity	≤90%	
Product size	175x130x185mm	147x130x200
Net weight	1.3	1.4

OPERATION GUIDE

Application

- » It is mainly used for microbiological experiments in biological safety cabinets, purification workbenches, mobile vehicles and other environments.

Applied area

- » Used for high temperature sterilization of small items such as inoculation loops and inoculation needles

Using Steps

- » Place the instrument on a stable workbench, file the switch in the OFF position
- » Plug in
- » Turn the switch to the required position



Instrument use environment

- » Indoor use
- » The working environment of the instrument is suitable for the temperature range of +5°C ... +40°C
- » The applicable temperature range of the instrument is ≤80%
- » There is no vibration and airflow that affect performance around
- » There is no conductive dust, explosive gas or corrosive gas in the surrounding air
- » Keep a safety distance of 10cm before and after the instrument is working



Safety Precautions

- » It is strictly prohibited to plug in the power connector and toggle the power button when the hands are wet with liquid!
- » It is strictly prohibited to unplug the power plug when the instrument is live!
- » It is strictly forbidden to maintain and clean the instrument when it is powered on!
- » It is strictly forbidden to install the instrument on an uneven, swaying and vibrating work surface!

DISPOSAL

For the disposal of batteries in the EU, the (EU) 2023/1542 directive of the European Parliament applies. Due to the contained pollutants, batteries must not be disposed of as household waste. They must be given to collection points designed for that purpose. In order to comply with the EU directive 2012/19/EU we take our devices back. We either re-use them or give them to a recycling company which disposes of the devices in line with law. For countries outside the EU, batteries and devices should be disposed of in accordance with your local waste regulations. If you have any questions, please contact PCE Instruments.

PCE INSTRUMENTS CONTACT INFORMATION

Germany

PCE Deutschland GmbH
Im Langel 26
D-59872 Meschede
Deutschland
Tel.: +49 (0) 2903 976 99 0
Fax: +49 (0) 2903 976 99 29
info@pce-instruments.com
www.pce-instruments.com/deutsch

France

PCE Instruments France EURL
2, rue Georges Kuhnmmunch
67250 Soultz-sous-Forêts
France
Tel.: +33 (0) 972 35 37 17
Fax: +33 (0) 972 35 37 18
info@pce-france.fr
www.pce-instruments.com/french

Spain

PCE Ibérica S.L.
Calle Mula, 8
02500 Tobarra (Albacete)
España
Tel.: +34 967 543 548
info@pce-iberica.es
www.pce-instruments.com/espanol

United Kingdom

PCE Instruments UK Ltd
Trafford House
Chester Rd, Old Trafford
Manchester M32 0RS
United Kingdom
Tel: +44 (0) 161 464902 0
Fax: +44 (0) 161 464902 9
info@pce-instruments.co.uk
www.pce-instruments.com/english

Italy

PCE Italia s.r.l.
Via Pesciatina 878 / B-Interno 6
55010 Loc. Gragnano
Capannori (Lucca)
Italia
Telefono: +39 0583 975 114
Fax: +39 0583 974 824
info@pce-italia.it
www.pce-instruments.com/italiano

Turkey

PCE Teknik Cihazları Ltd.Şti.
Halkalı Merkez Mah.
Pehlivan Sok. No.6/C
34303 Küçükçekmece - İstanbul
Türkiye
Tel: 0212 471 11 47
Faks: 0212 705 53 93
info@pce-cihazlari.com.tr
www.pce-instruments.com/turkish

The Netherlands

PCE Brookhuis B.V.
Twentepoort West 17
7609 RD Almelo
Nederland
Telefoon: +31 (0)53 737 01 92
info@pcebenelux.nl
www.pce-instruments.com/dutch

United States of America

PCE Americas Inc.
1201 Jupiter Park Drive, Suite 8
Jupiter / Palm Beach
33458 FL
USA
Tel: +1 (561) 320-9162
Fax: +1 (561) 320-9176
info@pce-americas.com
www.pce-instruments.com/us

Denmark

PCE Instruments Denmark ApS
Birk Centerpark 40
7400 Herning
Denmark
Tel.: +45 70 30 53 08
kontakt@pce-instruments.com
www.pce-instruments.com/dansk



Subject to change without notice