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Ziegelei 1

D-72336 Balingen

E-Mail: [info@kern-sohn.com](mailto:info@kern-sohn.com)

Tel: +49-[0]7433- 9933-0

Fax: +49-[0]7433-9933-149

Internet: [www.kern-sohn.com](http://www.kern-sohn.com)

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Density determination set for precision balance KERN EMB 2000-2V

## 1 Scope of delivery

- ⇒ Check packaging and density determination set immediately when unpacking for possible visible damage.
- ⇒ Make sure that all parts are completely present.

### 1 Weighing plate "density set"


### 2 Platform

3 Beaker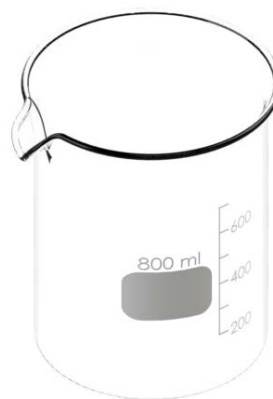

4 Immersion basket for descending solid matter (density  $> 1 \text{ g/cm}^3$ )

5 Immersion basket for floating solid matter (density  $< 1 \text{ g/cm}^3$ )

6 Sinker  
200 g stainless steel weight





7 Thermometer


8 Manual



## 2 Dimension [mm]

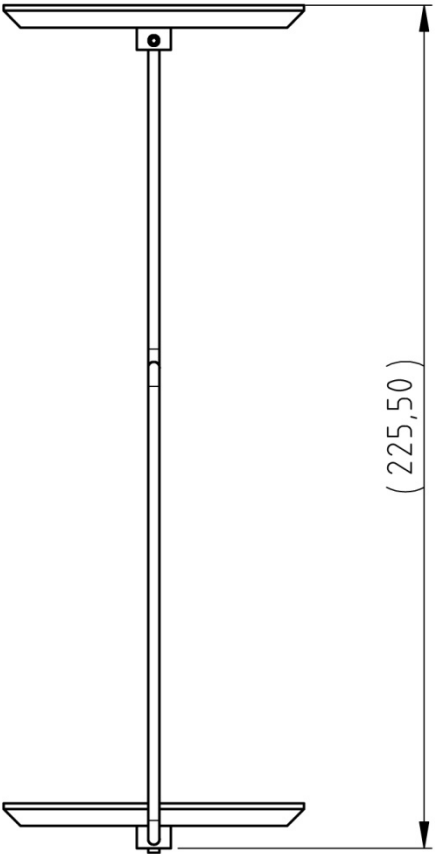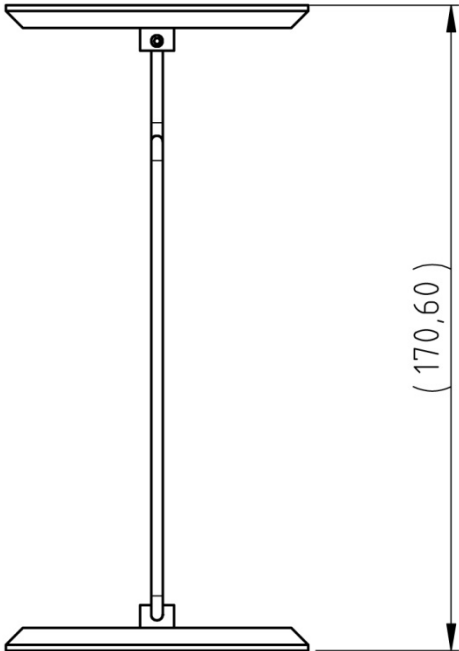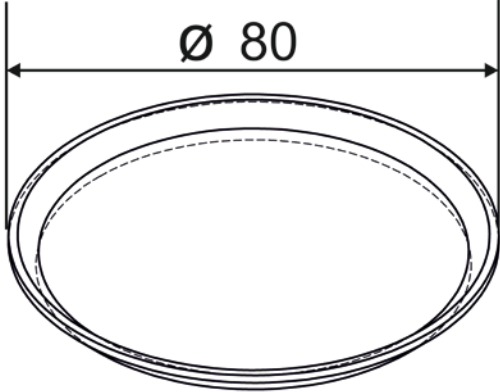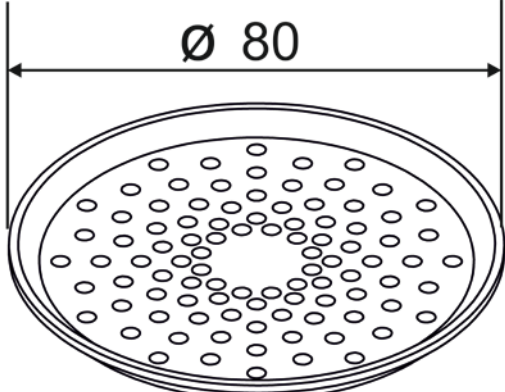Immersion basket for descending solid matter (density > 1 g/cm <sup>3</sup> )	Immersion basket for floating solid matter (density < 1 g/cm <sup>3</sup> )
 <p data-bbox="646 779 694 913">( 225,50 )</p>	 <p data-bbox="1268 683 1316 817">( 170,60 )</p>
Sample dish	
 <p data-bbox="411 1422 534 1467">Ø 80</p>	 <p data-bbox="1013 1422 1136 1467">Ø 80</p>

Beaker

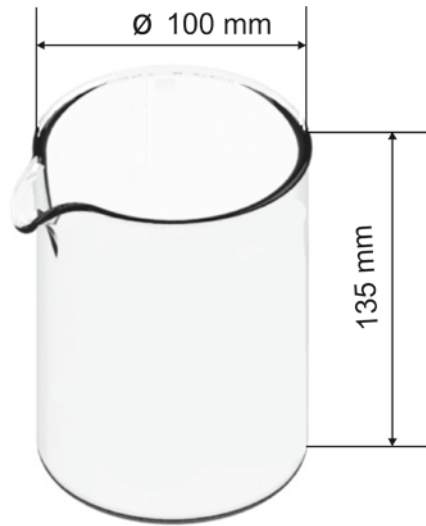Platform
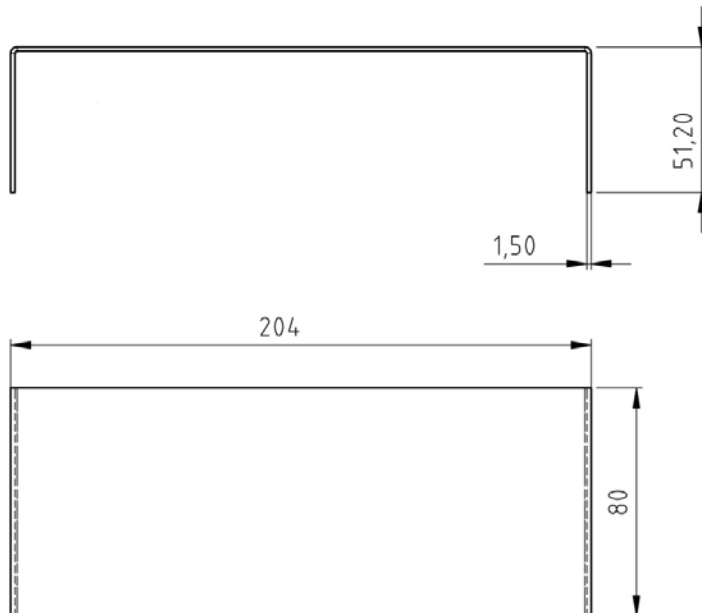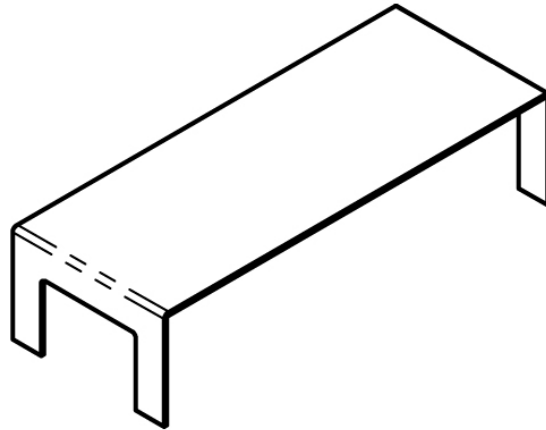
### 3 Installation with KERN EMB 2000-2V

Fig.1: KERN EMB 2000-2V with installed density sets KERN YDB-02

- ⇒ Disconnect scale from power supply.
- ⇒ Remove the standard weighing plate and replace it with the density set.
- ⇒ Place the platform for glass containers in a way that it does not touch the weighing plate.
- ⇒ Place beaker in the centre of the platform. Make sure that it has no contact with the frame.
- ⇒ Hang the immersion basket on the rack. Make sure that it is centred in the recess.
- ⇒ Pour the liquid into the glass beaker. Filling height should be approx.  $\frac{3}{4}$  of the capacity. Immerse thermometer
- ⇒ Heat the liquid, the instruments or the sinker until the temperature is constant. Observe the warm-up time of the balance.

For further information, as well as the procedure of density determination please see the operating instructions added to the density balance KERN EMB-V or on the KERN homepage ([www.kern-sohn.com](http://www.kern-sohn.com)).

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