-ebro[®]

MEASUREMENTS FOR LIFE

Winlog.light

Quick Reference Guide

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Content

Introduction	16
Survey	16
System requirements	16
Hardware requirements	16
Software requirements.....	16
Installation of Software	16
Inserting the CD-ROM	17
Starting Setup-Assistant	17
Enter user group	18
Choose installation folder	18
Start installation	19
Connecting Interface to computer	19
Solving installation problems	21
First steps with „Winlog.light“	22
Survey	22
Adjusting Programme.....	23
Preparation of validation	23
Evaluation of validation	24
Looking into the validation report.....	24

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Introduction

Survey

Winlog.light is a reporting and evaluation software. It meets the strict requirements of pharmaceutical validation, qualification and measuring tasks. The software has a programme surface which is easy to operate and offers numerous possibilities for generating meaningful reports.

System requirements

In order to operate the software on your computer without any problems, your computer has to fulfil the following requirements.

Hardware requirements

- Processor speed at least 1 GHz
- 1 GB working storage
- 300 MB free hard disk storage
- USB (Universal Serial Bus)

Software requirements

Operating system Microsoft Windows XP, VISTA or 7

Installation of Software

Winlog.light has its own installation programme, the Setup-Assistant, which guides you in several steps through the installation of the software. Therefore the installation of the software is very easy:

- insert the CD-ROM
- start Setup-Assistant
- enter user group
- choose installation folder
- start installation

Installation of Software

In the following the separate installation steps are briefly described. In case of problems the Setup-Assistant gives you corresponding hints with suggested solutions. Be sure to observe the messages of the Setup-Assistant on the monitor!

After the installation of the software you just have to connect the interface (write-reading device).

Inserting the CD-ROM

- Insert the CD-ROM into the respective drive of your computer.

Starting Setup-Assistant

If the autostart function is switched on on your computer, the Setup-Assistant will start automatically after a few moments. If the autostart function is switched off on your computer, you have to start the Setup-Assistant manually.

- In this case select „**Start: Execute**” from the Windows tool bar
- Now select the drive where the CD-ROM is (e.g.,D:“) and directly behind it type „**setup.exe**“.

*In case it is necessary, you can cancel the installation any time. Just click on the button „**Cancel**“ and confirm the query of the Setup-Assistant with „**Yes**“.*

- In order to continue with the installation, click on the button „**Next**“. After clicking on „**OK**“ the Setup-Assistant starts. Now the Setup-Assistant will guide you through the installation of the software.

Installation of Software

In case of problems, the Setup-Assistant gives you corresponding hints with suggested solutions. Be sure to observe the messages of the Setup-Assistant on the monitor!

Enter user group

You can install the software in such a way that only you have an access to the software. If you are system administrator, you can also indicate in installation that all users of the computer are permitted to have access to the software. After this decision, the Setup-Assistant asks you.

Presetting: All users

Choose installation folder

You can install the software in any folder you want. The plant has already asserted the folder

C:\Program Files\Winlog.light

- **In order to choose another, already existing folder**, click on „**Browse**“ and then choose the desired folder.
- **In order to indicate a not yet existing folder** and have it created by the Setup-Assistant, enter the folder to be created manually into the line „**Directory**“.

Now all necessary preparations are made.

- **In order to start the installation**, click on „**Next**“.

Installation of Software

Start installation

When all installation requirements have been fulfilled, the Setup-Assistant now installs the software completely.

In case of problems, the Setup-Assistant gives you corresponding hints with suggested solutions. Be sure to observe the messages of the Setup-Assistant on the monitor!

If everything has functioned without problems, then after some time the Setup-Assistant will give the message that the software was successfully installed and that the installation is now completed.

After the installation of the software you only have to connect the interface (write-/reading device) (see next page).

Connecting Interface to computer

Write-/reading devices (Interfaces) are connected with the computer over the USB (Universal Serial Bus). In order to connect your interface with the computer, proceed as follows:

- Make sure that the installation CD is in the computer drive. This contains control programmes (so called „drivers“) which are necessary for the operation of the interfaces.
- In case you are using an fourfold/eightfold interface, connect this over the included net adapter to the power supply and switch it on (switch lights up).
- Connect the interface with the help of the included USB cable to the computer.

The operating system automatically recognizes whether there are modifications to the USB and gives corresponding messages on the monitor screen.

Connecting Interface to computer

*The operating system treats the fourfold/eightfold interface like four/eight separate hardware devices. The message „**New hardware found**“ therefore appears 8/16 times subsequently for the fourfold/eightfold interface.*

Shortly after the message „**New hardware found**“ the „**Assistant for searching for new hardware**“ appears on the monitor screen automatically.

- Select „**Automatically install Software**“ and click on „**Next**“.
- Ignore the message „**Hardware installation**“ and click on „**Continue Installation**“.

Now the „**Assistant for searching for new hardware**“ installs the driver.

- If the installation was successful, click on „**Complete**“.

*In case you have connected the fourfold/eightfold interface, the procedure „**New hardware found**“ runs through another 7/15 times up to „**Complete**“.*

To solve installation problems

In case problems should arise with the installation of the drivers, a warning message appears.

In this case proceed as follows:

Select **Start** → **Settings** → **Control panel** → **Administrative tools** → **Computer management**.

The window „**Computer management**“ will open.

- Click there on „**Device Manager**“.

Now a list of the hardware devices contained in the computer appears on the right. All hardware devices which pose problems for the operating system are marked with a yellow symbol.

- If you also discover a COM port with a yellow symbol on your computer, double click on the respective COM port name.

The window „**Characteristics of USB Serial Port**“ will open.

- Click on the button „**Install Driver anew**“.

Now the „**Hardware-Update-Assistant**“ opens and offers to install the driver.

- Proceed as described with the „**Assistant for searching for new hardware**“. The now necessary steps are identical with those described there.

First steps with „Winlog.light“

Survey

After the double click on the programme symbol, the programme main window will open. This has two main areas: the area for the function selection (Navigator) and the working space.

The procedures of the programmes are subdivided into three areas: „**Validations**“, „**Preferences**“ and „**Reports**“.

In the **area „Validations“** validation plans can be drawn up, process data can be evaluated according to a validation procedure or validation reports can be opened for viewing.

The **area „Preferences“** offers entry possibilities for

- general information (e.g. folder for the source database, existent interfaces [write-/reading devices] for the data loggers or physical units to be used),
- the security (e. g. password, notification email addresses or electronic signature),
- the user administration (who may use the programme and with which rights?) as well as
- the devices used (e.g. type of devices, manufacturer, device programme).

In the **area „Reports“** the recorded programme tasks can be browsed, filtered and shown as a report according to various features under „**Audit Trail**“.

First steps with „Winlog.light“

Adjusting Programme

After the installation the programme needs certain basic information in order to work in accordance with the regulations. But also afterwards in the current use of the programme, it may be necessary to change certain preferences.

Adjusted can be

- general information as for example the folder of the database, the interface (write-/reading device), the units of measurements to be used or information on the users firm.
- Information on security such as passwords, protection from misuse, email addresses for automatic notifications or electronic signature;
- for users with access to the programme and their user rights;
- the hardware devices used (hardware device manufacturer, type of device etc.).

Please obtain further details from the included Help on the monitor.

Preparation of Validation

The programme helps you to prepare a validation respective plan one. It guides you in six subsequent masks (steps) through the preparation. In these masks it demands all necessary details to be given one after the other. After having finished the details you can view the validation plan shown and have it printed out.

Please obtain further details from the included Help on the monitor.

First steps with „Winlog.light“

Evaluation of validation

The programme helps you with the evaluation of a validation. After expiry of a validation cycle, it confirms whether a validation was successful or not according to the validation plan and the data from the data loggers. For this the programme selects the data from the data loggers and judges this according to the details given from the validation plan.

Each validation only lets itself be evaluated once. A validation report is made from the validation plan by the evaluation. This is available any time after the evaluation and can be looked into.

The programme guides you in three subsequent masks (steps) through the evaluation.

Please obtain further details from the included Help on the monitor.

Looking into the validation report

With the help of a report over a validation, each validation also allows itself to be judged at a later time and for example in the case of a dispute to be proven (Quality Assurance).

The programme directly files all data and details which are created in preparing a validation or in the evaluation of a validation, into the corresponding database. In calling up a validation plan or a validation report, the programme pick outs the necessary data from the database and puts them into a prefabricated report form.

Reports of already evaluated validations are saved in databases belonging to the programme secure from manipulation. They can be viewed at any time. In doing so the data is not changed.

You can print the report directly or e.g. export it into Portable Document Format (PDF).

Please obtain further details from the included Help on the monitor.

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ebro Electronic GmbH & Co. KG