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Ziegelei 1
D-72336 Balingen
E-Mail: info@kern-sohn.com

Tel: +49-[0]7433- 9933-0
Fax: +49-[0]7433-9933-149
Internet: www.kern-sohn.com

Zusatzbeschreibung Supplementary instructions Instructions supplémentaires

KERN BFB_IFB_KFx-TM_SFB_UFx

Version 1.0
02/2010

Deutsch
English
Français

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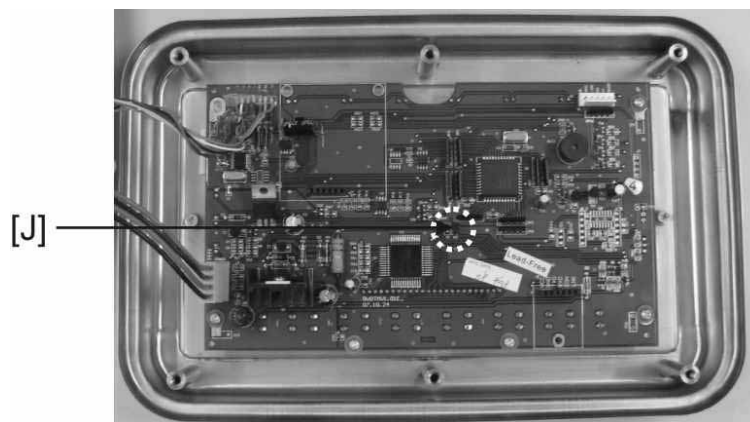
Supplementary instructions

1 Adjustment

As the acceleration value due to gravity is not the same at every location on earth, each display unit with connected weighing plate must be coordinated - in compliance with the underlying physical weighing principle - to the existing acceleration due to gravity at its place of location (only if the weighing system has not already been adjusted to the location in the factory). This adjustment process must be carried out for the first commissioning, after each change of location as well as in case of fluctuating environment temperature. To receive accurate measuring values it is also recommended to adjust the display unit periodically in weighing operation.

- Access to menu block "P2 CAL" / "P3CAL" is locked in verified weighing systems.

KERN BFB, IFB, KFB-TM, UFB

Access to menu requires breaking of seal and activation of adjustment switch (See chpt 2).

KERN SFB, KFN-TM, UFN

In order to access the menu you will have to short-circuit the two contacts of the circuit board with a jumper (See chpt 2).

Attention:

The instrument will have to be reverified by a bureau of standards and a new seal attached if a seal was broken before applications subject to verification can be recommenced.

Please follow verification instructions (See chpt 2.).

- The applied adjustment weight depends on the weighing system's capacity. Carry out adjustment as near as possible to the maximum weight of the weighing system. Info about test weights can be found on the Internet at: <http://www.kern-sohn.com>
- Observe stable environmental conditions. Warm-up time required for stabilisation.

1.1 KERN BFB, IFB, KF_x-TM, UF_x, SFB verified models:

- ⇒ Switch on instrument and press  during self-test. Pn
- ⇒ Press , ,  in turn and the first menu block “PO
CHK” will appear. POCHK
- ⇒ Press  repeatedly until “P2 CAL” appears. P2CAL
- ⇒ Press  and the first menu item “COUNT” will appear.
KERN BFB, IFB, KFB-TM, UFB:
At verified scales press first the adjustment switch. Count
- ⇒ Press  repeatedly until “CAL” appears. CAL
- ⇒ Press . UnLd
- ⇒ Wait for stability display, then press . Ensure that there are no objects on the weighing plate. STABLE
UnLd
- ⇒ The currently set adjustment weight will be displayed. 15.000 kg
- ⇒ Change by pressing the navigation keys (See manual chpt 2.1.1); select desired setting; currently enabled digit will be flashing. STABLE
LoAd
- ⇒ Acknowledge by . PASS
- ⇒ Carefully place adjusting weight in the centre of the weighing plate. Wait for stability display, then press . PASS
- ⇒ After the adjustment the balance will carry out a self-test. Remove adjusting weight **during** self test, balance will return into weighing mode automatically. In case of an adjustment error or incorrect adjusting weight the display will show an error message, repeat adjustment process. STABLE
ZERO
GROSS
0.000 kg

1.2 KERN SFB non-verifiable models:

- ⇒ Switch on instrument and press  during self-test. Pn
- ⇒ Press , ,  in turn and the first menu block “POCHK” will appear. POCHK
- ⇒ Press  repeatedly until “P3 CAL” appears. P3CAL
- ⇒ Press  and the first menu item “COUNT” will appear. CoUnt
- ⇒ Press  repeatedly until “CAL” appears. CAL
- ⇒ Press . If necessary press  repeatedly until “nonLin” appears. noLin
- ⇒ Press . Ensure that there are no objects on the weighing plate. UnLd
- ⇒ Wait for stability display, then press . STABLE
UnLd
- ⇒ The currently set adjustment weight will be displayed. ,15.000 kg
- ⇒ Change by pressing the navigation keys (See manual chpt 2.1.1); select desired setting; currently enabled digit will be flashing. STABLE
LoAd
- ⇒ Acknowledge by . PASS
- ⇒ Carefully place adjusting weight in the centre of the weighing plate. Wait for stability display, then press . STABLE
GROSS
0.000 kg
- ⇒ After the adjustment the balance will carry out a self-test. Remove adjusting weight **during** self test, balance will return into weighing mode automatically. In case of an adjustment error or incorrect adjusting weight the display will show an error message, repeat adjustment process.

2 Verification

General introduction:

According to EU directive 90/384/EEC balances must be verified if they are used as follows (legally controlled area):

- e) For commercial transactions if the price of goods is determined by weighing.
- f) For the production of medicines in pharmacies as well as for analyses in the medical and pharmaceutical laboratory.
- g) For official purposes.
- h) For manufacturing final packages.

In cases of doubt, please contact your local trade in standard.

Verification notes:

An EU Qualification Approval is in existence for verified weighing systems. If a balance is used where obligation to verify exists as described above, it must be verified and re-verified in regular intervals.

Reverification is carried out according to relevant national statutory regulations. The validity for verification of balances in Germany is e.g. 2 years.

The legal regulation of the country where the balance is used must be observed!

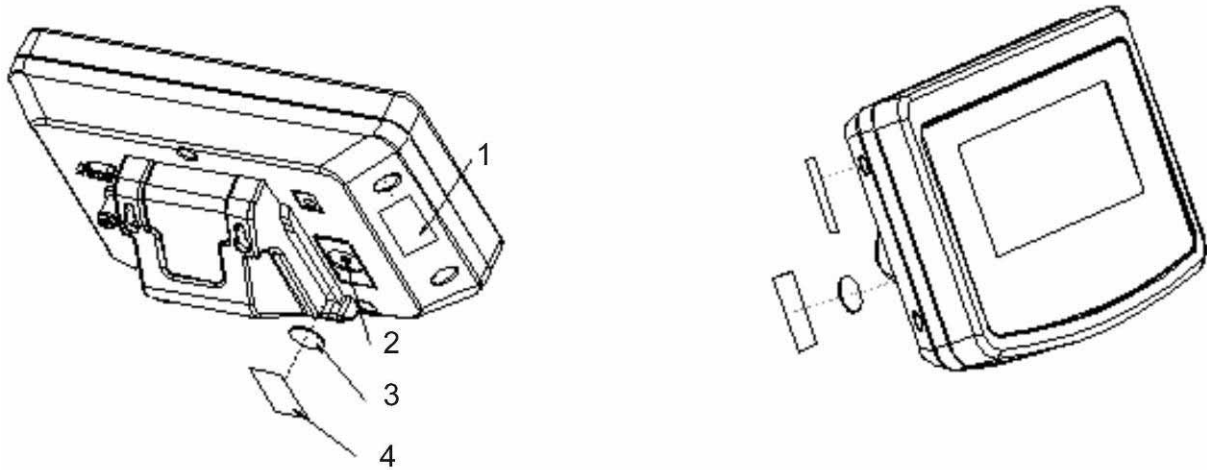
- Verification of the weighing system is invalid without the "seal".

Notes on verified weighing systems

KERN BFB, IFB, KFB-TM, UFB:

Position for seals and adjustment switch

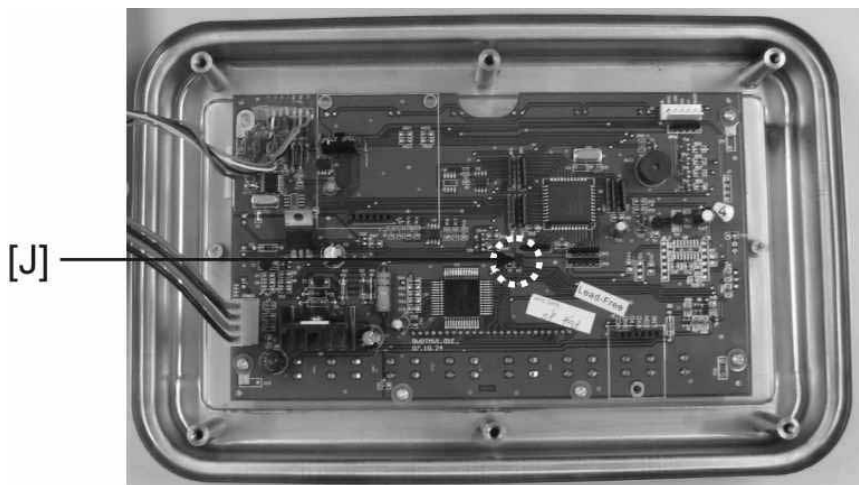
- 5. Self-destructing seal
- 6. Adjustment switch
- 7. Cover adjustment switch
- 8. Self-destructing seal

KERN SFB, KFN-TM, UFN:

Access to conductor plate:

- Remove seal
- How to open the display unit
- For adjustment / access the menu, the jumper [J] has to be set on both pins as shown in the picture.

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