MIKROSKOP MICROSCOPE 40X-1024X

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EN Operating instructions

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General Information

About This instruction Manual

Please read the safety instructions in this manual carefully. To avoid damage to the unit and the risk of injury, only use this product as described in the manual. Keep the instruction manual handy so that you can easily look up information on all the functions.

CAREFUL!

You will find this symbol before every section of text that deals with the risk of minor to severe injuries resulting from improper use.

Intended Use

This product is intended only for private use.

It was developed for magnified viewing of natural and man-made objects.

General Warning

DANGER!

Tools with sharp edges and points are often used when working with this device. Because there is a risk of injury from such tools, store this device and all tools and accessories in a location that is out of the reach of children.

DANGER!

This device contains electronic components which operate via a power source (power supply and/or batteries). Only use the device as described in the manual, otherwise you run the risk of an electric shock.

DANGER!

Do not expose the device to high temperatures. Use only the supplied power supply or the recommended batteries. Do not short-circuit the device or batteries or throw them into a fire! Excessive heat or improper handling could trigger a short-circuit, a fire or an explosion. Never bend, pinch or pull the power and connecting cables, extensions and adapters. Protect the cables from sharp edges and heat. Before operating, check the device, cables and connections for damage. Never use a damaged unit or a unit with damaged power cables. Damaged parts must be exchanged immediately by an authorised service centre.

DANGER!

Children should only use the device under adult supervision. Keep packaging material, like plastic bags and rubber bands, out of the reach of children, as they pose a risk of choking.

CAREFUL!

Children must not have access to the included chemicals and liquids. Do not drink the chemicals. Wash hands thoroughly with running water after using the chemicals. In the event that the chemicals come into contact with your eyes or mouth, rinse thoroughly with water. If you are in pain after exposure, contact a doctor immediately and show him the substances.

NOTE!

Do not disassemble the device. In the event of a defect, please contact your dealer. The dealer will contact the Service Centre and can send the device in to be repaired, if necessary.

Do not expose the device to temperatures above 45 °C.

Operating instructions

Parts overview:

- ❶ 10X WF eyepiece
- ❷ 16X WF eyepiece
- ❸ Barlow lens
- ❹ Electronic eyepiece (MicroOcular)
- ❺ Microscope head
- ❻ Objective revolver
- ❼ Microscope stage
- ❽ Focus wheel
- ❾ LED lighting (transmitted light)
- ❿ Electricity supply
- ⓫ Microscope base
- ⓬ Main plug
- ⓭ 5 slides, 10 covering glasses and 5 preparations in a plastic box
- ⓮ Colour filter disc
- ⓯ LED lighting (reflected light)
- ⓰ Direct light/transmitted light switch
- ⓱ Microscope tools: a) pipette; b) tweezers
- ⓲ Hatchery
- ⓳ MicroCut
- ⓴ Specimens: a) yeast; b) gum media (specimen inclusion medium); c) sea salt; d) brine shrimp eggs
- ⓵ Locking screw
- ⓶ Photomizer Software

1. General/Location

Make sure you position your microscope on a stable, solid surface. An electricity supply is required for observation with the electric illuminator.

Position your device so that it can be disconnected from the power supply at any time. The wall you use socket should be located near the device and easily accessible, since the plug on the power cord serves as a disconnecting device for the power supply. Always pull on the plug to separate the device from the power supply. Never pull on the cord.

2. Electric LED lighting

Before use, make sure the light switch (16) is set to 'off'.

The microscope has two light sources. Lighting can be of three types. Set the switch (16) to 'II' to light the specimen from above (reflected light) or 'I' to light it from below (transmitted light). Use setting 'III' to light the specimen simultaneously by transmitted and reflected light. The transmitted light unit (9) is used for transparent specimens (those on glass slides). To view solid, non-transparent specimens, use the reflected light unit (15). Use of both forms of lighting simultaneously is only recommended for semitransparent specimens. This operating mode is not recommended for transmitted light specimens on slides as it may cause reflection on the slide.

3. Colour filter disc

The colour filter (14) under the microscope table (7) aids in viewing very bright and transparent objects. Just select the right colour for the specimen in question. The components of colourless or transparent objects (e.g. starch particles, single-cell specimens) can thus be better recognised.

4. Microscope setup

The microscope head (5) will now be prepared for your first observation. First, loosen the screw (21) and rotate the head into a convenient position. Begin every observation with the lowest magnification. Place the microscope's table (7) with the focus knob (8) into the lowest position and rotate the objective revolver (6) until it locks on the lowest magnification (4X).

**NOTE:**

Make sure to place the microscope's table (7) in its lowest position before changing the objective in order to prevent damage to the microscope.

Insert the 10X eyepiece (No. 1, 1) in the Barlow lens (No. 1, 3).

Take care that the Barlow lens is inserted completely into the monocular head (No. 1, 4).

5. Observation

After you have set up the microscope with the proper illumination, the following principles are important:

Begin each observation at the lowest magnification, so that the centre and position of the object to be viewed is in focus. The higher the magnification, the more light is required for good picture quality.

Place a permanent slide culture directly under the microscope lens on the plate (7). The specimen to be examined must be directly over the lighting.

Look through the eyepiece (1 and 2) and carefully turn the focus wheel (8) until you can see a sharp picture.

Now you can progress to a higher magnification. Slowly pull the Barlow lens (No. 2, 3) out of the monocular barrel (No. 2, 4). When the Barlow lens is nearly entirely pulled out, the magnification is raised to 2X.

For even higher magnification, you can put the 16X eyepiece (2) into the objective revolver (6) and rotate the objective revolver to a higher magnification (10X or 40X).

**TIP:**

Depending on the preparation, higher magnifications do not always lead to better pictures.

When changing the magnification of your microscope by changing or adjusting the eyepiece, objective lens or Barlow lens, you must readjust the focus wheel (8) to sharpen the image.

**NOTE:**

Please be very careful when doing this. If you move the mechanical plate upward too fast, the objective lens and the slide can touch and become damaged.

6. Condition and prepare viewed objects

6.1. Condition

With the Barlow lens nearly fully extended, your microscope's magnification can be doubled. Both transparent and non-transparent specimens can be examined with this microscope, which features both direct and transmitted light. If opaque specimens are being examined, such as small animals, plant parts, tissues, stones and the like, the light is reflected from the specimen, through the lens and eyepiece, where it is magnified, to the eye (reflected light principle, switch position I). If opaque specimens are being examined, the light from below goes through the specimen, lens and eyepiece to the eye and is magnified en route (direct light principle, switch position II).

Some small water organisms, plant parts and animal components are transparent by nature, but many others require pretreatment — that is, you need to make a thinnest possible slice of the object by hand cutting or using a microtome, and then examine this sample.

6.2. Creation of thin preparation cuts

Specimens should be sliced as thin as possible. A little wax or paraffin is needed to achieve the best results. Put the wax into a heat-safe bowl and heat it over a flame until the wax is melted. You can use a candle flame to melt the wax.

**DANGER!**

Be extremely careful when dealing with hot wax, as there is a danger of being burned.

Then, dip the specimen several times in the liquid wax. Allow the wax that encases the specimen to harden. Use a MicroCut (19) or other small knife or scalpel to make very thin slices of the object in its wax casing.

**DANGER!**

Be extremely careful when using the MicroCut, knife or scalpel. These instruments are very sharp and pose a risk of injury.

Place the slices on a glass slide and cover them with another slide before attempting to view them with the microscope.

6.3. Creation of your own preparation

Put the object to be observed on a glass slide and cover the object with a drop of distilled water (No. 3) using the pipette (No. 3, 17a).

Set a cover glass (available at a well-stocked hobby shop) perpendicular to the edge of the water drop, so that the water runs along the edge of the cover glass (No. 4). Now lower now the cover glass slowly over the water drop.

**TIP:**

The gum medium (20b) supplied is used to make permanent slide cultures. Add it instead of distilled water.

The gum medium hardens so that the specimen is permanently affixed to its slide.

7. Experiments

Now that you're familiar with your microscope's functions and how to prepare slides, you can complete the following experiments and observe the results under your microscope.

7.1. Newspaper print

Objects:

1. A small piece of paper from a newspaper with parts of a picture and some letters
 2. A similar piece of paper from an illustrated magazine:
- Use your microscope at the lowest magnification and make a slide preparation from each object. Place the slide with the newspaper on the microscope table and observe the slide. The letters in the newspaper appear broken because the newspaper is printed on raw, inferior paper. Now observe the slide with the magazine preparation. Letters of the magazine appear smoother and more complete. The picture from the newspaper consists of many small points, which appear somewhat dirty. The pixels (raster points) of the magazine image appear sharper.

7.2. Textile fibres

Objects and accessories:

1. Threads of different textiles: Cotton, linen, wool, silk, Celanese, nylon and any others you can find.
2. Two needles:

Put each thread on a glass slide and fray each with the help of the two needles. Put a drop of water over each thread with the pipette and cover each with a cover glass. Adjust the microscope to a low magnification. Cotton fibres are of plant origin and look, under the microscope, like a flat, twisted band. The fibres are thicker and rounder at the edges than in the centre. Cotton fibres consist primarily of long, collapsed tubes. Linen fibres are also of plant origin; they are round and run in straight lines. The fibres shine like silk and exhibit numerous swellings along the shaft of the fibre. Silk is of animal origin and consists of solid fibres of smaller diameter than the hollow vegetable fibres. Each silk fibre is smooth and even and has the appearance of a small glass rod. Wool fibres are also of animal origin; the surface consists of overlapping scales, which appear broken and wavy. If possible, compare wool fibres from different weaving mills, and note the differences in the appearance of the fibres. Experts can determine the country of origin of wool based on its appearance under a microscope. Celanese is artificially manufactured by a long chemical process. All Celanese fibres show hard, dark lines on a smooth, shining surface.

The fibres crinkle in the same way after drying. Observe the similarities and differences between the different fibres.

7.3. Saltwater brine shrimps

Accessories:

1. Brine shrimp eggs (20d)
2. Sea salt (20c)
3. Hatchery (18)
4. Yeast (20a)

CAREFUL!

These eggs are not fit for human consumption.

7.3.1. Winter eggs of *Artemia salina*

Artemia salina are species of brine shrimp typically found in salt lakes – bodies of water with a higher salinity than even the ocean. During a drought, a salt lake can become a hostile habitat for organisms, and entire populations of *Artemia salina* sometimes die off. During drought conditions, to ensure that the species will repopulate the salt lake when the drought ends, *Artemia salina* lay thick-shelled eggs called winter eggs that can survive for up to ten years in a dormant state. Winter eggs can withstand heat, cold and chemicals. These eggs hatch when favourable conditions return to their ambient environment. The eggs provided (20d) are of this type.

7.3.2. Hatching winter eggs

To hatch the brine shrimps, create a solution with an appropriate salinity and temperature. First, fill two containers with a half litre of freshwater each, and let them both stand for about thirty hours. Next, pour half of the provided salt (20c) into one container and stir the solution until the salt dissolves. Pour some of this solution into the hatchery (18). Place a few eggs close to the lid. Position the hatchery somewhere with plenty of light but not in direct sunlight. The ambient temperature should ideally hover around 25 °C. As water in the hatchery evaporates, gradually

add freshwater from the second container. After two to three days, the eggs will hatch brine shrimp larvae, called nauplii.

7.3.3. Observing *Artemia salina* under a microscope

Using the pipette (17a), move some larvae from the container to a microscope slide for examination. When viewing the larvae, you'll notice that they swim through the solution using hairlike limbs! Each day, examine a few more. You can even view the entire hatchery under the microscope if you remove its lid. The larvae will mature in six to ten weeks, depending on the temperature of the water. Soon, you will have an entire generation of saltwater brine shrimps that reproduce frequently!

7.3.4. Feeding your *Artemia salina*

Feed your brine shrimps often to keep them alive. The best food is dry powdered yeast (20a). Give them some every other day. Be careful not to overfeed them, as doing so can cause the water to stagnate and poison the brine shrimps. If the water does begin to stagnate (you'll see it darkening), transfer the brine shrimps to the fresh saline solution you have prepared earlier (see 7.3.2).

8. MicroOcular setup

NOTE:

The MicroOcular only works without the Barlow lens supplied. Magnification setting is changed by using the MicroOcular and must be reset by re-focussing.

Remove the Barlow lens (Fig 5, 3) and eyepiece currently in use from the eyepiece holder (Fig 5, 5) and replace them with the MicroOcular (PC-Ocular) (Fig 6, 4) and reducer lens (Fig 6, F) as in illustration 6 in the holders (Fig 6, 5).

**NOTE:**

Please do not yet connect MicrOcular and PC. Please go through the items below in sequence.

9. Installation and use of the software

9.1. Software and installation information

A software CD is included with your microscope. The software and drivers on this CD must be installed on your computer in order to use your MicrOcular. Once you connect the MicOcular to your computer after installation, you can view pictures on and save them to your computer using Photomizer. To install the software and driver correctly, simply follow the installation steps.

9.2. Software installation

1. **Important:** Before inserting the CD, first plug the USB cable into the USB port on your PC. Windows will now recognize the new device, and will indicate this in a notification window. Now please click on "Cancel".
2. Now insert the CD-ROM that came in your package into the CD/DVD drive of your computer. The installation menu starts automatically. If it does not, go to the Windows explorer and select the CD/DVD drive (most of the time, it's the "D" drive, but it can have another letter). From there, start the file "autorun.exe" by double-clicking with the left mouse button.

9.2.1. Driver installation

To install the driver software, click on the menu point "Install Driver" with the mouse cursor. Then follow the installation program instructions. During the software installation the correct driver for your operating system will be installed automatically. No manual input is needed. In rare cases the device may not be recognised by your computer. As a rule you need then only install the driver manually from the CD. If this fails please refer to the troubleshooting chapter that follows.

9.2.2. Installation of the Image Editing Software Photomizer

The image editing software "Photomizer" is located on the software CD. You can edit your pictures here.

1. To install the software, click once on the menu point "Install PHOTOMIZER" with the left mouse button.
2. The Photomizer Software requires Microsoft .NET Framework 4.0, which it will install if it is not already on your system. If it is already on your system, you may skip to step 5.
3. In the welcome window, please accept the Microsoft license agreement, and then click "Install". The installation can take a few minutes.
4. As soon as everything is installed, click "Finish".
5. Now you will be presented with a choice, in which you can choose your language.
Make your selection and confirm it by clicking „OK“.
6. When you see "Welcome", click on „Next“.
7. In the next window, you will be asked for the „Destination Folder“. Here, just click on „Next“.
8. Now, the window with the Setup-Status will appear – here, a progress bar will inform you about the ongoing installation. This process can take a few minutes.
9. The window "Photomizer is being installed" appears. Click on "Finish". The installation ends.

**TIP:**

To use the MicOcular long-term we recommend to always connect it to the same USB port.

10. Using the MicOcular

10.1. Preparation

1. Slide a specimen under your microscope and focus on it.
2. Remove the eyepiece and Barlow lens from the eyepiece support and the dust cap from your MicOcular and install same in the eyepiece supports instead of the Barlow lens.
3. Start your PC if you haven't yet and connect your MicOcular to the USB port of your computer.

10.2. Showing and saving MicOcular images on your PC

1. Start the Photomizer Software.
2. Click „Open camera“
3. In the event that you have connected more than one unit, you can choose the desired unit in the subsequent selection. Click on "USB 2.0 Webcam". If only one device is connected, this step is omitted.
4. The camera image should now be visible on your screen. Focus the microscope image.
5. Click „Capture“ to record an image. It will then be shown on the right.
6. Click on it to select it and then click „Transfer image“.
7. This takes you to the Photomizer software.
8. File - Save as

10.3. The Photomizer Software

If you need help in the „Photomizer“ programme click „?“ and then „Open help“. If you need further assistance please visit the maker's home page at www.photomizer.net/bresser

Care and maintenance

Before cleaning, separate the device from the power supply by removing the plug. Only use a dry cloth to clean the exterior of the device.

**NOTE:**

Do not use any cleaning fluid to avoid damaging the electronics.

Clean the lenses eyepieces and lenses only with a soft, lint-free cloth, like a microfibre cloth.

**NOTE:**

Do not apply excess pressure to the cloth so as to avoid scratching the lenses.

To remove more stubborn dirt, moisten the cleaning cloth with an eyeglass-cleaning solution and wipe the lenses gently. Protect the device from dust and moisture. After use, particularly in high humidity, let the device acclimatize for a short period of time, so that the residual moisture can dissipate before storing.

Troubleshooting

Problem

No picture visible

Picture flickers while viewing with PC eyepiece

Solution

- switch light on
- adjust focus

- if necessary, adjust resolution of the video graphics board

Software installation reports
 “not XP approved”

- confirm by clicking “OK”

System requirements for PC eyepiece

Minimum system requirements: PC with an Intel Pentium IV processor or higher; Windows XP with Service Pack 3*, Windows Vista (32/64Bit) with Service Pack 2* or Windows 7 (32/64Bit) with Service Pack 1*; .NET Framework 4.0*; min. 1024 MB RAM (64Bit = 2048 MB); min. 500 MB free hard drive space; free USB port; CD/DVD/BD drive.

*available for free via Windows Update (Internet connection needed)

Magnification table

<i>Eyepiece</i>	<i>Objective</i>	<i>Magnification</i>	<i>with Barlow lens</i>
10X	4X	40X	64X
10X	10X	100X	160X
10X	40X	400X	640X
16X	4X	64X	102X
16X	10X	160X	256X
16X	40X	640X	1024X

DISPOSAL

 Dispose of the packaging materials properly, according to their type, such as paper or cardboard. Contact your local waste-disposal service or environmental authority for information on the proper disposal.

 Do not dispose of electronic devices in the household garbage!
 As per Directive 2002/96/EC of the European Parliament on waste electrical and electronic equipment and its adaptation into German law, used electronic devices must be collected separately and recycled in an environmentally friendly manner.

 In accordance with the regulations concerning batteries and rechargeable batteries, disposing of them in the normal household waste is explicitly forbidden. Please make sure to dispose of your used batteries as required by law — at a local collection point or in the retail market. Disposal in domestic waste violates the Battery Directive.

-  Cd¹
 -  Hg²
 -  Pb³
- ¹ battery contains cadmium
² battery contains mercury
³ battery contains lead

Declaration of Conformity

Product Type: Microscope
 Product Name: Microscope 40X-1024X
 Article No.: 90-39100

Bresser GmbH has issued a 'Declaration of Conformity' in accordance with applicable guidelines and corresponding standards. This can be viewed any time upon request.

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MIKROSKOP MICROSCOPE 40X-1280X

National Geographic's net proceeds support vital exploration, conservation, research, and education programs.

Visit our website:

www.nationalgeographic.com

© 2014 National Geographic Society
NATIONAL GEOGRAPHIC and Yellow Border Design
are trademarks of the National Geographic Society,
used under license. All rights reserved.

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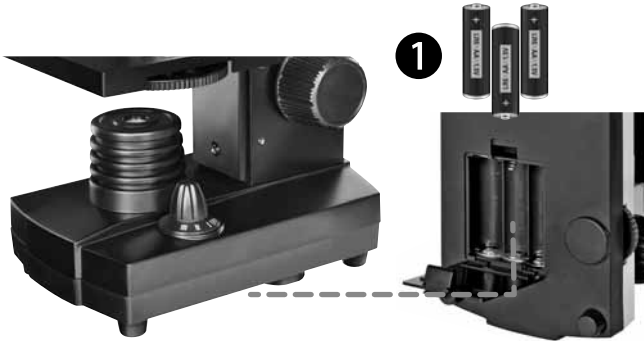
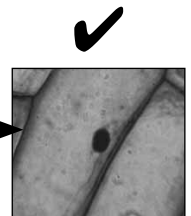
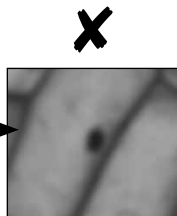
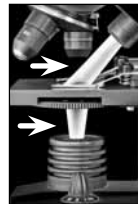
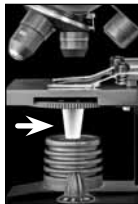
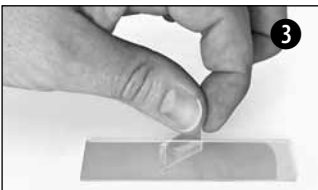
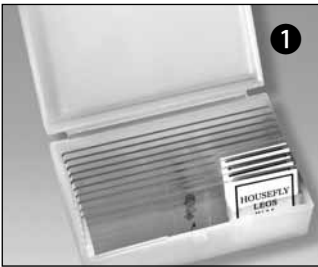
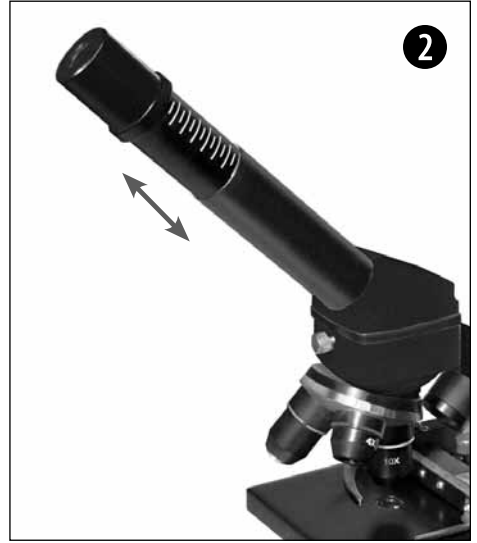
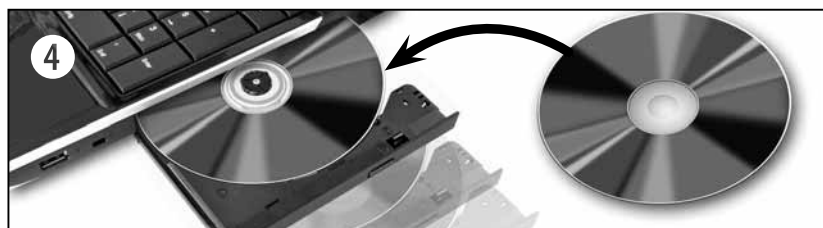
EN About this manual

This quick-start guide will help you put the device into operation quickly, but it is not intended to replace the more detailed user manual.

Please read the safety instructions and the user manual carefully before using the device. This quick-start manual contains the most important steps to start using your new device. Comprehensive operating instructions and further information, including troubleshooting advice, is available on our website: www.bresser.de (Download: <http://www.bresser.de/9039100>)

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Name: Telefon:

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General Warning

- Tools with sharp edges and points are often used when working with this device. Because there is a risk of injury from such tools, store this device and all tools and accessories in a location that is out of the reach of children.
- This device contains electronic components which operate via a power source (batteries). Only use the device as described in the manual, or you run the risk of an electric shock.
- Do not expose the device to high temperatures. Use only the power adapter supplied. Do not short-circuit the device or batteries, or throw them into a fire. Excessive heat or improper handling could trigger a short-circuit, a fire or even an explosion.
- Never bend, pinch, pull or run over the power and connecting cables, extensions and adapters. Protect the cables from sharp edges and heat. Before operating, check the device, cables and connections for damage. Never use a damaged unit or a unit with damaged power cables. Damaged parts must be exchanged immediately by an authorised service centre.
- Children should only use the device under adult supervision. Keep packaging material, like plastic bags and rubber bands, out of the reach of children, as they pose a choking hazard.
- Position your device so that it can be disconnected from the power supply at any time. The wall socket you use should be located near the device and easily accessible, since the plug on the power cable serves as a disconnecting device for the power supply.
- Do not disassemble the device. In the event of a defect, please contact your dealer. The dealer will contact the Service Centre and can send the device in to be repaired, if necessary.
- Do not expose the device to high temperatures.

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Using the Software CD

Packaged with your new microscope, you'll find a software CD that contains the driver and image-processing software. You can install the software on your computer. Then when you connect your microscope to your computer with the USB cable, you can look at the pictures from the microscope on the computer monitor, and save them using the CamApp software. You can edit the pictures with the Photomizer software. We recommend that you always connect your microscope to the same USB port.

Important: Before inserting the CD into your computer, first plug the microscope's USB cable into the USB port on your PC. Windows will now recognize the new device and will indicate this in a notification window. Click on 'Cancel', since the driver for the device will be installed automatically later. Now insert the CD that came packaged with your microscope into the CD/DVD drive of your computer. The installation menu should start automatically. If it does not, go to the Windows Explorer and select the CD/DVD drive (most of the time, it's the 'D' drive, but it can have another letter). From there, start the file 'menu.exe' by double-clicking with the left mouse button.

Driver installation

To install the driver software, click 'Install Driver' on the menu with the left mouse button. Then follow the installation program instructions. During the software installation the correct driver for your operating system will be installed automatically. No manual input is needed. Once the computer restarts, the driver is installed. In rare cases, the device may not be recognised by your computer. You should then install the driver manually from the CD. Select 'Repair Program' during the installation of the driver. After the driver has been installed, restart the computer.

Software installation

Note: The Photomizer software requires Microsoft .NET Framework 4. This must be installed first, if it has not been installed on your operating system already.

1. To install the software, click 'Install Software' on the menu once with the left mouse button.
2. A selection window will appear in which you can select the preferred language. Select your preferred language and confirm by clicking 'OK'.
3. In the welcome window, click on 'Next'.
4. In the next window, you will be asked for the destination folder. Here, just click on 'Next'.
5. The options 'Create Desktop Shortcut' and 'Link in the Quick Launch Bar' set up program shortcuts on your desktop and taskbar. Click on 'Install'.
6. Now, the window with the setup status will appear. Here, a progress bar will inform you about the ongoing installation. This process can take a few minutes.
7. The window informing you that Photomizer is being installed will appear. Click on 'Finish' to complete the installation.

Warranty and warranty term extension

The warranty term is two years from the date of purchase. Please keep your proof of purchase. Register at www.bresser.de/warranty and fill out a brief questionnaire to get your warranty term extended to **five years**. Registration must be completed within three months of purchase (date of receipt) to validate the warranty. If you register thereafter, the warranty term will not be extended. If you have problems with your device, please contact our customer service first. Do not send any products without consulting us first by telephone. Many problems with your device can be solved over the phone. If the problem cannot be resolved by phone, we will take care of transporting your device to be repaired. If the problem occurred after the warranty ended or it is not covered by our warranty terms, you will receive an estimate of repair costs, free of charge.

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NATIONAL GEOGRAPHIC and Yellow Border Design are trademarks of the National Geographic Society, used under license. All rights reserved.