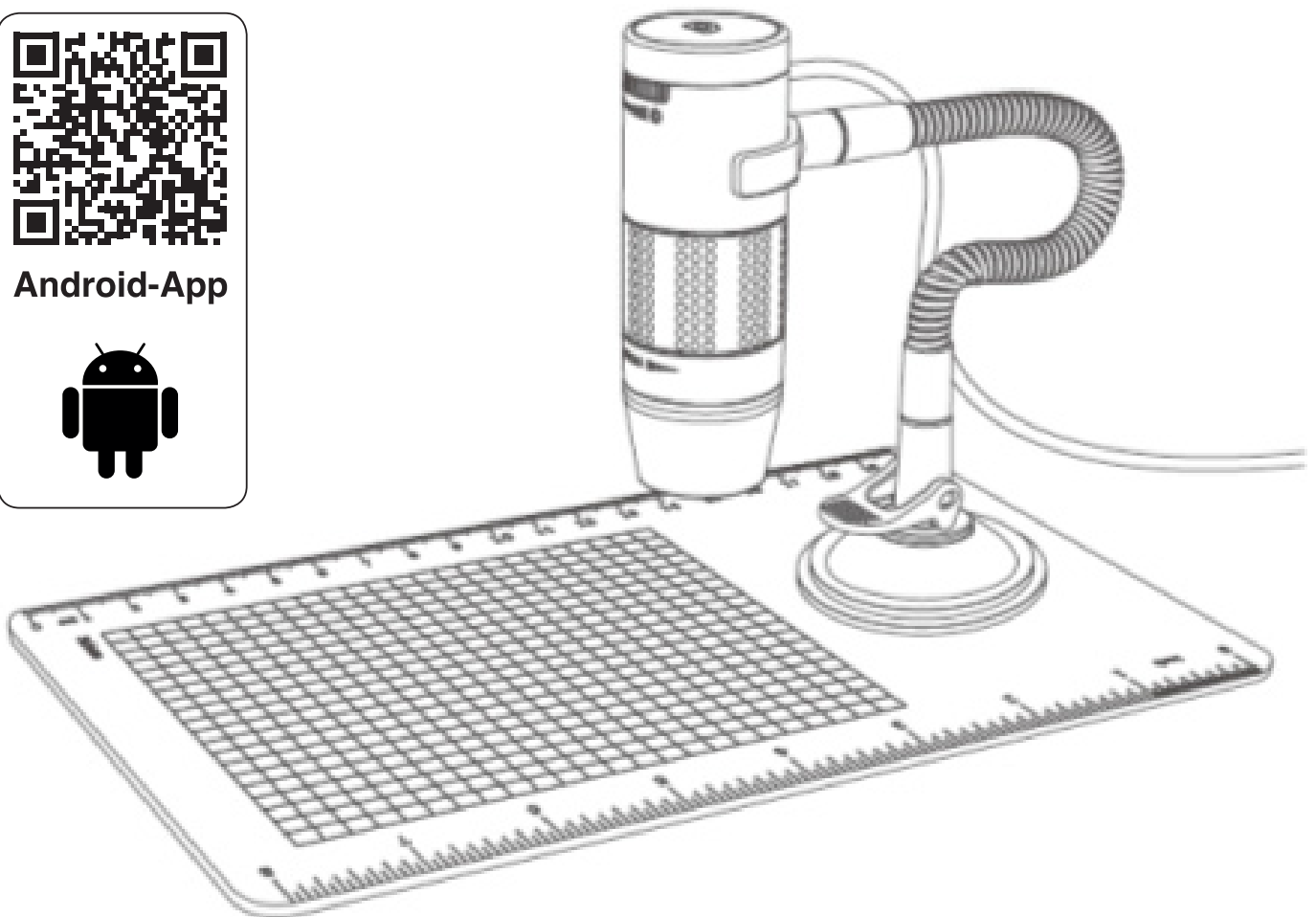
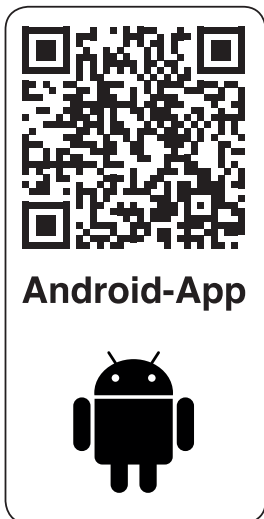
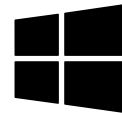
reflecta[®]

DigiMicroscope Flex

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
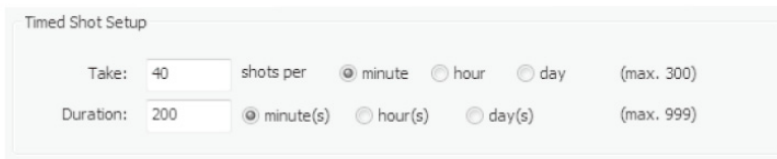
The 'Timed Shot Setup' dialog box contains two rows of settings. The first row is labeled 'Take:' and has a text input field with the value '40', followed by the text 'shots per'. To the right are three radio buttons: 'minute' (which is selected), 'hour', and 'day'. Further right is the text '(max. 300)'. The second row is labeled 'Duration:' and has a text input field with the value '200', followed by three radio buttons: 'minute(s)' (selected), 'hour(s)', and 'day(s)'. Further right is the text '(max. 999)'.

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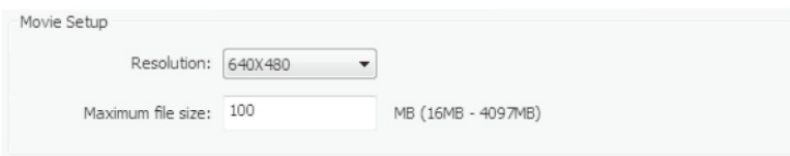
The 'Movie Setup' dialog box has two rows. The first row is labeled 'Resolution:' and has a dropdown menu showing '640X480'. The second row is labeled 'Maximum file size:' and has a text input field with the value '100', followed by the text 'MB (16MB - 4097MB)'.

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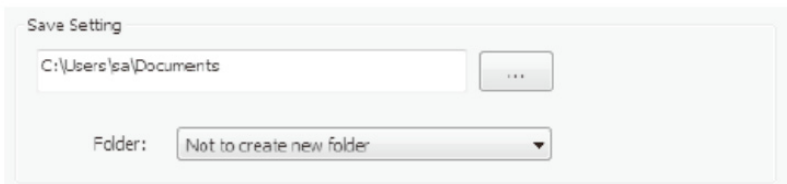
The 'Save Setting' dialog box has two rows. The first row shows a text input field with the path 'C:\Users\sas\Documents' and a button with three dots '...'. The second row is labeled 'Folder:' and has a dropdown menu showing 'Not to create new folder'.

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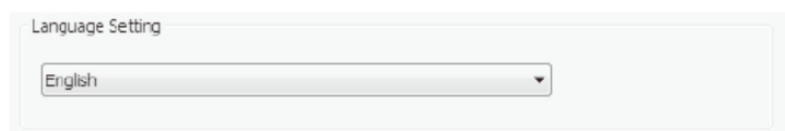The 'Language Setting' dialog box has one row with a dropdown menu showing 'English'.

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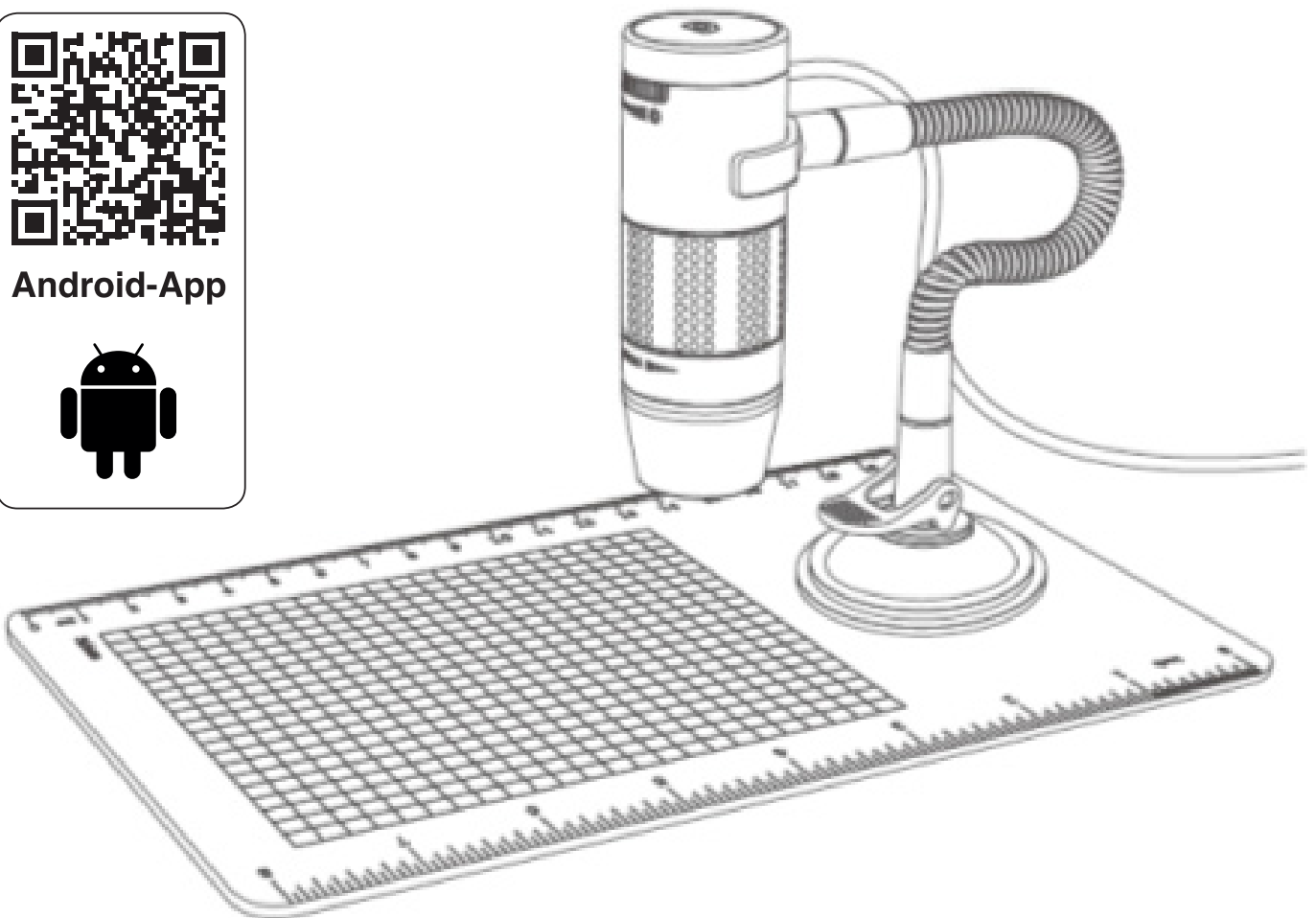
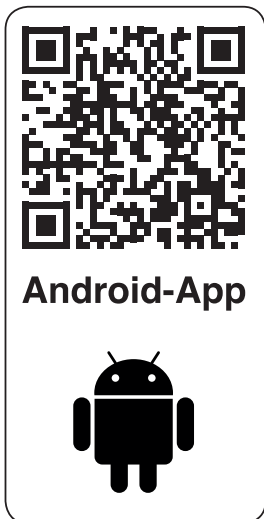
EN

CONTENTS

BEFORE USE	4	USING THE XPLOVIEW SOFTWARE	17
Important information	4	Button menu	17
Care and maintenance	4	Full screen viewing	18
Warning	4	Image rotation / flip	19
Product description	5	System settings menu	20
Computer requirements	5	Device setup	20
		Timed shot setup	21
		Movie setup	21
		Save setting	21
		Language setting	21
		Advanced settings	22
		Saved files	23
		Uninstalling the xploview software	23
PRODUCT AT A GLANCE	6		
Package contents	6		
Product overview	7		
Product specification	8		
GETTING STARTED	9		
Software installation	9		
Connecting the device	10		
Starting the xploview software	10		
Assembling the stand	11		
Attaching the microscope to the stand	12		
BASICS	12		
Focusing	12		
Preparation position	12		
Viewing using the microscope stand	13		
Viewing directly on an object	14		
For 60x magnification	14		
For 250x magnification	15		
Adjusting the LED light intensity	16		
Taking a snapshot	16		

BEFORE USE

Important information

Please read this instruction manual carefully before using this product, and retain this instruction for future reference.

Improvements and changes to this text necessitated by typographical errors, or improvements to the software and / or equipments may be made at any time without notice.

Care and maintenance

- Avoid vibration, shock and pressure e.g. dropping the microscope.
- Keep the device dry and protect it from water or vapour.
- Do not leave your device in a place with an extreme high or low temperature.
- Do not touch the device with a wet hand as it may damage the device, or cause an electric shock to the user.
- Do not use or store the device in dusty, dirty areas as its moving parts may be damaged.
- Do not use harsh chemicals, cleaning solvents or strong detergents to clean the device. Wipe it with a soft cloth slightly dampened in a mild soap-and-water solution.

Warning

- Do not place lighted digital microscope upon the eye, doing so may cause permanent eye damage.
- Do not attempt to open or dismantle the digital microscope.

Product description

- This digital microscope differs from a traditional optical microscope, where, instead of looking through an eyepiece, magnified image or live video is displayed on a computer monitor via USB connection.
- This device has 2 magnification powers, it can magnify images up to 60x and 250x when placed directly on top of the object and through a 17" monitor. Magnification power depends on the size of the monitor used - the larger the monitor, the higher the magnification.
- This device has 4 LED lights, where the intensity can be adjusted.

Computer requirements

Windows based PC

Compatible operating systems

Windows 10 (32 bit or 64 bit)
Windows 8 (32 bit or 64 bit)
Windows 7 (32 bit or 64 bit)
Windows Vista (32 bit or 64 bit)
Windows XP SP2, SP3

USB

USB 2.0

Mac OS based PC

Compatible operating systems

Mac OS X 10.5.6 - Mac OS X 10.11.x


USB

USB 2.0


PRODUCT AT A GLANCE

Package contents

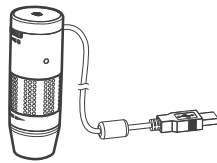Digital microscope
with USB cable

Installation CD

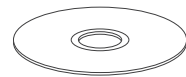
Microscope stand
with suction cup


User manual

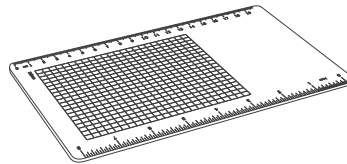Gridded
observation board

Product overview

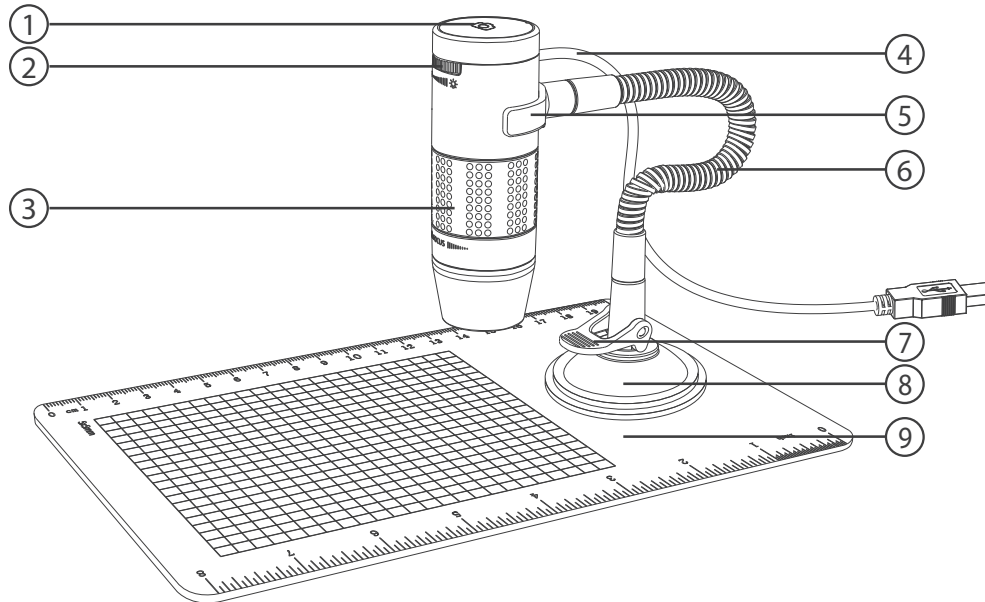1. Snapshot button
2. Light intensity knob
3. Focusing knob
4. USB cable
5. Stand bracket

6. Flexible tube
7. Locking lever
8. Suction cup
9. Observation board

Product specification

Connection type	USB 2.0
Effective magnifications (on a 17" monitor)	60x, 250x
Effective viewing area	5.6 x 4.2mm (60x), 1.4 x 1.1mm (250x),
Illumination	Four LEDs
Power supply	5V=250mA (via USB)
Sensor	CMOS
Maximum snapshot resolution	1600 x 1200 pixels (UXGA)
Maximum video capturing resolution	640 x 480 pixels (VGA)
Size	Ø 30 x 86mm
Weight	207 grams

GETTING STARTED

Software installation

Windows based PC

- Insert the supplied application CD to the CD-ROM of the computer.
- Double click the "xploview v3.2.xx.exe" icon <  > located on the driver CD.
- Follow the xploview setup wizard to install the application software for the digital microscope.

Mac OS based PC

- Insert the supplied application CD to the CD-ROM of the computer.
- Double click the "xploview v3.2.xx.dmg" icon <  > located on the driver CD.
- Drag the xploview icon <  > into the Applications folder.

Connecting the device

Connect the device to the computer using the USB cable provided.

Starting the xploview software

Windows based PC

The xploview software can be launched by double clicking the xploview icon <  > from the desktop, or from the start menu.

Mac OS based PC

The xploview software can be launched by double clicking the xploview icon <  > from the Applications menu.

Note

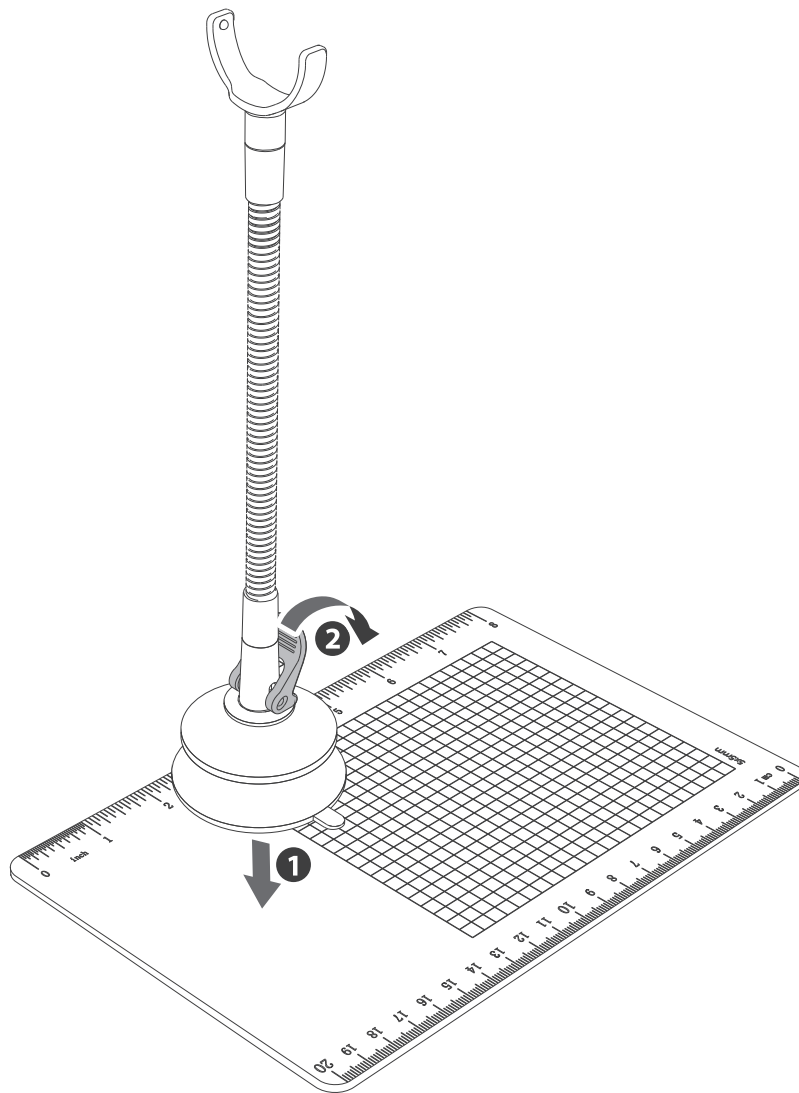
When the device is connected to the computer for the first time, a driver will automatically be installed by Windows or Mac OS. This process could take up to a few minutes.

Each time when the xploview software is started, the device should start up automatically. Otherwise, the device can be selected from the device setup section of the system settings menu.

It is normal that the LED will flash once when the device is plugged in. When xploview is running, the LED will switch on again automatically.


Assembling the stand

- Make sure the mounting surface is free from oil and dirt, and ensure the locking lever is released.
- Press the suction cup down and secure it into place by pushing the locking lever down.

Attaching the microscope to the stand

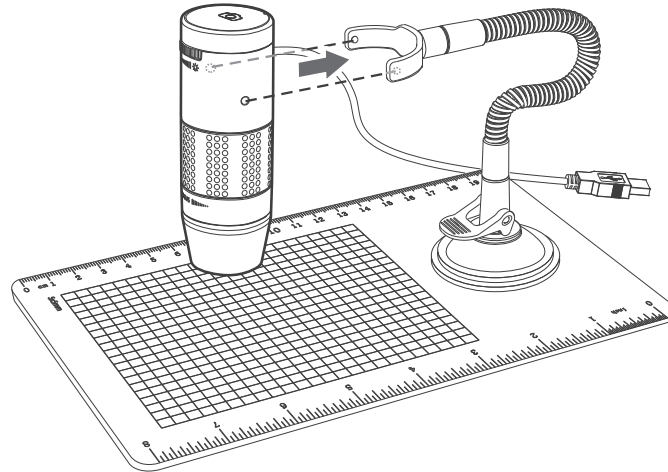Secure the microscope to the stand as shown below.


BASICS

Focusing

Preparation position



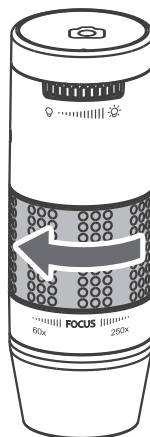
Turn the focusing knob clockwise to the furthest position.

Viewing using the microscope stand

Place the object on the observation board.

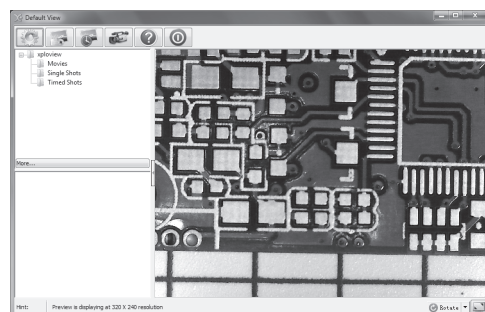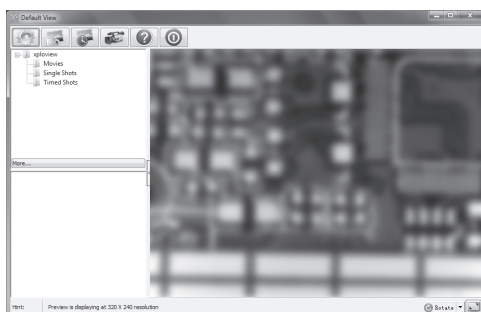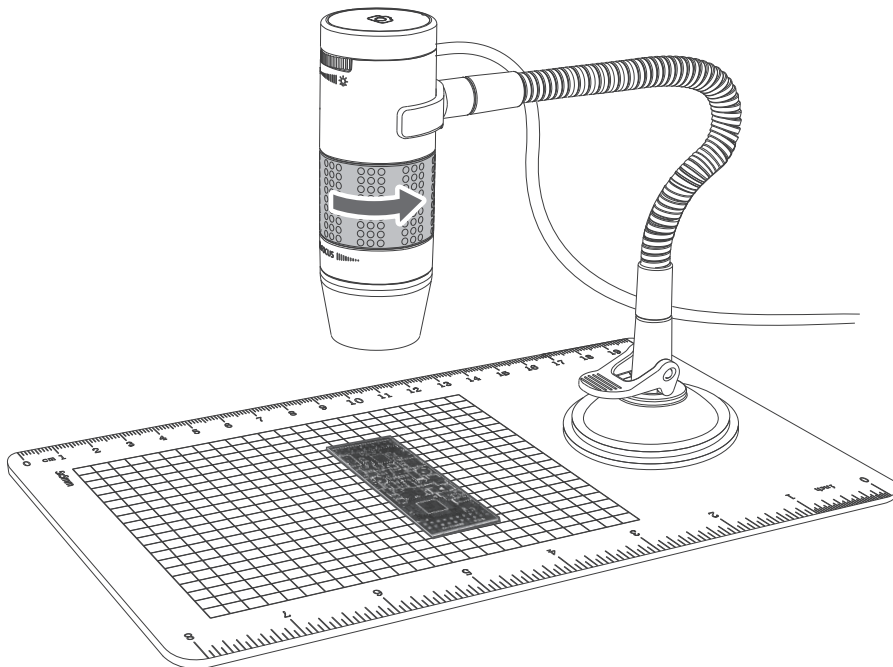From the preparation position, turn the focusing knob anticlockwise until the image is sharp and clear. The closer the microscope to the object, the higher the magnification power.



Viewing directly on an object

For 60x magnification (17" monitor)



Place the microscope directly on top of an object .

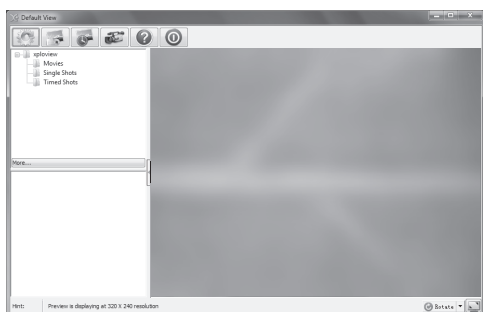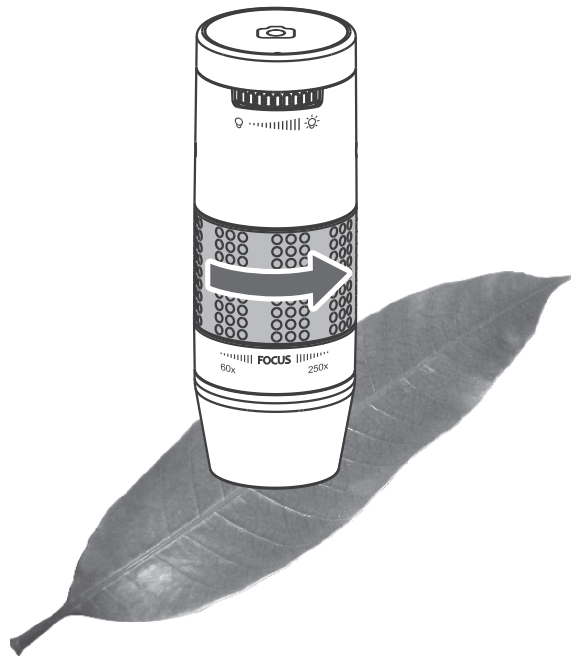From the preparation position, turn the focusing knob anticlockwise until the first focal point has been reached where the image is sharp and clear.
Magnification of 60x can be achieved when viewing on a 17" monitor.


For 250x magnification (17" monitor)

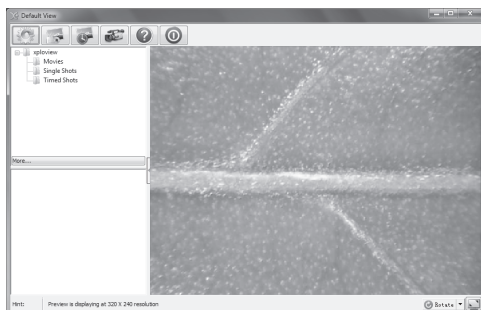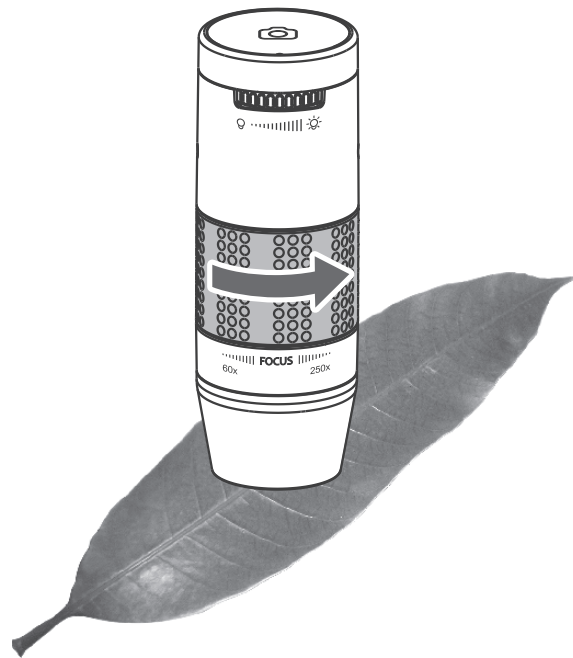At 60x magnification, turn the focusing knob further in anticlockwise direction until the second focal point has been reached where the image is sharp and clear. Magnification of 250x can be achieved when viewing on a 17" monitor.


Adjusting the LED light intensity

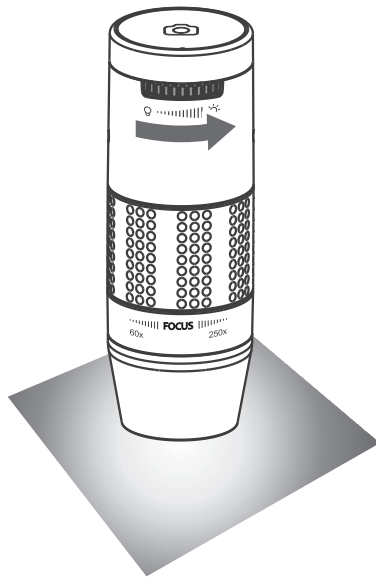Turn the light intensity knob clockwise or anticlockwise to increase or decrease the LED light intensity.


Increasing the LED light intensity

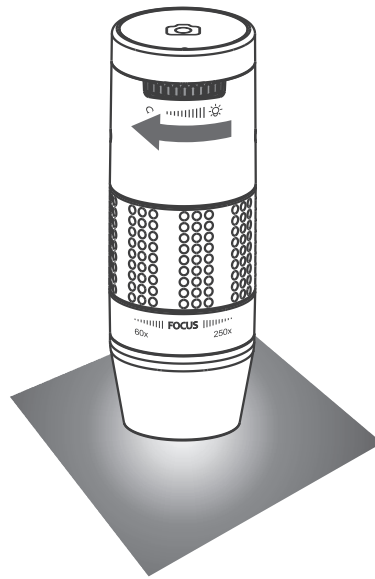Reducing the LED light intensity

Taking a snapshot

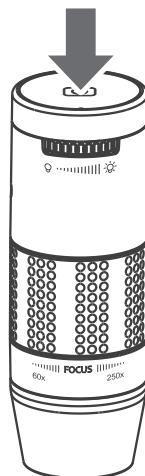The snapshot button is touch-sensitive. Tap the snapshot button located on top of the device to take a snapshot.

USING THE XPLOVIEW SOFTWARE

Button menu

The icons on the button menu :

Open **System Settings Menu** (see system settings menu on page 20).

Capture on screen image.

To start and to stop **Timed Shot**. Images will be captured at a regular interval (see system settings menu on page 21 to adjust frequency and duration).

To start and to stop **Video Recording**.

Information of the application software. This information maybe helpful when updating software.

Shut down the application software.

Full screen viewing



To activate the full screen mode, click the full screen button <  > located on the bottom right corner of the xpoview application software window.

To exit full screen mode, either double click on the screen, or press the "Esc" button on the keyboard.


Image rotation / flip

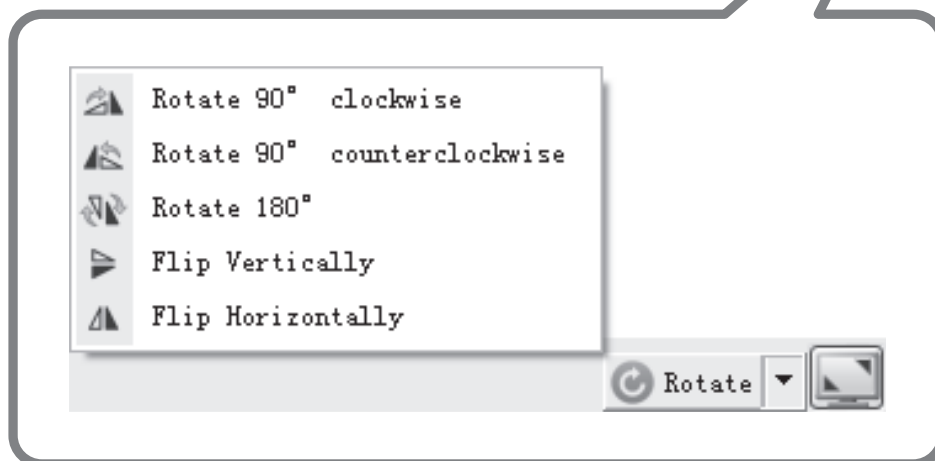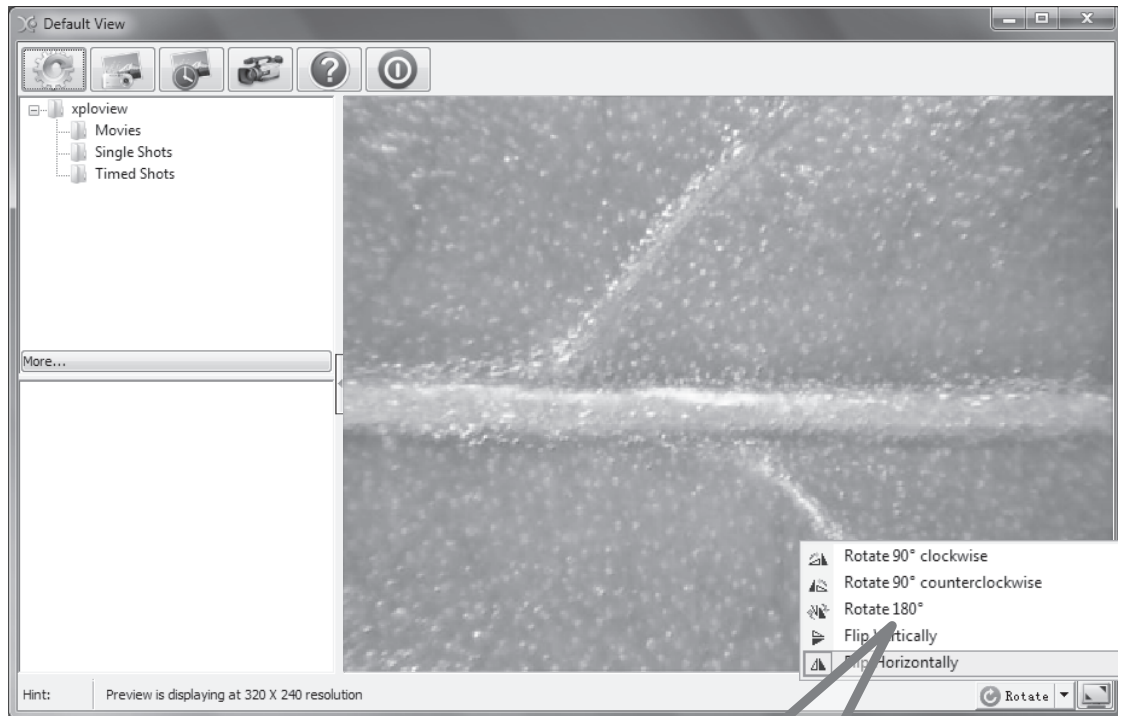Click <  Rotate > button to rotate or flip the image.

System settings menu

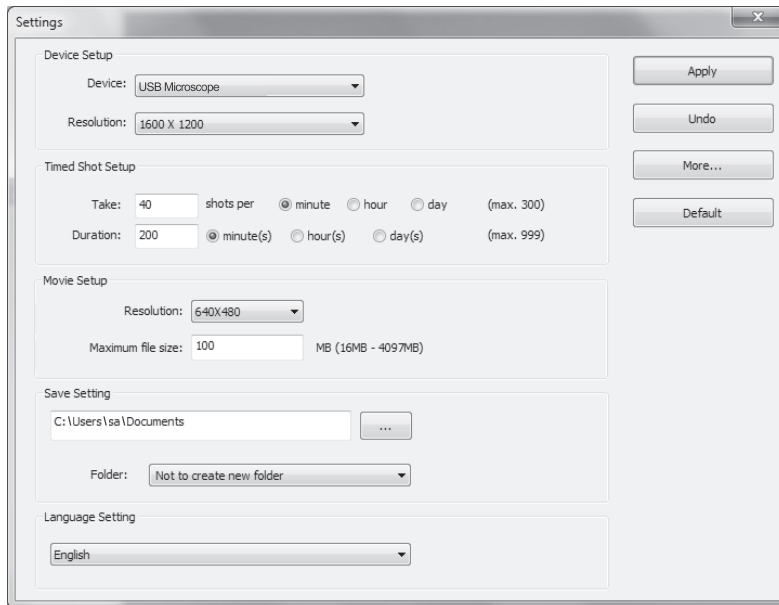
The first time the xploview software is started, the default settings will be loaded, you may change these settings manually in the system settings menu.

Device setup


If the image captured by the digital microscope was not displayed by default, you can change this by selecting it from the "Device" drop-down menu.

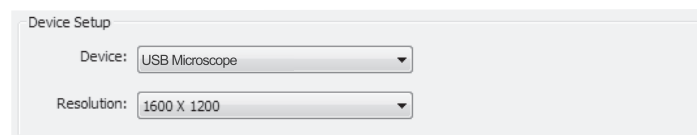The resolution of images you capture can be changed from the "Resolution" drop-down menu.

Timed shot setup

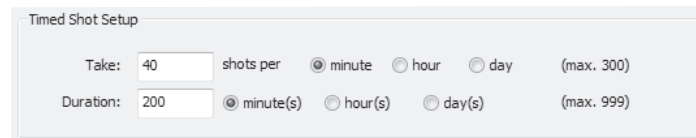The frequency and duration of automatic images capture can be adjusted under this option.

The screenshot shows the 'Timed Shot Setup' dialog box. It contains two rows of controls. The first row has a 'Take:' label, a text input field with '40', the text 'shots per', and three radio buttons labeled 'minute', 'hour', and 'day'. The 'minute' radio button is selected. To the right of these controls is the text '(max. 300)'. The second row has a 'Duration:' label, a text input field with '200', and three radio buttons labeled 'minute(s)', 'hour(s)', and 'day(s)'. The 'minute(s)' radio button is selected. To the right of these controls is the text '(max. 999)'.

Movie setup


The resolution of videos you record can be changed from the "Resolution" menu. You can also set a maximum file size for each video.

The screenshot shows the 'Movie Setup' dialog box. It contains two rows of controls. The first row has a 'Resolution:' label, a dropdown menu showing '640X480', and a small downward arrow. The second row has a 'Maximum file size:' label, a text input field with '100', and the text 'MB (16MB - 4097MB)'.

Save setting

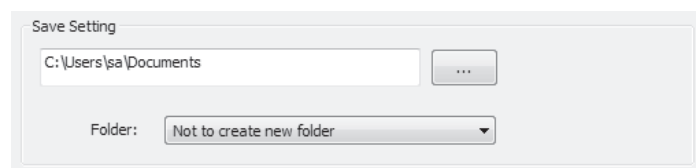
The default location for captured images or videos can be changed under this option.

The screenshot shows the 'Save Setting' dialog box. It contains two rows of controls. The first row has a text input field with 'C:\Users\sa\Documents' and a button with three dots. The second row has a 'Folder:' label and a dropdown menu showing 'Not to create new folder'.

Language setting

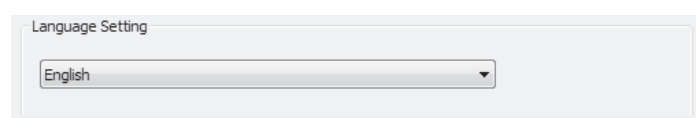The language of the xpoview software can be changed under this option.


The screenshot shows the 'Language Setting' dialog box. It contains one row of controls with a dropdown menu showing 'English'.


Advanced settings

By clicking the "More..." button on the right of the system settings menu, you will be able to manually adjust all of the image settings. Note that the settings available may be different, depending on your operating system.

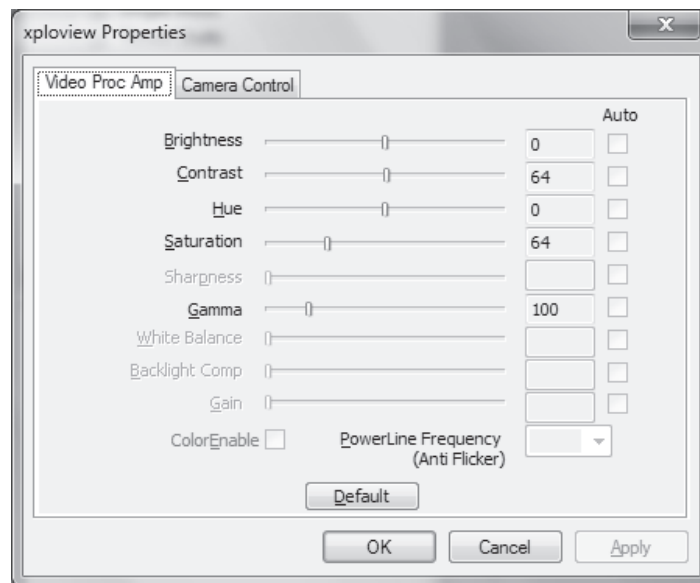Windows based PC


Mac OS based PC

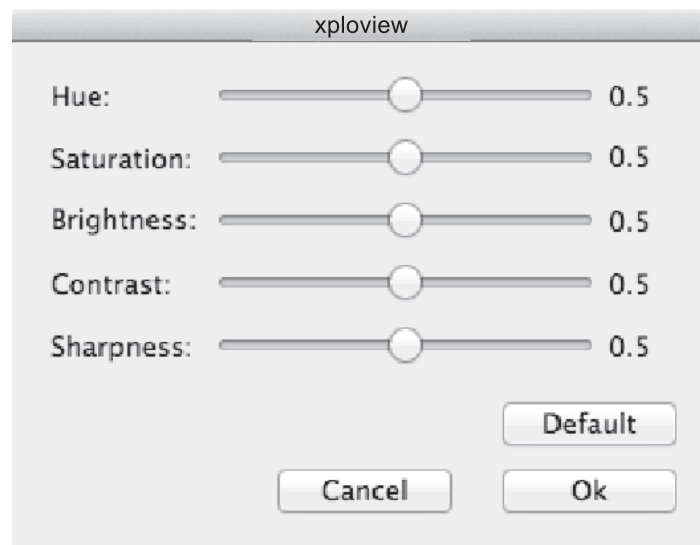Saved files

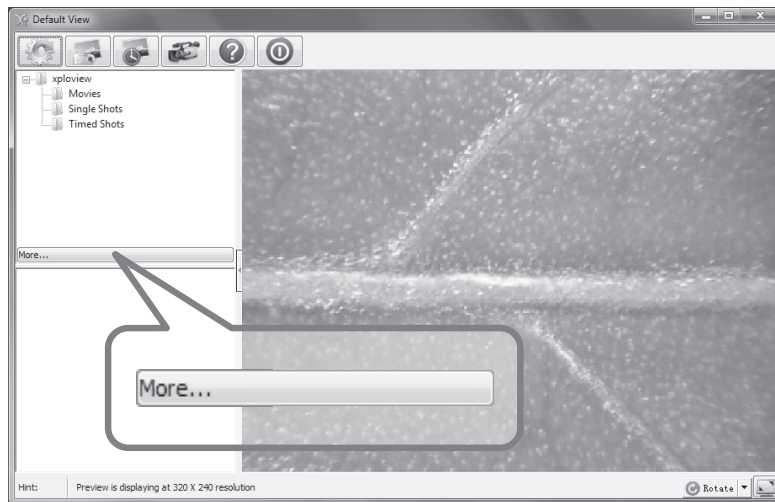With the xploview application software opened, you can locate the saved files folder by clicking the "More..." button located on the left of the main software window.

Uninstalling the xploview software

Windows based PC

Select "Uninstall" from the start menu (Start > All Programs > xploview > Uninstall).

Mac OS based PC

Drag the xploview application icon from the "Applications" folder to "Trash".

FCC compliance statement (United States only)

This device complies with part 15 of the FCC Rules. Operation is subject to the following two conditions: (1) This device may not cause harmful interference, and (2) this device must accept any interference received, including interference that may cause undesired operation.

Changes or modifications not expressly approved by the party responsible for compliance could void the user's authority to operate this device.

This device has been tested and found to comply with the limits for a Class B digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This device generates, uses and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this device does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- Reorient or relocate the receiving antenna.
- Increase the separation between the equipment and receiver.
- Connect the device into an outlet on a circuit different from that to which the receiver is connected.
- Consult the dealer or an experienced radio/TV technician for help.

Legal information

This document is published without any warranty. While the information provided is believed to be accurate, it may include errors or inaccuracies. In no event shall the manufacturer or its distributors be liable for incidental or consequential damages of any nature, including but not limited to loss of profits or commercial loss, arising out of the use of the information in this document.

All rights reserved. Mac, Mac OS and OS X are trademarks of Apple Inc., registered in the U.S. and other countries. Windows is a registered trademark of Microsoft Corporation in the United States and other countries. All other trademarks and brands are property of their respective owners.

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
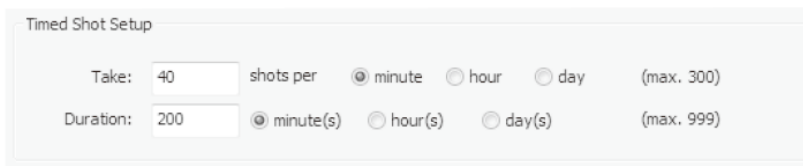
The screenshot shows the 'Timed Shot Setup' window. It contains two rows of settings. The first row is 'Take: 40 shots per' with radio buttons for 'minute' (selected), 'hour', and 'day', and a '(max. 300)' label. The second row is 'Duration: 200' with radio buttons for 'minute(s)' (selected), 'hour(s)', and 'day(s)', and a '(max. 999)' label.

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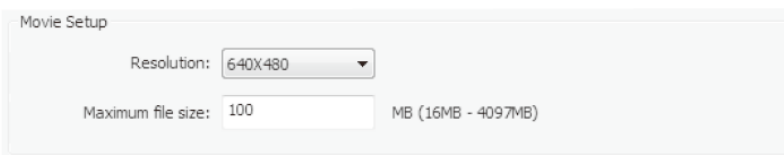
The screenshot shows the 'Movie Setup' window. It has two main settings: 'Resolution: 640x480' with a dropdown arrow, and 'Maximum file size: 100 MB (16MB - 4097MB)' with a text input field.

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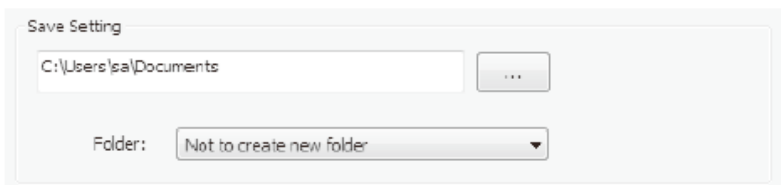
The screenshot shows the 'Save Setting' window. It features a text input field containing 'C:\Users\sa\Documents' with a browse button ('...') to its right. Below this is a 'Folder:' dropdown menu currently set to 'Not to create new folder'.

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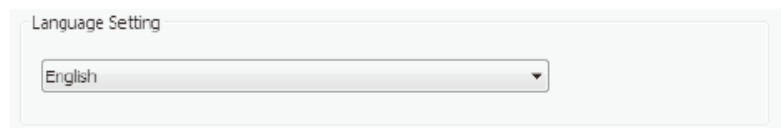The screenshot shows the 'Language Setting' window. It contains a single dropdown menu with 'English' selected.

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
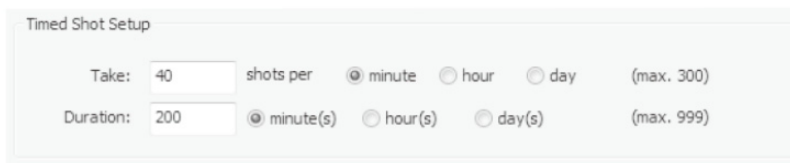
The screenshot shows the 'Timed Shot Setup' dialog box. It has two rows of controls. The first row is labeled 'Take:' and contains a text input field with the value '40', followed by the text 'shots per'. To the right are three radio buttons labeled 'minute', 'hour', and 'day'. The 'minute' radio button is selected. To the right of these is the text '(max. 300)'. The second row is labeled 'Duration:' and contains a text input field with the value '200'. To the right are three radio buttons labeled 'minute(s)', 'hour(s)', and 'day(s)'. The 'minute(s)' radio button is selected. To the right of these is the text '(max. 999)'.

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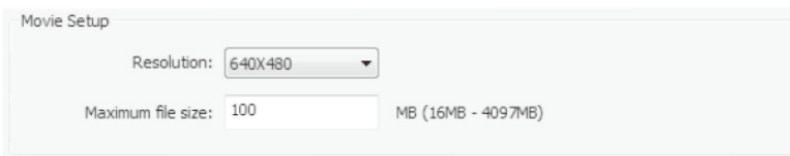
The screenshot shows the 'Movie Setup' dialog box. It has two rows of controls. The first row is labeled 'Resolution:' and contains a dropdown menu with the value '640X480'. The second row is labeled 'Maximum file size:' and contains a text input field with the value '100', followed by the text 'MB (16MB - 4097MB)'.

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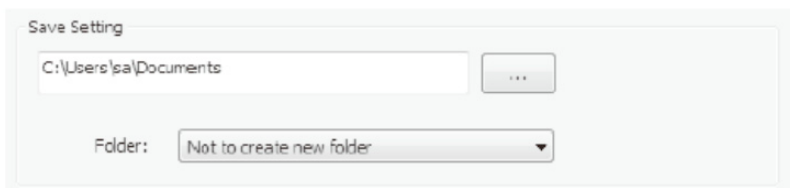
The screenshot shows the 'Save Setting' dialog box. It has two rows of controls. The first row contains a text input field with the value 'C:\Users\sa\Documents' and a button with three dots '...'. The second row is labeled 'Folder:' and contains a dropdown menu with the value 'Not to create new folder'.

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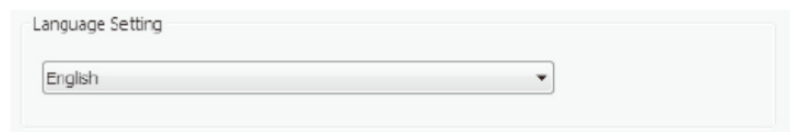The screenshot shows the 'Language Setting' dialog box. It has one row of controls containing a dropdown menu with the value 'English'.

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