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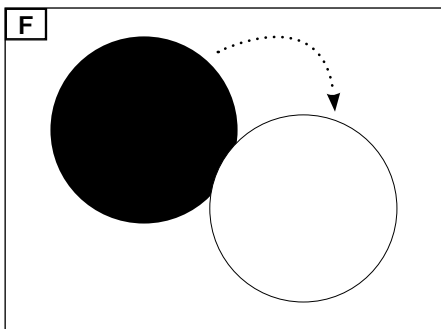
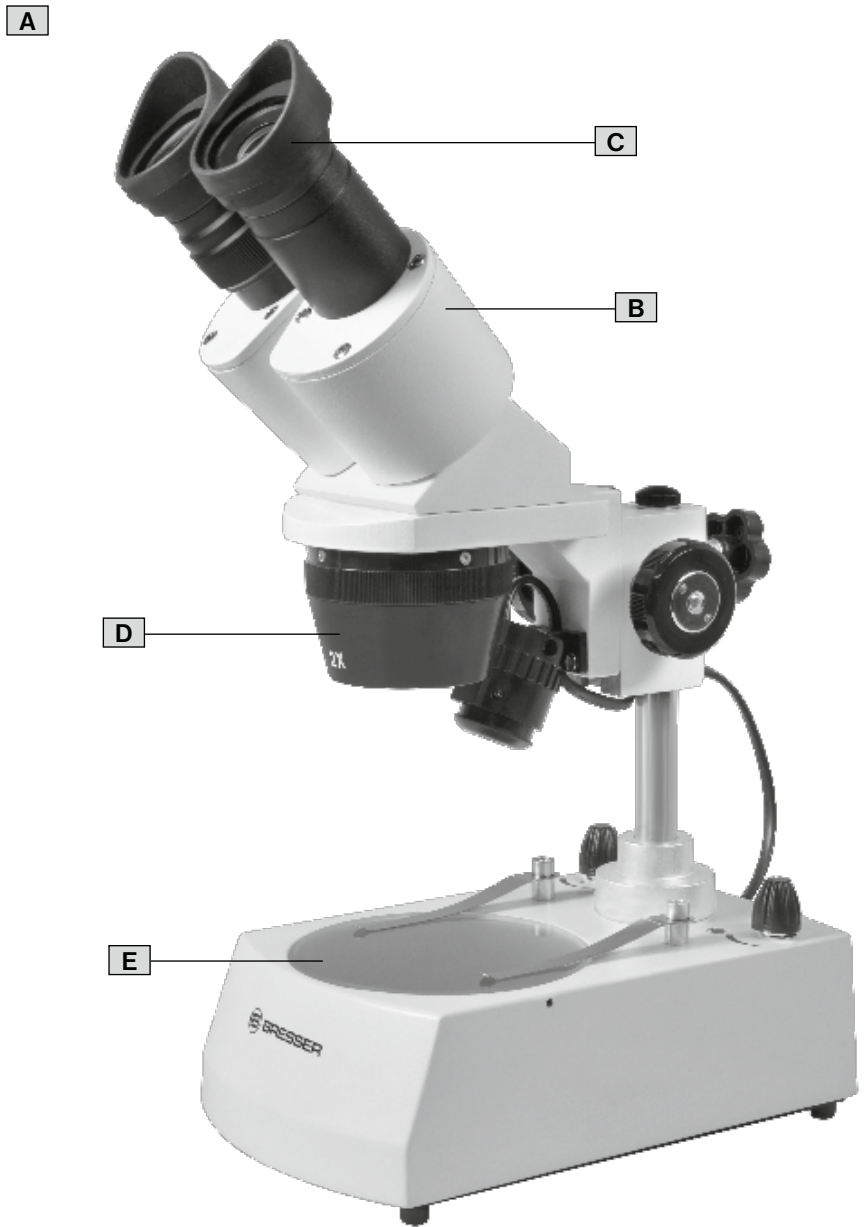
GB

ABOUT THIS INSTRUCTION MANUAL

These operating instructions are to be considered a component of the device. Please read the safety instructions and the operating instructions carefully before use. Keep these instructions for renewed use at a later date.

When the device is sold or given to someone else, the instruction manual must be provided to the new owner/user of the product.

INTENDED USE

This product is intended only for private use. It was developed for the magnified display of things in nature.

This device is not intended for use by individuals (including children) with limited physical, sensory or mental capabilities or those lacking in experience and/or knowledge, unless they are supervised by an individual responsible for their safety or have received instructions from them regarding the use of the device.

Children must be supervised to ensure that they do not play with the device.

GENERAL WARNING

CHOKING HAZARD!

Keep packaging material, like plastic bags and rubber bands, out of the reach of children, as these materials pose a choking hazard. There is a RISK OF CHOKING.

RISK of BODILY INJURY!

Tools with sharp edges and points are often used when working with this device. Because there is a risk of injury from such tools, store this device and all tools and accessories in a location that is out of the reach of children.

RISK of ELECTRIC SHOCK!

This device contains electronic components that operate via a power source (batteries). Children should only use the device under adult supervision. Only use the device as described in the manual; otherwise, you run the risk of an electric shock.

Never bend, pinch or pull the power and connecting cables, extensions and adapters. Protect the cables from sharp edges and heat. Before operating, check the device, cables and connections for damage. Never use a damaged unit or a unit with damaged power cables. Damaged parts must be exchanged immediately by an authorised service centre.

RISK of FIRE/EXPLOSION!

Do not expose the device to high temperatures. Use only the recommended batteries. Do not short-circuit the device or batteries, or throw them into a fire. Excessive heat or improper handling could trigger a short-circuit, a fire or an explosion.

RISK of property damage!

Protect the device from severe shocks!

The manufacturer is not liable for damage related to improperly installed batteries! Do not disassemble the device. In the event of a defect, please contact your dealer. The dealer will contact the Service Centre and can send the device in to be repaired, if necessary.

1. PACKAGE CONTENTS (FIG. 1)

- (A) Microscope with
 - (B) Microscope head
 - (C) 1 Pair WF10x eyepieces
 - (D) 2x/4x objectives
 - (E) Object plate, transparent and one-sided matte
 - (F) Object plate, one-sided black and one-sided white
- (G) Smartphone holder
- (H) Transport case

2. PARTS OVERVIEW (FIG. 2)

- 1 Eyepiece
- 2 Rubber eyecups
- 3 Diopter setting
- 4 Binocular head
- 5 Fixing knob
- 6 Focusing wheel
- 7 Turnable interchangeable objective
- 8 Rotary knob (On/Off switch with dimmer for transmitted light)
- 9 Rotary knob (On/Off switch with dimmer for incident light)
- 10 Slide clamp
- 11 Object plate
- 12 Battery compartment
- 13 Locking screw for battery compartment
- 14 clamp screw for the eyepiece
- 15 Smartphone holder

3. LOCATION

Before you set up the microscope, you must choose a suitable location. First you must make sure that your microscope is on a stable and solid surface.

4. INSERTING BATTERIES (FIG. 3)

The battery compartment (12) is located at the bottom of the microscope. Loosen the locking screw (13) of the battery compartment cover. Remove the cover. Insert three AA type batteries (Mignon LR6). Make sure that the poles of the batteries are in the correct position as marked on the inner side of the battery compartment. Re-insert the battery compartment cover and tighten the locking screw (13) only hand-tight to avoid any damage to the cover.

You can also use accumulators of the same type instead of normal batteries. Due to the lower voltage of accumulators, a lower illumination can occur.

5. ELECTRIC ILLUMINATION (FIG. 4)

The microscope is equipped with a battery-powered LED illumination. Therefore the usage with an off-grid power supply (e.g. outside) is possible.

On/Off switch and dimmer are combined in two different rotary switches (8/9) for each light source (transmitted light and incident light). To power on the device turn the rotary switch (8/9) until you hear a 'click'. Starting at this point, the brightness can be adjusted. Always use a light intensity that allows for a good illumination of the object. The transmitted light source and incident light source can be controlled independently.

NOTE!

Do not operate the microscope with maximum brightness over a longer time. This will reduce the LED lifetime.

The battery runtime is up to 30 hours (at minimum illumination power).

6. OBSERVATION

Place the desired object in the middle of the object plate (11) and secure it with the clamps (10) if necessary. The objective's magnification (2x or 4x) can be adjusted by turning the objective holder (7). It is recommended to start observations always with the 2x objective. Now look through the eyepieces (1) at the binocular head (4). Adjust the sharpness of the image with the focusing wheel (6).

The glass plate can be replaced by a black/white plate. Depending on the colour, the image contrast can be enhanced.

7. FOCUSING AND DIOPTRIC COMPENSATION

Look at an object. Close your left eye and adjust the focus with the focusing wheel (6) for your right eye. Close your right eye and adjust the focus for your left eye with the diopter setting (3). Now the microscope is adjusted for your eyes.

8. INTERPUPILLAR DISTANCE

This microscope is equipped with an adjustable binocular head. To set up the microscope to your individual eye distance, look through the eyepieces. Move the eyepieces against each other until you find the ideal position.

9. RUBBER EYECUPS

The rubber eyecups (2) can be used to minimize the entry of disturbing sidelight. For this put the rubber eyecups on the eyepieces. Turn the eyepieces so that the higher edges of the eyepieces are pointing outwards.

! NOTE!

The LED illumination has a lifespan of approx. 20,000 hours. In the case of a technical error please contact the service center.

10. SMARTPHONE HOLDER (FIG. 5)

Put the Smartphone holder (15) over the eyepiece and fix it with the knurled screw.

When using a device with binocular head, install the holder to the left or right eyepiece.

! NOTE!

The suction cups must be free of dust and dirt. It may help to slightly dampen them.

Make sure that the Smartphone holder is mounted securely and the Smartphone is firmly fixed on the plate.

Smartphones with a rough surface can not be fixed as good as that one with flat surfaces. Make sure that it will not slip down.

Start the camera App on your Smartphone. The camera lens must be placed directly over the eyepiece to get the image centered in the middle of the LCD.

It can be necessary to enlarge the image to the full LCD size by using the camera zoom. A shadowing at the edges is possible.

Remove the Smartphone from the holder after use.

11. CLEANING AND MAINTENANCE

Your microscope is a high-quality optical instrument. Please take care that neither dust nor fingerprints take place on the optical surfaces. If necessary, clean the microscope body with a slightly moistened soft cloth. Use a separate clean cloth with some drops of alcohol to clean the optical surfaces. After using, the microscope should be protected with its dust cover.

Remove the batteries from the unit if it is not used for a longer time.

Remember:

A well maintained microscope will keep its optical quality for years and thus maintain its value.

12. DISPOSAL

 Dispose of the packaging materials properly, according to their type, such as paper or cardboard. Please take the current legal regulations into account when disposing of your device. You can get more information on the proper disposal from your local waste-disposal service or environmental authority.

 Do not dispose of electronic devices in the household garbage! As per Directive 2002/96/EC of the European Parliament on waste electrical and electronic equipment and its adaptation into German law, used electronic devices must be collected separately and recycled in an environmentally friendly manner.

 In accordance with the regulations concerning batteries and rechargeable batteries, disposing of them in the normal household waste is explicitly forbidden. Please make sure to dispose of your used batteries as required by law – at a local collection point or in the retail market. Disposal in domestic waste violates the Battery Directive. Batteries that contain toxins are marked with a sign and a chemical symbol. „Cd“ = cadmium, „Hg“ = mercury, „Pb“ = lead.

13. EC DECLARATION OF CONFORMITY

Product type: Microscope with transmitted and incident light

Product description: Erudit ICD 20x-40x

Art. No.: 5803600

Bresser GmbH has issued a „Declaration of Conformity“ in accordance with applicable guidelines and corresponding standards. This can be viewed any time upon request.

Bresser GmbH

DE-46414 Rhede/Westf. · Germany

14. TECHNICAL DATA

	Modell Erudit ICD (Art. No. 5803600)
Microscope head	binocular
Eyepiece(s)	1 piece DIN WF 10x
Objectives	2x / 4x
Magnification	20x / 40x
Power supply	4.5V (3 batteries, AA/LR6 type)
Illumination	LED* (battery-powered)

* The LED illumination used in this device has a very long lifespan and does not need to be changed. It can not be changed due to technical reasons!

15. WARRANTY & SERVICE

The regular guarantee period is 2 years and begins on the day of purchase. To benefit from an extended voluntary guarantee period as stated on the gift box, registration on our website is required.

You can consult the full guarantee terms as well as information on extending the guarantee period and details of our services at www.bresser.de/warranty_terms.

Would you like detailed instructions for this product in a particular language? Then visit our website via the link below (QR code) for available versions.

Alternatively you can also send an email to manuals@bresser.de or leave a message on +49 (0) 28 72 – 80 74-220*. Please always state your name, precise address, a valid phone number and email address, as well as the article number and name.

*Number charged at local rates in Germany (the amount you will be charged per phone call will depend on the tariff of your phone provider); calls from abroad will involve higher costs.

www.bresser.de/download/5803600

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Irrefiner und technische Änderungen vorbehalten. · Errors and technical changes reserved.
Verzigtungen en technische veranderingen voorbehouden. · Sous réserve d'erreurs et de modifications techniques.
Queda reservada la posibilidad de incluir modificaciones o de que el texto contenga errores.
Con riserva di errori e modifiche tecniche. · Остативем за собой право на ошибки и технические изменения.
Manual_5803600_ErudiNCI_de-en_BRESSER_v01817a

Bresser GmbH

Gutenbergstr. 2 · DE-46414 Rhede
Germany

www.bresser.de · service@bresser.de