



**KERN & Sohn GmbH**

Ziegelei 1  
72336 Balingen-Frommern  
Germany

+0049-[0]7433-9933-0  
+0049-[0]7433-9933-149  
info@kern-sohn.com

# Installation Instructions

## Interface Adapter with Cable for USB\_Device/Slave

### KERN KUP-03

Type TYKUP-03-A  
Version 1.1  
2023-06  
GB



You will find the current version of these instructions also online under:  
<https://www.kern-sohn.com/shop/de/DOWNLOADS/>  
Under the column Operating instructions

TYKUP-03-A-IA-e-2311



# Interface Adapter with Cable for USB\_Device/ Slave

Version 1.1 2023-06

## Installation Instructions

### Contents

1	<b>General hints</b> .....	2
1.1	Installation.....	2

## 1 General hints

<b>i</b>	<ul style="list-style-type: none"><li>• <b>Only KERN KUP-adapters may be connected to the 15-pol-sub-D-connection of the balance!</b></li><li>• Use the KUP adapter to transfer data from the device to a PC and vice-versa.</li></ul>
----------	--

### 1.1 Installation

1. Switch off the appliance
2. Plug in the interface adapter at the KUP connection
3. Switch on the appliance. The interface adapter is automatically recognized by the appliance.

#### Communication parameters:

Communication parameters (baud rate, bits and parity) of balance and printer / PC must match.

Interface: USB (factory setting)	
Baud rate	9600
Data bits	8
Stop bits	1
Parity	none

## USB driver:

The USB driver required for the scale is already pre-installed on devices with an up-to-date operating system.

If the system is not recognized automatically, you can obtain the correct driver from the Downloads area at [www.kern-sohn.com](http://www.kern-sohn.com).

It will be recognized as a virtual COM port and will be assigned a corresponding numbered connection.

Example (Microsoft Windows Device Manager):

